USER GUIDE

Chromium Next GEM Single Cell 3' Reagent Kits v3.1

with Feature Barcoding technology for CRISPR Screening



FOR USE WITH

Chromium Next GEM Single Cell 3' GEM, Library & Gel Bead Kit v3.1, 16 rxns PN-1000121 Chromium Next GEM Single Cell 3' GEM, Library & Gel Bead Kit v3.1, 4 rxns PN-1000128 Chromium Single Cell 3' Feature Barcode Library Kit, 16 rxns PN-1000079 Chromium Next GEM Chip G Single Cell Kit, 48 rxns PN-1000120 Chromium Next GEM Chip G Single Cell Kit, 16 rxns PN-1000127 Single Index Kit T Set A, 96 rxns PN-1000213



Notices

Document Number

CG000205 • Rev D

Legal Notices

© 2019 10X Genomics, Inc (10x Genomics). All rights reserved. Duplication and/or reproduction of all or any portion of this document without the express written consent of 10x Genomics, is strictly forbidden. Nothing contained herein shall constitute any warranty, express or implied, as to the performance of any products described herein. Any and all warranties applicable to any products are set forth in the applicable terms and conditions of sale accompanying the purchase of such product. 10x Genomics provides no warranty and hereby disclaims any and all warranties as to the use of any third-party products or protocols described herein. The use of products described herein is subject to certain restrictions as set forth in the applicable terms and conditions of sale accompanying the purchase of such product. A non-exhaustive list of 10x Genomics' marks, many of which are registered in the United States and other countries can be viewed at: www.10xgenomics.com/trademarks. 10x Genomics may refer to the products or services offered by other companies by their brand name or company name solely for clarity, and does not claim any rights in those third party marks or names. 10x Genomics products may be covered by one or more of the patents as indicated at: www.10xgenomics.com/patents. The use of products described herein is subject to 10x Genomics Terms and Conditions of Sale, available at www.10xgenomics.com/legal-notices, or such other terms that have been agreed to in writing between 10x Genomics and user. All products and services described herein are intended FOR RESEARCH USE ONLY and NOT FOR USE IN DIAGNOSTIC PROCEDURES.

Instrument & Licensed Software Updates Warranties

Updates to existing Instruments and Licensed Software may be required to enable customers to use new or existing products. In the event of an Instrument failure resulting from an update, such failed Instrument will be replaced or repaired in accordance with the 10x Limited Warranty, Assurance Plan or service agreement, only if such Instrument is covered by any of the foregoing at the time of such failure. Instruments not covered under a current 10x Limited Warranty, Assurance Plan or service agreement will not be replaced or repaired.

Support

Email: support@10xgenomics.com

10x Genomics

6230 Stoneridge Mall Road Pleasanton, CA 94588 USA

Document Revision **Summary**

Document Number CG000205

Chromium Next GEM Single Cell 3' Reagent Kits v3.1

Title User Guide with Feature Barcoding technology for CRISPR

Screening

Revision Rev C to Rev D

Revision Date November 2019

Specific Changes:

• Updated index plate kit name and associated part numbers.

General Changes:

• Updated for general minor consistency of language and terms throughout.

Table of Contents

Introduction	6
Chromium Next GEM Single Cell 3' Reagent Kits v3.1	7
Chromium Accessories	11
Recommended Thermal Cyclers	11
Additional Kits, Reagents & Equipment	12
Protocol Steps & Timing	14
Stepwise Objectives	15
CRISPR Screening Overview	19
Tips & Best Practices	20
Step 1	26
GEM Generation & Barcoding	27
1.1 Prepare Master Mix	28
1.2 Load Chromium Next GEM Chip G	30
1.3 Run the Chromium Controller	31
1.4 Transfer GEMs	31
1.5 GEM-RT Incubation	32
Step 2	33
Post GEM-RT Cleanup & cDNA Amplification	34
2.1 Post GEM-RT Cleanup — Dynabeads	35
2.2 cDNA Amplification	37
2.3 cDNA Cleanup – SPRIselect	39
2.3A Pellet Cleanup	39
2.3B Transferred Supernatant Cleanup	39
2.4 Post cDNA Amplification QC & Quantification	40
Step 3	41
3' Gene Expression Library Construction	42
3.1 Fragmentation, End Repair & A-tailing	44
3.2 Post Fragmentation, End Repair & A-tailing Double Sided Size Selection – SPRIselect	45
3.3 Adaptor Ligation	46
3.4 Post Ligation Cleanup – SPRIselect	47
3.5 Sample Index PCR	48
3.6 Post Sample Index PCR Double Sided Size Selection – SPRIselect	49
3.7 Post Library Construction OC	50

Step 4	51
CRISPR Screening Library Construction	52
4.1 Guide RNA cDNA Cleanup – SPRIselect	53
4.2 Feature PCR	54
4.3 Post Feature PCR Cleanup – SPRIselect	55
4.4 Sample Index PCR	56
4.5 Post Sample Index PCR Double Sided Size Selection – SPRIselect	57
4.6 Post Library Construction QC	58
Sequencing	59
Troubleshooting	62
GEMs	63
Chromium Controller Errors	65
Appendix	66
Post Library Construction Quantification	67
Agilent TapeStation Traces	68
Compatible sgRNA Specifications	69
Oligonucleotide Sequences	70

Introduction

Chromium Next GEM Single Cell 3' Reagent Kits v3.1 Chromium Accessories Recommended Thermal Cyclers Additional Kits, Reagents & Equipment Protocol Steps & Timing Stepwise Objectives CRISPR Screening Overview

Chromium Next GEM Single Cell 3' Reagent Kits v3.1

Chromium Next GEM Single Cell 3' GEM, Library & Gel Bead Kit v3.1, 16 rxns PN-1000121

Chromium Next GEM Single Cell 3' GEM Kit v3.1 16 rxns PN-1000123 (store at -20°C)

Chromium Next GEM Single Cell 3'		
GEM Kit v3.1	#	PN
RT Reagent B	1	2000165
RT Enzyme C	1	2000085
Template Switch Oligo	1	3000228
Reducing Agent B	1	2000087
Cleanup Buffer	2	2000088
cDNA Primers	1	2000089
O Amp Mix	1	2000047
10xGenomics.com		10x GENOMICS

Chromium Next GEM Single Cell 3' Library Kit v3.1 16 rxns PN-1000157 (store at -20°C)

Single Cell 3' Library Kit v3.1	#	PN
Fragmentation Enzyme	1	2000090
Fragmentation Buffer	1	2000091
 Ligation Buffer 	1	2000092
DNA Ligase	1	220110
Adaptor Oligos	1	2000094
SI Primer	1	2000095
O Amp Mix	1	2000047

Chromium Next GEM Single Cell 3' Gel Bead Kit v3.1, 16 rxns PN-1000122 (store at -80°C)



Dynabeads[™] MyOne[™] SILANE PN-2000048 (store at 4°C)


Dynabeads MyOne 1 2000048 SILANE		#	PN
	Dynabeads MyOne SILANE	1	2000048

Chromium Next GEM Single Cell 3' GEM, Library & Gel Bead Kit v3.1, 4 rxns PN-1000128

Chromium Next GEM Single Cell 3'GEM Kit v3.1 4 rxns PN-1000130 (store at -20°C)

Chromium Next GEM Single Cell 3' GEM Kit v3.1 PΝ RT Reagent B 1 2000165 RT Enzyme C 1 2000102 Template Switch Oligo 1 3000228 Reducing Agent B 2000087 1 2000088 Cleanup Buffer cDNA Primers 1 2000089 O Amp Mix 1 2000103 10xGenomics.com 10x

Chromium Next GEM Single Cell 3' Library Kit v3.1 4 rxns PN-1000158 (store at -20°C)

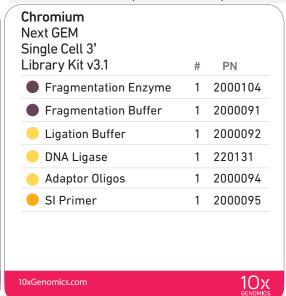

Chromium Next GEM Single Cell 3' Gel Bead Kit v3.1, 4 rxns PN-1000129 (store at -80°C)



Dynabeads[™] MyOne[™] SILANE PN-2000048 (store at 4°C)



Chromium Single Cell 3' Feature Barcode Library Kit, 16 rxns PN-1000079 (store at -20°C)



Chromium Next GEM Chip G Single Cell Kit, 48 rxns PN-1000120 (store at ambient temperature)



Chromium Next GEM Chip G Single Cell Kit, 16 rxns PN-1000127 (store at ambient temperature)



Single Index Kit T Set A, 96 rxns PN-1000213 (store at -20°C)



Chromium Accessories

Product	Part Number (Kit)	Part Number (Item)
10x Vortex Adapter	120251	330002
10x Magnetic Separator	120250	230003
Chromium Next GEM Secondary Holder	1000195	3000332

Recommended Thermal Cyclers

Thermal cyclers used must support uniform heating of 100 μ l emulsion volumes.

Supplier	Description	Part Number
BioRad	C1000 Touch Thermal Cycler with 96-Deep Well Reaction Module	1851197
Eppendorf	MasterCycler Pro	North America 950030010 International 6321 000.019
Thermo Fisher Scientific	Veriti 96-Well Thermal Cycler	4375786

Additional Kits, Reagents & Equipment

The items in the table below have been validated by 10x Genomics and are highly recommended for the Single Cell 3' protocols. Substituting materials may adversely affect system performance. This list does not include standard laboratory equipment, such as water baths, centrifuges, vortex mixers, pH meters, freezers, etc.

Supplier	Description		Part Number (US)	
Plastics				
Eppendorf	PCR Tubes 0.2 ml 8-tube strips DNA LoBind Tubes, 1.5 ml DNA LoBind Tubes, 2.0 ml	Choose either Eppendorf, USA Scientific or	951010022 022431021 022431048	
USA Scientific	TempAssure PCR 8-tube strip	Thermo Fisher Scientific PCR	1402-4700	
Thermo Fisher Scientific	MicroAmp 8-Tube Strip, 0.2 ml MicroAmp 8 -Cap Strip, clear	8-tube strips.	N8010580 N8010535	
Rainin	Tips LTS 200UL Filter RT-L200FLR Tips LTS 1ML Filter RT-L1000FLR Tips LTS 20UL Filter RT-L10FLR	Tips LTS 1ML Filter RT-L1000FLR		
Kits & Reagents				
Thermo Fisher Scientific	Nuclease-free Water Low TE Buffer (10 mM Tris-HCl pH 8.0, 0.1 mM E	DTA)	AM9937 12090-015	
Millipore Sigma	Ethanol, Pure (200 Proof, anhydrous)		E7023-500ML	
Beckman Coulter	SPRIselect Reagent Kit	B23318		
Bio-Rad	10% Tween 20	1662404		
Ricca Chemical Company	Glycerin (glycerol), 50% (v/v) Aqueous Solution		3290-32	
Qiagen	Qiagen Buffer EB		19086	
Equipment				
VWR	Vortex Mixer		10153-838	
Thermo Fisher Scientific	MYFUGE 12 Mini Centrifuge (alternatively, use any equivalent mini centrifuge)		C1012	
Eppendorf	Eppendorf ThermoMixer C Eppendorf SmartBlock 1.5 ml, Thermoblock for 24 reaction vessel (alternatively, use a temperature-controlled Heat Block)		5382000023 5360000038	
Rainin	Pipet-Lite Multi Pipette L8-50XLS+ Pipet-Lite Multi Pipette L8-200XLS+ Pipet-Lite Multi Pipette L8-10XLS+ Pipet-Lite Multi Pipette L8-20XLS+ Pipet-Lite LTS Pipette L-2XLS+ Pipet-Lite LTS Pipette L-10XLS+ Pipet-Lite LTS Pipette L-20XLS+ Pipet-Lite LTS Pipette L-100XLS+ Pipet-Lite LTS Pipette L-200XLS+ Pipet-Lite LTS Pipette L-1000XLS+ Pipet-Lite LTS Pipette L-1000XLS+		17013804 17013805 17013802 17013803 17014393 17014388 17014392 17014384 17014391	

Additional Kits, Reagents & Equipment

The items in the table below have been validated by 10x Genomics and are highly recommended for the Single Cell 3' protocols. Substituting materials may adversely affect system performance. This list does not include standard laboratory equipment, such as water baths, centrifuges, vortex mixers, pH meters, freezers, etc.

Supplier	Description		Part Number (US)
Quantification & Quality Contro	pl		
Agilent	2100 Bioanalyzer Laptop Bundle High Sensitivity DNA Kit 4200 TapeStation	Choose Bioanalyzer,	G2943CA 5067-4626 G2991AA
Thermo Fisher Scientific	High Sensitivity D1000 ScreenTape/Reagents High Sensitivity D5000 ScreenTape/Reagents	TapeStation, or Qubit based on availability & preference.	5067-5592/ 5067-5593 5067-5584/ 5067-5585
Thermo Fisher Scientific	Qubit 4.0 Flourometer Qubit dsDNA HS Assay Kit		Q33226 Q32854
Advanced Analytical	Fragment Analyzer Automated CE System - 12 cap Fragment Analyzer Automated CE System - 48/96 cap High Sensitivity NGS Fragment Analysis Kit		FSv2-CE2F FSv2-CE10F DNF-474
KAPA Biosystems	KAPA Library Quantification Kit for Illumina Platforms		KK4824

Protocol Steps & Timing

Day	Steps	Timing Stop & Store
	Cell Preparation	
2 h	Dependent on Cell Type	~1-1.5 h
211	Step 1 – GEM Generation & Barcoding	
	 1.1 Prepare Reaction Mix 1.2 Load Chromium Next GEM Chip G 1.3 Run the Chromium Controller 1.4 Transfer GEMs 1.5 GEM-RT Incubation 	20 min 10 min 18 min 3 min 55 min 550 4°C ≤72 h or −20°C ≤1 week
4 h	Step 2 – Post GEM-RT Cleanup & cDNA Amplification	_
	 2.1 Post GEM RT-Cleanup – Dynabead 2.2 cDNA Amplification 2.3 cDNA Cleanup – SPRIselect 	45 min 40 min
	2.3A Pellet Cleanup 2.3B Transferred Supernatant Cleanup 2.4 cDNA QC & Quantification	20 min 4° C ≤72 h or −20°C ≤4 weeks 30 min 4° C ≤72 h or −20°C ≤4 weeks 50 min
6 h	Step 3 – 3' Gene Expression Library Construction	
	 3.1 Fragmentation, End Repair & A-tailing 3.2 Post Fragmentation, End Repair & A-tailing Double Sided Size Selection – SPRIselect 	50 min 30 min
	3.3 Adaptor Ligation3.4 Post Ligation Cleanup – SPRIselect	25 min 20 min
	3.5 Sample Index PCR	40 min 5 ⁵ 10 4°C ≤72 h
	3.6 Post Sample Index PCR Double Sided Size Selection – SPRIselect	30 min 4°C ≤72 h or -20°C long term
	3.7 Post Library Construction QC	50 min
8 h Plus	Step 4 – CRISPR Screening Library Construction	
	4.1 Guide RNA cDNA Cleanup	20 min
	4.2 Feature PCR	50 min
	4.3 Post Feature PCR Cleanup – SPRIselect	20 min
	4.4 Sample Index PCR	30 min
	4.5 Post Sample Index PCR Size Selection – SPRIselect4.6 Post Library Construction QC	20 min 4°C ≤72 h or -20°C long term 50 min

Stepwise Objectives




The Chromium Single Cell Gene Expression Solution with Feature Barcoding technology upgrades short read sequencers to deliver a scalable microfluidic platform for assessing CRISPR-mediated perturbations of gene expression by profiling 500-10,000 individual cells per sample. A pool of ~3,500,000 10x Barcodes to separately index each cell's transcriptome along with the CRISPR-mediated perturbations. It was done by partitioning thousands of cells into nanoliter-scale Gel Beads-in-emulsion (GEMs), where all generated cDNA (from poly-adenylated mRNAs and single-guide RNAs/sgRNAs) share a common 10x Barcode. Libraries are generated and sequenced from the cDNA and 10x Barcodes are used to associate individual reads back to the individual partitions.

This document outlines the protocol for generating Single Cell 3' Gene Expression and CRISPR Screening libraries from the same cells.

Single Cell 3' v3 Gel Beads


In addition to the poly(dT) primer that enables the production of barcoded, full-length cDNA from poly-adenylated mRNA, the Single Cell 3' v3.1 Gel Beads also include two additional primer sequences (Capture Sequence 1 and Capture Sequence 2), that enable capture and priming of Feature Barcoding technology compatible targets or analytes of interest.

The poly(dT) primers along with one of the capture sequence primers are used in this protocol for generating Single Cell 3' Gene Expression and CRISPR Screening libraries.

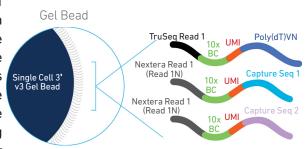

Step 1 **GEM Generation &** Barcoding

GEMs are generated by combining barcoded Single Cell 3' v3.1 Gel Beads, a Master Mix containing cells, and Partitioning Oil onto Chromium Next GEM Chip G. To achieve single cell resolution, cells are delivered at a limiting dilution, such that the majority (~90-99%) of generated GEMs contain no cell, while the remainder largely contain a single cell.

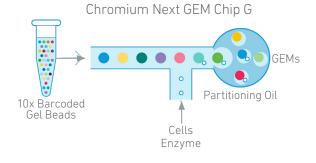

Step 1 **GEM Generation &** Barcoding

Immediately following **GEM** generation, the Gel Bead is dissolved releasing the three types of primers and any co-partitioned cell is lysed. The poly(dT) and the capture sequence primers in the gel bead are engaged simultaneously in two different reactions inside individual GEMs (primer with Capture Sequence 2 is not shown in the illustrated example).

A. Primers containing:

- an Illumina TruSeq Read 1 (read 1 sequencing primer)
- 16 nt 10x Barcode
- 12 nt unique molecular identifier (UMI)
- 30 nt poly(dT) sequence


are mixed with the cell lysate and a Master Mix containing reverse transcription (RT) reagents. Incubation of the GEMs produces barcoded, full-length cDNA from poly-adenylated mRNA.

B. In the same partition, primers containing:

- an Illumina Nextera Read 1 (Read 1N; read 1 sequencing primer)
- 16 nt 10x Barcode
- 12 nt unique molecular identifier (UMI)
- Capture Sequence 1

are mixed with the cell lysate and the Master Mix containing RT reagents. Incubation of the GEMs produces barcoded, cDNA from the sgRNA.


See Appendix for Compatible sqRNA Specifications

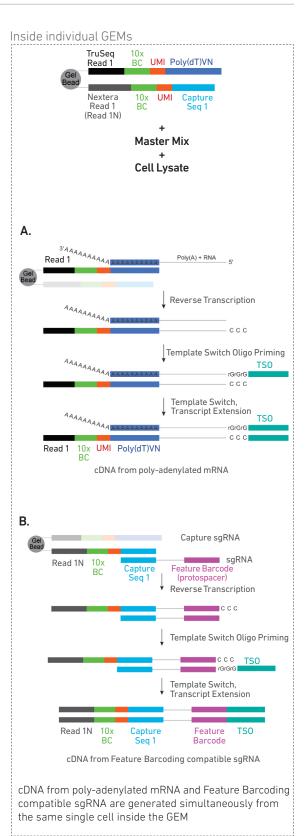

Step 2 Post GEM-RT Cleanup & cDNA Amplification




After incubation, GEMs are broken and pooled fractions are recovered. Silane magnetic beads are used to purify the cell barcoded products from the post GEM-RT reaction mixture, which includes leftover biochemical reagents and primers. The cell barcoded cDNA molecules are amplified via PCR to generate sufficient mass for library constructions. Size selection is used to separate the amplified cDNA molecules for 3' Gene Expression and and CRISPR Screening library construction.

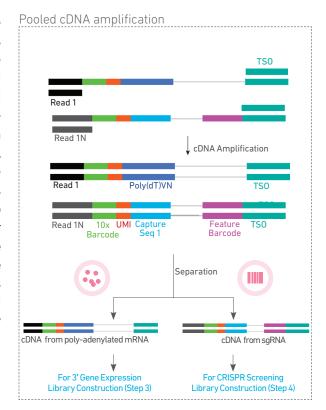

Step 3 3' Gene Expression **Library Construction**




cDNA Full-length from poly-adenylated mRNA is used to generate 3' Gene Expression libraries. Enzymatic fragmentation and size selection are used to optimize the cDNA amplicon size. TruSeq Read 1 (read 1 primer sequence) is added to the molecules during GEM incubation. P5, P7, a sample index, and TruSeq Read 2 (read 2 primer sequence) are added via End Repair, A-tailing, Adaptor Ligation, and PCR. The final libraries contain the P5 and P7 primers used in Illumina bridge amplification.

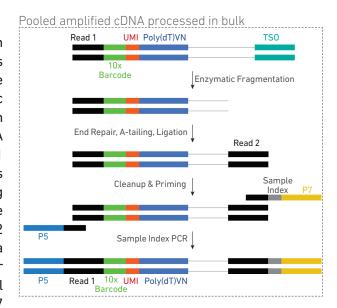

Step 4 **CRISPR Screening Library Construction**



Amplified cDNA from sgRNA molecules is used to generate CRISPR Screening libraries. P5, P7, a sample index, and TruSeq Read 2 (read 2 primer sequence) are added via PCR. The final libraries contain the P5 and P7 primers used in Illumina bridge amplification.

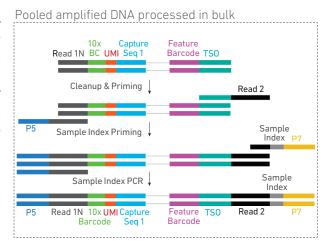


Step 5 Sequencing

The Single Cell 3' libraries comprise standard Illumina paired-end constructs which begin and end with P5 and P7. The 16 bp 10x Barcode and 12 bp UMI are encoded in Read 1, while Read 2 is used to sequence the cDNA fragment in 3' Gene Expression libraries and the Feature Barcode (sqRNA protospacer) in the CRISPR Screening libraries. Sample index sequences are incorporated as the sample index read. Standard Illumina sequencing primer sites TruSeg Read 1 and TruSeg Read 2 in the 3' Gene Expression libraries and Nextera Read 1 and Truseq Read 2 in the CRISPR Screening libraries are used in paired-end sequencing.

Illumina sequencer compatibility, sample indices, library loading and pooling, recommended read depths & run parameters for sequencing are summarized in step 5.

Chromium Single Cell 3' Gene Expression Library

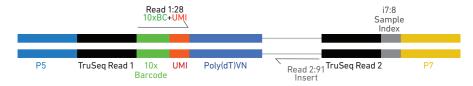

Chromium Single Cell 3' CRISPR Screening Library




*Minimum required Read 2 length for CRISPR Screening libraries is 70 bp

See Appendix for Oligonucleotide Sequences

CRISPR Screening Overview

The Chromium Single Cell Gene Expression Solution with Feature Barcoding technology provides a high-throughput and scalable approach to obtain gene expression profiles along with perturbation phenotypes via direct capture of poly-adentylated mRNAs and single-guide RNAs (sgRNAs) from the same single cell (see Stepwise Objectives). For compatibility with Feature Barcoding technology, sqRNAs should be engineered containing either Capture Sequence 1 or Capture Sequence 2. Two possible locations for integrating the capture sequence in the sgRNA include (1) within the sgRNA hairpin structure, or (2) immediately before the sgRNA termination signal, elongating the 3'-end of the sgRNA. However, alternate sgRNA integration locations for either of the two capture sequences may be possible depending on the specific application, type of construct used etc.

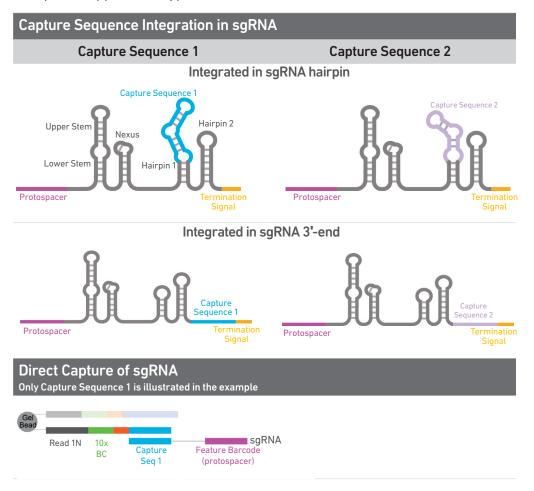



Performing sgRNA QC by qPCR, NGS, or other methods is recommended prior to proceeding with the Single Cell Gene Expression and CRISPR Screening Solution.

See Appendix for Compatible sgRNA Specifications

Tips & **Best Practices**



Icons









Emulsion-safe **Plastics**

• Use validated emulsion-safe plastic consumables when handling GEMs as some plastics can destabilize GEMs.

Cell Concentration

- Recommended starting point is to load ~1,600 cells per reaction, resulting in recovery of ~1000 cells, and a multiplet rate of ~0.8%. The optimal input cell concentration is $700-1,200 \text{ cells/}\mu\text{l}.$
- The presence of dead cells in the suspension may also reduce the recovery rate. Consult the 10x Genomics Single Cell Protocols Cell Preparation Guide and the Guidelines for Optimal Sample Preparation flowchart (Documents CG00053 and CG000126 respectively) for more information on preparing cells.

Multiplet Rate (%)	# of Cells Loaded	# of Cells Recovered
~0.4%	~800	~500
~0.8%	~1,600	~1,000
~1.6%	~3,200	~2,000
~2.3%	~4,800	~3,000
~3.1%	~6,400	~4,000
~3.9%	~8,000	~5,000
~4.6%	~9,600	~6,000
~5.4%	~11,200	~7,000
~6.1%	~12,800	~8,000
~6.9%	~14,400	~9,000
~7.6%	~16,000	~10,000

General Reagent Handling

- · Fully thaw and thoroughly mix reagents before use.
- Keep all enzymes and Master Mixes on ice during setup and use. Promptly move reagents back to the recommended storage.
- Calculate reagent volumes with 10% excess of 1 reaction values.
- Cover Partitioning Oil tubes and reservoirs to minimize evaporation.
- If using multiple chips, use separate reagent reservoirs for each chip during loading.
- Thoroughly mix samples with the beads during bead-based cleanup steps.

50% Glycerol Solution

- Purchase 50% glycerol solution from Ricca Chemical Company, Glycerin (glycerol), 50% (v/v) Aqueous Solution, PN-3290-32.
- Prepare 50% glycerol solution:
 - i. Mix an equal volume of water and 99% Glycerol, Molecular Biology Grade.
 - ii. Filter through a 0.2 µm filter.
 - iii. Store at –20°C in 1-ml LoBind tubes. 50% glycerol solution should be equilibrated to room temperature before use.

Pipette Calibration

- Follow manufacturer's calibration and maintenance schedules.
- Pipette accuracy is particularly important when using SPRIselect reagents.

Chromium **Next GEM Chip** Handling



- Minimize exposure of reagents, chips, and gaskets to sources of particles and fibers, laboratory wipes, frequently opened flip-cap tubes, clothing that sheds fibers, and dusty surfaces.
- After removing the chip from the sealed bag, use in ≤ 24 h.
- · Execute steps without pause or delay, unless indicated. When multiple chips are to be used, load, run, and collect the content from one chip before loading the next.
- Fill all unused input wells in rows labeled 1, 2, and 3 on a chip with an appropriate volume of 50% glycerol solution before loading the used wells. DO NOT add glycerol to the wells in the bottom NO FILL row.
- Avoid contacting the bottom surface of the chip with gloved hands and other surfaces. Frictional charging can lead to inadequate priming of the channels, potentially leading to either clogs or wetting failures.
- Minimize the distance that a loaded chip is moved to reach the Chromium Controller.
- Keep the chip horizontal to prevent wetting the gasket with oil, which depletes the input volume and may adversely affect the quality of the resulting emulsion.

Chromium **Next GEM** Secondary Holders







Chromium **Next GEM Chip** Loading





- The holder lid flips over to become a stand, holding the chip at 45 degrees for optimal recovery well content removal.
- Squeeze the black sliders on the back side of the holder together to unlock the lid and return the holder to a flat position.



- Align notch on the chip (upper left corner) and the holder.
- Insert the left-hand side of the chip under the guide. Depress the right-hand side of the Next GEM Chip chip until the spring-loaded clip engages.
- Close the lid before dispensing reagents into the wells.



- Place the assembled chip and holder flat on the bench with the lid closed.
- Dispense at the bottom of the wells without introducing bubbles.
- When dispensing Gel Beads into the chip, wait for the remainder to drain into the bottom of the pipette tips and dispense again to ensure complete transfer.
- Refer to Load Chromium Next GEM Chip G for specific instructions.



Gel Bead Handling



- Use one tube of Gel Beads per sample. DO NOT puncture the foil seals of tubes not used at the time.
- Equilibrate the Gel Beads strip to room temperature before use.
- Store unused Gel Beads at –80°C and avoid more than 12 freeze-thaw cycles. DO NOT store Gel Beads at -20°C.



- Snap the tube strip holder with the Gel Bead strip into a 10x Vortex Adapter. Vortex 30 sec.
- Centrifuge the Gel Bead strip for ~5 sec after removing from the holder. Confirm there are no bubbles at the bottom of the tubes and the liquid levels look even. Place the Gel Bead strip back in the holder and secure the holder lid.
- If the required volume of beads cannot be recovered, place the pipette tips against the sidewalls and slowly dispense the Gel Beads back into the tubes. DO NOT introduce bubbles into the tubes and verify that the pipette tips contain no leftover Gel Beads. Withdraw the full volume of beads again by pipetting slowly.

10x Gasket **Attachment**

- · After reagents are loaded, attach the gasket by holding the tongue (curved end, to the right) and hook it on the left-hand tabs of the holder. Gently pull the gasket Notched Cut toward the right and hook it on the two right-hand tabs.
- DO NOT touch the smooth side of the gasket. DO NOT press down on the top of the gasket after attachment.
- Keep the assembly horizontal to avoid wetting the gasket with Partitioning Oil.




10x Magnetic Separator

- Offers two positions of the magnets (high and low) relative to a tube, depending on its orientation. Flip the magnetic separator over to switch between high (magnet.High) or low (magnet.Low) positions.
- If using MicroAmp 8-Tube Strips, use the high position (magnet•High) only throughout the protocol.



Magnetic Bead Cleanup Steps

- During magnetic bead based cleanup steps that specify waiting "until the solution clears", visually confirm clearing of solution before proceeding to the next step. See adjacent panel for an example.
- The time needed for the solution to clear may vary based on specific step, reagents, volume of reagents etc.

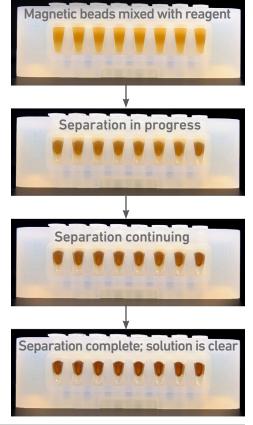


SPRIselect Cleanup & Size Selection

- · After aspirating the desired volume of SPRIselect reagent, examine the pipette tips before dispensing to ensure the correct volume is transferred.
- Pipette mix thoroughly as insufficient mixing of sample and SPRIselect reagent will lead to inconsistent results.
- Use fresh preparations of 80% Ethanol.

Tutorial — SPRIselect Reagent: DNA Sample Ratios

SPRI beads selectively bind DNA according to the ratio of SPRIselect reagent (beads).

Example: Ratio = Volume of SPRIselect reagent added to the sample = $50 \mu l$ = 0.5X Volume of DNA sample 100 µl

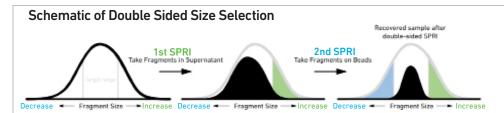

After the first SPRI, supernatant is transferred for a second SPRI while larger fragments are discarded (green). After the second SPRI, fragments on beads are eluted and kept while smaller fragments are discarded (blue). Final sample has a tight fragment size distribution with reduced overall amount (black).

Tutorial — Double Sided Size Selection

Step a – First SPRIselect: Add 50 μl SPRIselect reagent to 100 μl sample (0.5X).

Ratio = Volume of SPRIselect reagent added to the sample = $50 \mu l$ = 0.5X Volume of DNA sample

Step b – Second SPRIselect: Add 30 µl SPRIselect reagent to supernatant from step a (0.8X).

Ratio = $\underline{\text{Total Volume of SPRIselect reagent added to the sample (step a + b)}}$ = $\underline{50 \, \mu l + 30 \, \mu l}$ = 0.8X Original Volume of DNA sample 100 µl

Enzymatic Fragmentation

 Ensure enzymatic fragmentation reactions are prepared on ice and then loaded into a thermal cycler pre-cooled to 4°C prior to initiating the Fragmentation, End Repair, and A-tailing incubation steps.

Sample Indices in Sample Index PCR

- Choose the appropriate sample index sets to ensure that no sample indices overlap in a multiplexed sequencing run.
- Each well in the Single Index plate contains a unique mix of 4 oligos.
- The sample indexes can therefore be used in any combination.
- · Each sample index set is base-balanced to avoid monochromatic signal issues when it is the sole sample loaded on an Illumina sequencer.

Step 1

GEM Generation & Barcoding

- 1.1 Prepare Single Cell Master Mix
- 1.2 Load Chromium Next GEM Chip G
- 1.3 Run the Chromium Controller
- 1.4 Transfer GEMs
- 1.5 **GEM-RT Incubation**

1.0 **GEM Generation &** Barcoding



GET STARTED!						
Action		Item	10x PN	Preparation & Handling	Storage	
Equilibrate to Room Temperature		Single Cell 3' v3.1 Gel Beads	2000164	Equilibrate to room temperature 30 min before loading the chip.	-80°C	
	•	RT Reagent B	2000165	Vortex, verify no precipitate, centrifuge briefly.	−20°C	
		Template Switch Oligo	3000228	Centrifuge briefly, resuspend in 80 µl Low TE Buffer. Vortex 15 sec at maximum speed, centrifuge briefly, leave at room temperature for ≥ 30 min. After resuspension, store at -80°C.	-20°C	
	0	Reducing Agent B	2000087	Vortex, verify no precipitate, centrifuge briefly.	-20°C	
Place on Ice		RT Enzyme C	2000085/ 2000102	Centrifuge briefly before adding to the mix.	-20°C	
		Cell Suspension				
Obtain		Partitioning Oil	2000190	-	Ambient	
		Chromium Next GEM Chip G	2000177	-	Ambient	
		10x Gasket	370017/ 3000072	See Tips & Best Practices.	Ambient	
		Chromium Next GEM Secondary Holder	3000332	See Tips & Best Practices.	Ambient	
		10x Vortex Adapter	330002	See Tips & Best Practices.	Ambient	
		50% glycerol solution If using <8 reactions	-	See Tips & Best Practices.	-	



Firmware Version 4.0 or higher is required in the Chromium Controller or the Chromium Single Cell Controller used for this Single Cell 3' v3.1 protocol.

1.1 Prepare Master Mix



a. Prepare Master Mix on ice. Pipette mix 15x and centrifuge briefly.

Master Mix Add reagents in the order listed	PN	1X (µl)	4Χ + 10% (μl)	8X + 10% (µl)
RT Reagent B	2000165	18.8	82.2	165.0
Template Switch Oligo	3000228	2.4	10.4	20.8
Reducing Agent B	2000087	2.0	8.6	17.3
RT Enzyme C	2000085/ 2000102	8.7	38.4	76.8
Total	-	31.8	139.9	279.8

b. Add 31.8 µl Master Mix into each tube of a PCR 8-tube strip on ice.


Assemble Chromium Next GEM Chip G



After removing the chip from the sealed bag, use the chip in ≤ 24 h.



See Tips & Best Practices for chip handling instructions.

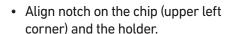

- · Insert the left-hand side of the chip under the guide. Depress the righthand side of the chip until the springloaded clip engages.
- · Close the lid before dispensing reagents into the wells.
- The assembled chip is ready for loading the indicated reagents. Refer to step 1.2 for reagent volumes and loading order.



Chromium Next GEM Chip G

Notch



For GEM generation, load the indicated reagents only in the specified rows, starting from row labeled 1, followed by rows labeled 2 and 3. DO NOT load reagents in the bottom row labeled NO FILL. See step 1.2 for details.





Cell Suspension Volume Calculator Table

(for step 1.2 of Chromium Next GEM Single Cell 3' v3.1 protocol)

Volume of Cell Suspension Stock per reaction (µl) | Volume of Nuclease-free Water per reaction (µl)

Cell Stock	Targeted Cell Recovery										
Concentration (Cells/µl)	500	1000	2000	3000	4000	5000	6000	7000	8000	9000	10000
100	8.3 35.0	16.5 26.7	33.0 10.2	n/a	n/a	n/a	n/a	n/a	n/a	n/a	n/a
200	4.1 39.1	8.3 35.0	16.5 26.7	24.8 18.5	33.0 10.2	41.3 2.0	n/a	n/a	n/a	n/a	n/a
300	2.8 40.5	5.5 37.7	11.0 32.2	16.5 26.7	22.0 21.2	27.5 15.7	33.0 10.2	38.5 4.7	n/a	n/a	n/a
400	2.1	4.1	8.3	12.4	16.5	20.6	24.8	28.9	33.0	37.1	41.3
	41.1	39.1	35.0	30.8	26.7	22.6	18.5	14.3	10.2	6.1	2.0
500	1.7	3.3	6.6	9.9	13.2	16.5	19.8	23.1	26.4	29.7	33.0
	41.6	39.9	36.6	33.3	30.0	26.7	23.4	20.1	16.8	13.5	10.2
600	1.4	2.8	5.5	8.3	11.0	13.8	16.5	19.3	22.0	24.8	27.5
	41.8	40.5	37.7	35.0	32.2	29.5	26.7	24.0	21.2	18.5	15.7
700	1.2 42.0	2.4 40.8	4.7 38.5	7.1 36.1	9.4 33.8	11.8 31.4	14.1 29.1	16.5 26.7	18.9 24.3	21.2 22.0	23.6 19.6
800	1.0	2.1	4.1	6.2	8.3	10.3	12.4	14.4	16.5	18.6	20.6
	42.2	41.1	39.1	37.0	35.0	32.9	30.8	28.8	26.7	24.6	22.6
900	0.9	1.8	3.7	5.5	7.3	9.2	11.0	12.8	14.7	16.5	18.3
	42.3	41.4	39.5	37.7	35.9	34.0	32.2	30.4	28.5	26.7	24.9
1000	0.8	1.7	3.3	5.0	6.6	8.3	9.9	11.6	13.2	14.9	16.5
	42.4	41.6	39.9	38.3	36.6	35.0	33.3	31.7	30.0	28.4	26.7
1100	0.8	1.5	3.0	4.5	6.0	7.5	9.0	10.5	12.0	13.5	15.0
	42.5	41.7	40.2	38.7	37.2	35.7	34.2	32.7	31.2	29.7	28.2
1200	0.7	1.4	2.8	4.1	5.5	6.9	8.3	9.6	11.0	12.4	13.8
	42.5	41.8	40.5	39.1	37.7	36.3	35.0	33.6	32.2	30.8	29.5
1300	0.6	1.3	2.5	3.8	5.1	6.3	7.6	8.9	10.2	11.4	12.7
	42.6	41.9	40.7	39.4	38.1	36.9	35.6	34.3	33.0	31.8	30.5
1400	0.6	1.2	2.4	3.5	4.7	5.9	7.1	8.3	9.4	10.6	11.8
	42.6	42.0	40.8	39.7	38.5	37.3	36.1	35.0	33.8	32.6	31.4
1500	0.6	1.1	2.2	3.3	4.4	5.5	6.6	7.7	8.8	9.9	11.0
	42.7	42.1	41.0	39.9	38.8	37.7	36.6	35.5	34.4	33.3	32.2
1600	0.5	1.0	2.1	3.1	4.1	5.2	6.2	7.2	8.3	9.3	10.3
	42.7	42.2	41.1	40.1	39.1	38.0	37.0	36.0	35.0	33.9	32.9
1700	0.5	1.0	1.9	2.9	3.9	4.9	5.8	6.8	7.8	8.7	9.7
	42.7	42.2	41.3	40.3	39.3	38.3	37.4	36.4	35.4	34.5	33.5
1800	0.5	0.9	1.8	2.8	3.7	4.6	5.5	6.4	7.3	8.3	9.2
	42.7	42.3	41.4	40.5	39.5	38.6	37.7	36.8	35.9	35.0	34.0
1900	0.4	0.9	1.7	2.6	3.5	4.3	5.2	6.1	6.9	7.8	8.7
	42.8	42.3	41.5	40.6	39.7	38.9	38.0	37.1	36.3	35.4	34.5
2000	0.4	0.8	1.7	2.5	3.3	4.1	5.0	5.8	6.6	7.4	8.3
	42.8	42.4	41.6	40.7	39.9	39.1	38.3	37.4	36.6	35.8	35.0

Grey boxes: Yellow boxes:

Volumes that would exceed the allowable water volume in each reaction

Indicate a low transfer volume that may result in higher cell load variability

Blue boxes: Optimal range of cell stock concentration to maximize the likelihood of achieving the desired cell recovery target

1.2 **Load Chromium** Next GEM Chip G



After removing the chip from the sealed bag, use in ≤ 24 h. When loading the chip, raising and depressing the pipette plunger should each take ~5 sec. When dispensing, raise the pipette tips at the same rate as the liquid is rising, keeping the tips slightly submerged.

a. Dispense 50% Glycerol into Unused Chip Wells (if < 8 samples per chip)

i. 70 ul to unused wells in row labeled 1.

ii. 50 μ l to unused wells in row labeled 2.

iii. 45 µl to unused wells in row labeled 3.

DO NOT add 50% glycerol solution to the bottom row of NO FILL wells. DO NOT use any substitute for 50% glycerol solution.

b. Prepare Master Mix + Cell Suspension

Refer to the Cell Suspension Volume Calculator Table. Add the appropriate volume of nucleasefree water and corresponding volume of single cell suspension to Master Mix for a total of 75 µl in each tube. Gently pipette mix the cells suspension before adding to the Master Mix.

c. Load Row Labeled 1

Gently pipette mix the Master Mix + Cell Suspension and using the same pipette tip, dispense 70 µl Master Mix + Cell Suspension into the bottom center of each well in row labeled 1 without introducing bubbles.





d. Prepare Gel Beads

Snap the tube strip holder with the Gel Bead strip into a 10x Vortex Adapter. Vortex 30 sec. Centrifuge the Gel Bead strip for ~5 sec. Confirm there are no bubbles at the bottom of the tubes and the liquid levels are even. Place the Gel Bead strip back in the holder. Secure the holder lid.



e. Load Row Labeled 2

Puncture the foil seal of the Gel Bead tubes. Slowly aspirate 50 µl Gel Beads. Dispense into the wells in row labeled 2 without introducing bubbles. Wait 30 sec.





f. Load Row Labeled 3

Dispense 45 µl Partitioning Oil into the wells in row labeled 3 from a reagent reservoir. Failure to add Partitioning Oil to the top row labeled 3 will prevent GEM generation and can damage the Chromium Controller.





Attach the gasket and run the chip in the Chromium Controller immediately after loading the Partitioning Oil.

q. Attach 10x Gasket

Align the notch with the top left-hand corner. Ensure the gasket holes are aligned with the wells. Avoid touching the smooth surface.



Keep horizontal to avoid wetting the gasket. DO NOT press down on the gasket.

1.3 Run the Chromium Controller



- a. Press the eject button on the Controller to eject the tray.
- **b.** Place the assembled chip with the gasket in the tray, ensuring that the chip stays horizontal. Press the button to retract the tray.
- c. Confirm the Chromium Chip G program on screen. Press the play button.



d. At completion of the run (~18 min), the Controller will chime. Immediately proceed to the next step.



Firmware Version 4.0 or higher is required in the Chromium Controller or the Chromium Single Cell Controller used for the Single Cell 3' v3.1 protocol.





1.4 Transfer GEMs



- a. Place a tube strip on ice.
- b. Press the eject button of the Controller and remove the chip.
- c. Discard the gasket. Open the chip holder. Fold the lid back until it clicks to expose the wells at 45 degrees.



- d. Check the volume in rows labeled 1-2. Abnormally high volume in any well indicates a clog.
- e. Slowly aspirate 100 µl GEMs from the lowest points of the recovery wells in the top row labeled 3 without creating a seal between the tips and the bottom of the wells.




- f. Withdraw pipette tips from the wells. GEMs should appear opaque and uniform across all channels. Excess Partitioning Oil (clear) in the pipette tips indicates a potential clog.
- g. Over the course of ~20 sec, dispense GEMs into the tube strip on ice with the pipette tips against the sidewalls of the tubes.
- h. If multiple chips are run back-to-back, cap/ cover the GEM-containing tube strip and place on ice for no more than 1 h.

Expose Wells at 45 Degrees





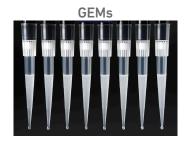

1.5 **GEM-RT Incubation**

Use a thermal cycler that can accommodate at least 100 μ l volume. A volume of $125 \,\mu l$ is the preferred setting on the Bio-Rad C1000 Touch. In alternate thermal cyclers, use highest reaction volume setting.

a. Incubate in a thermal cycler with the following protocol.

Lid Temperature	Reaction Volume	Run Time
53°C	125 µl	~55 min
Step	Temperature	Time
1	53°C	00:45:00
2	85°C	00:05:00
3	4°C	Hold



b. Store at 4°C for up to 72 h or at -20°C for up to a week, or proceed to the next step.

Step 2

Post GEM–RT Cleanup & cDNA Amplification

- Post GEM-RT Cleanup Dynabeads 2.1
- 2.2 cDNA Amplification
- 2.3 cDNA Cleanup – SPRIselect
- 2.4 cDNA QC & Quantification

2.0 Post GEM-RT Cleanup & cDNA Amplification



GET STARTE	D!			
Action	Item	10x PN	Preparation & Handling	Storage
Equilibrate to Room Temperature	Reducing Agent B	2000087	Thaw, vortex, verify no precipitate, centrifuge.	–20°C
	Feature cDNA Primers 1 Verify name & PN	2000096	Vortex, centrifuge briefly.	-20°C
	Dynabeads MyOne SILANE	2000048	Vortex thoroughly (≥30 sec) immediately before adding to the mix.	4°C
	Beckman Coulter SPRIselect Reagent	-	Manufacturer's recommendations.	-
	Agilent Bioanalyzer High Sensitivity Kit If used for QC and quantification	-	Manufacturer's recommendations.	-
	Agilent TapeStation ScreenTape and Reagents If used for QC and quantification	-	Manufacturer's recommendations.	-
	Qubit dsDNA HS Assay Kit If used for QC and quantification	-	Manufacturer's recommendations.	-
Place on ice	○ Amp Mix	2000047/ 2000103	Vortex, centrifuge briefly.	-20°C
Thaw at 65°C	Cleanup Buffer	2000088	Thaw for 10 min at 65°C at max speed on a thermomixer. Verify no visible crystals. Cool to room temperature.	-20°C
Obtain	Recovery Agent	220016	-	Ambient
	Qiagen Buffer EB	-	Manufacturer's recommendations.	-
	Bio-Rad 10% Tween 20	-	Manufacturer's recommendations.	-
	10x Magnetic Separator	230003	-	Ambient
	Prepare 80% Ethanol Prepare 15 ml for 8 reactions.	-	-	-

2.1 Post GEM-RT Cleanup -**Dynabeads**



a. Add 125 µl Recovery Agent to each sample at room temperature. DO NOT pipette mix or vortex the biphasic mixture. Wait 2 min.

The resulting biphasic mixture contains Recovery Agent/Partitioning Oil (pink) and agueous phase (clear), with no persisting emulsion (opaque).



If biphasic separation is incomplete:

Firmly secure the cap on the tube strip, ensuring that no liquid is trapped between the cap and the tube rim. Mix by inverting the capped tube strip 5x, centrifuge briefly, and proceed to step b. DO NOT invert without firmly securing the caps.



A smaller aqueous phase volume indicates a clog during GEM generation.



b. Slowly remove and discard 125 µl Recovery Agent/Partitioning Oil (pink) from the bottom of the tube. DO NOT aspirate any aqueous sample.

c. Prepare Dynabeads Cleanup Mix.



	Dynabeads Cleanup Mix Add reagents in the order listed	PN	1X (µl)	4X + 10% (µl)	8X + 10% (μl)
	Cleanup Buffer	2000088	182	801	1602
Resuspend clump	Dynabeads MyOne SILANE Vortex thoroughly (≥30 sec) immediately before adding to the mix. Aspirate the full liquid volume with a pipette tip to verify that the beads have not settled in the bottom of the tube. If clumps are present, pipette mix to resuspend completely. DO NOT centrifuge before use.	2000048	8	35	70
	Reducing Agent B	2000087	5	22	44
	Nuclease-free Water		5	22	44
	Total	-	200	880	1760



- d. Vortex and add 200 µl to each sample. Pipette mix 10x (pipette set to 200 μl).
- e. Incubate 10 min at room temperature. Pipette mix again at ~5 min after start of incubation to resuspend settled beads.

Add Dynabeads Cleanup Mix



f. Prepare Elution Solution I. Vortex and centrifuge briefly.

Elution Solution I Add reagents in the order listed	PN	1X (μl)	10Χ (μl)
Buffer EB	-	98	980
10% Tween 20	-	1	10
Reducing Agent B	2000084	1	10
Total	-	100	1000



g. At the end of 10 min incubation, place on a 10x Magnetic Separator•High position (magnet•High) until the solution clears.

A white interface between the aqueous phase and Recovery Agent is normal.

- h. Remove the supernatant.
- i. Add $300 \mu l$ 80% ethanol to the pellet while on the magnet. Wait 30 sec.
- i. Remove the ethanol.
- k. Add 200 µl 80% ethanol to pellet. Wait 30 sec.
- **l.** Remove the ethanol.
- m.Centrifuge briefly. Place on the magnet•Low.
- n. Remove remaining ethanol. Air dry for 1 min.
- o. Remove from the magnet. Immediately add 35.5 µl Elution Solution I.
- p. Pipette mix (pipette set to 30 μl) without introducing bubbles.
- q. Incubate 2 min at room temperature.
- r. Place on the magnet•Low until the solution clears.
- **s.** Transfer $35 \mu l$ sample to a new tube strip.

2.2 cDNA Amplification





a. Prepare cDNA Amplification Mix on ice. Vortex and centrifuge briefly.

cDNA Amplification Reaction Mix Add reagents in the order listed	PN	1Χ (μl)	4X + 10% (μl)	8X + 10% (μl)
○ Amp Mix	2000047/ 2000103	50	220	440
Feature cDNA Primers 1 Verify name & PN Use indicated primer only	2000096	15	66	132
Total	-	65	286	572

- b. Add 65 μl cDNA Amplification Reaction Mix to 35 μl sample.
- c. Pipette mix 15x (pipette set to 90 µl). Centrifuge briefly.
- d. Incubate in a thermal cycler with the following protocol.

Lid Temperature	Reaction Volume	Run Time
105°C	100 µl	~30-45 min
Step	Temperature	Time
1	98°C	00:03:00
2	98°C	00:00:15
3	63°C	00:00:20
4	72°C	00:01:00
5	Go to Step 2, see table	e below for total # of cycles
6	72°C	00:01:00
7	4°C	Hold

The optimal number of cycles is a trade-off between generating sufficient final mass for library construction and minimizing PCR amplification artifacts. The number of cDNA cycles should also be reduced if large numbers of cells are sampled.

Recommended starting point for cycle number optimization.

Targeted Cell Recovery	Total Cycles
<500	13
500-6,000	12
>6,000	11




e. Store at 4°C for up to 72 h or or −20°C for ≤1 week, or proceed to the next step.

Step Overview (steps 2.2 & 2.3)

Amplification Products Generated in Step 2.2 - cDNA Amplification Post GEM-RT Cleanup Products Feature cDNA Primers 1 PN-2000096 cDNA Amplification Amplification **Products** Poly(dT)VN cDNA from poly-adenylated mRNA Capture Seg 1 Feature Barcode cDNA from Feature Barcoding compatible sgRNA

Step 2.3 - cDNA Cleanup - SPRIselect Overview

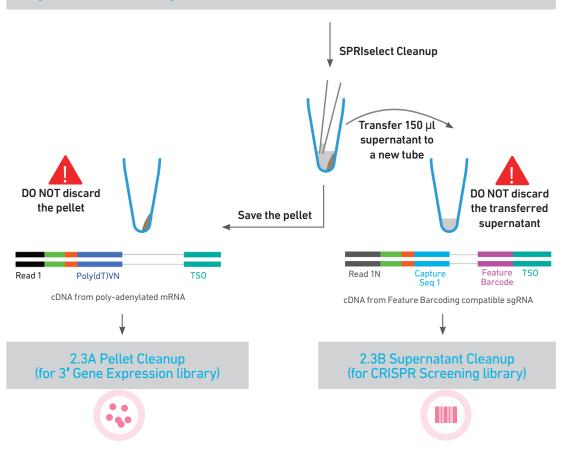

Step 2 cDNA Amplification & QC

2.3 cDNA Cleanup – SPRIselect

- a. Vortex to resuspend the SPRIselect reagent. Add 60 μ l SPRIselect reagent (0.6X) to each sample and pipette mix 15x (pipette set to 150 μ l).
- b. Incubate 5 min at room temperature.
- c. Place on the magnet•High until the solution clears.



- d. Transfer and save 150 μl supernatant in a new tube strip without disturbing the pellet. Maintain at room temperature. DO NOT discard the transferred supernatant (cleanup for CRISPR Screening library construction).
- **e.** DO NOT discard the pellet (cleanup for 3' Gene Expression library construction). **Immediately** proceed to Pellet Cleanup (step 2.3A).



2.3A Pellet Cleanup

(for 3' Gene Expression library)

- i. Add 200 µl 80% ethanol to the pellet. Wait 30 sec.
- ii. Remove the ethanol.
- iii. Repeat steps i and ii for a total of 2 washes.
- iv. Centrifuge briefly and place on the magnet•Low.
- v. Remove any remaining ethanol. Air dry for 2 min. DO NOT exceed 2 min as this will decrease elution efficiency.
- vi. Remove from the magnet. Add 40.5 μl Buffer EB. Pipette mix 15x.
- vii. Incubate 2 min at room temperature.
- viii. Place the tube strip on the magnet•Hlgh until the solution clears.
- ix. Transfer 40 µl sample to a new tube strip.
- X. STOP

Store at 4°C for up to 72 h or at -20°C for up to 4 weeks, or proceed to step 2.4 followed by step 3 for 3' Gene Expression Library Construction.



2.3B Transferred Supernatant Cleanup

(for CRISPR Screening library)


- i. Vortex to resuspend the SPRIselect reagent.
 Add 60 μl SPRIselect reagent (1.2X) to 150 μl of the transferred supernatant and pipette mix 15x (pipette set to 150 μl).
- ii. Incubate for 5 min at room temperature.
- iii. Place on the magnet•High until the solution clears.
- iv. Remove supernatant.
- v. Add 200 µl 80% ethanol to the pellet. Wait 30 sec.
- vi. Remove the ethanol.
- vii. Repeat steps v and vi for a total of 2 washes.
- viii. Centrifuge briefly and place on the magnet•Low.
- ix. Remove any remaining ethanol. Air dry for 2 min. DO NOT exceed 2 min as this will decrease elution efficiency.
- x. Remove from the magnet. Add 50.5 μ l Buffer EB. Pipette mix 15x.
- xi. Incubate 2 min at room temperature.
- xii. Place the tube strip on the magnet•**High** until the solution clears.
- xiii.Transfer **50 μl** sample to a new tube strip.
- Store at 4°C for up to 72 h or at -20°C for up to 4 weeks, or proceed directly to step 4 for CRISPR Screening Library Construction.

Step 2 cDNA Amplification & QC

2.4 Post cDNA Amplification QC & Quantification

a. Run 1 µl of sample from Pellet Cleanup (step 2.3A-x), diluted 1:10 on an Agilent Bioanalyzer High Sensitivity chip. DO NOT run sample from 2.3B Transferred Supernatant Cleanup step.


For input cells with low RNA content (<1pg total RNA/cell), 1 µl undiluted product may be run. Lower molecular weight product (35 – 150 bp) may be present. This is normal and does not affect sequencing or application performance.

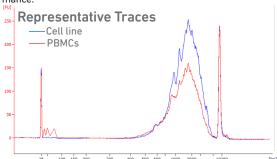


EXAMPLE CALCULATION

i. Select Region

Under the "Electropherogram" view choose the "Region Table". Manually select the region of ~200 - ~9000 bp

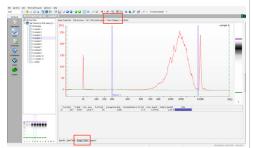

ii. Note Concentration [pg/µl]

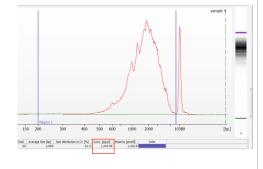

iii. Calculate

Multiply the cDNA concentration [pg/ μ l] reported via the Agilent 2100 Expert Software by the elution volume (40 µl) of the Post cDNA Amplification Reaction Clean Up sample (taking any dilution factors into account) and then divide by 1000 to obtain the total cDNA yield in ng.

Example Calculation of cDNA Total Yield

Concentration: 1244.06_pg/µl Elution Volume: 40

Dilution Factor: 10

Total cDNA Yield

- = Conc'n (pg/µl) x Elution Volume x Dilution Factor 1000 (pg/ng)
- $= 1244.06(pg/\mu l) \times 40 \times 10 = 497.62 ng$ 1000 (pg/ng)



Carry forward ONLY 25% of total cDNA yield into 3' Gene Expression Library Construction (step 3)

- = 0.25 x Total cDNA yield
- = 0.25 x 497.62= 124.4 ng

Refer to step 3.5 for appropriate number of Sample Index PCR cycles based on carry forward cDNA yield/input cDNA.

Alternate Quantification Methods:

Agilent TapeStation. See Appendix for representative traces

Agilent Bioanalyzer or Agilent TapeStation are the recommended methods for accurate quantification.

(If using Qubit Fluorometer and Qubit dsDNA HS Assay Kit, see Appendix)

Step 3

3' Gene Expression Library Construction

- 3.1 Fragmentation, End Repair & A-tailing
- 3.2 Post Fragmentation End Repair & A-tailing Double Sided Size Selection – SPRIselect
- 3.3 Adaptor Ligation
- 3.4 Post Ligation Cleanup – SPRIselect
- 3.5 Sample Index PCR
- 3.6 Post Sample Index PCR Double Sided Size Selection – SPRIselect
- 3.7 Post Library Construction QC

3.0 3' Gene Expression **Library Construction**




GET STARTED!						
Action	Item	10x PN	Preparation & Handling	Storage		
Equilibrate to Room Temperature	Fragmentation Buffer	2000091	Vortex, verify no precipitate, centrifuge briefly.	-20°C		
	Adaptor Oligos	2000094	Vortex, centrifuge briefly.	-20°C		
	Ligation Buffer	2000092	Vortex, verify no precipitate, centrifuge briefly.	-20°C		
	SI Primer	2000095	-	-20°C		
	Single Index Plate T Set A	2000240	-	-20°C		
	Beckman Coulter SPRIselect Reagent	-	Manufacturer's recommendations.	-		
	Agilent TapeStation Screen Tape and Reagents If used for QC		Manufacturer's recommendations.			
	Agilent Bioanalyzer High Sensitivity kit If used for QC	-	Manufacturer's recommendations.	-		
Place on Ice	Fragmentation Enzyme	2000090/ 2000104	Centrifuge briefly.	-20°C		
	DNA Ligase	220110/ 220131	Centrifuge briefly.	-20°C		
	○ Amp Mix	2000047/ 2000103	Centrifuge briefly.	-20°C		
	KAPA Library Quantification Kit for Illumina Platforms	-	Manufacturer's recommendations.	-		
Obtain	Qiagen Buffer EB	-	-	Ambient		
	10x Magnetic Separator	230003	See Tips & Best Practices.	Ambient		
	Prepare 80% Ethanol Prepare 20 ml for 8 reactions	-	Prepare fresh.	Ambient		

Step Overview (Step 3.1d)

Correlation between input & library complexity

A Single Cell 3' Gene Expression library is generated using a fixed proportion (10 μ l, 25%) of the total cDNA (40 μ l) obtained at step 2.3A-ix. The complexity of this library will be comparable to one generated using a higher proportion (>25%) of the cDNA. The remaining proportion (30 µl, 75%) of the cDNA may be stored at 4°C for up to 72 h or at -20°C for longer-term storage (up to 4 weeks).

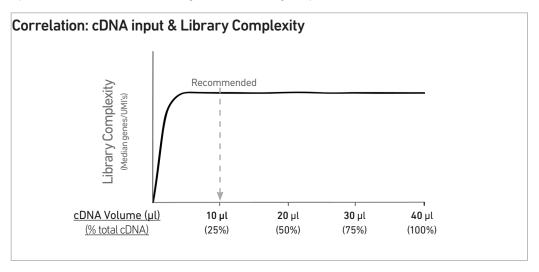

Note that irrespective of the total cDNA yield (ng), which may vary based on cell type, targeted cell recovery etc., this protocol has been optimized for a broad range of input mass (ng), as shown in the example below. The total number of SI PCR cycles (step 3.5e) should be optimized based on carrying forward a fixed proportion (10 µl, 25%) of the total cDNA yield calculated during Post cDNA Amplification QC & Quantification (step 2.4).

Example: Library Construction Input Mass & SI PCR Cycles						
Targeted Total cDNA Cell Recovery (ng)		cDNA Input into Fragmentation		SI PCR Cycle		
	Volume (μl)	Mass (ng)	Number			
Low 6	250 ng	10 μl	62.5 ng	13		
High	1900 ng	10 μl	475 ng	10		
Low	1 ng	10 μl	0.25 ng	16		
High	200 ng	10 μl	50 ng	12		
	Targeted Cell Recovery Low High	Targeted Cell Recovery Low 250 ng High 1900 ng Low 1 ng	Targeted Cell Recovery Total cDNA Fragme Volume (μl) Low 250 ng 10 μl High 1900 ng 10 μl Low 1 ng 10 μl High 1900 ng 10 μl	Targeted Cell Recovery Total cDNA Yield (ng) Volume (μl) Low 250 ng 10 μl 62.5 ng High Low 1 ng 1 ng 10 μl 0.25 ng High High		

3.1 Fragmentation, End Repair & A-tailing

a. Prepare a thermal cycler with the following incubation protocol.

Lid Temperature	Reaction Volume	Run Time
65°C	50 μl	~35 min
Step	Temperature	Time
Pre-cool block Pre-cool block prior to preparing the Fragmentation Mix	4°C	Hold
Fragmentation	32°C	00:05:00
End Repair & A-tailing	65°C	00:30:00
Hold	4°C	Hold



- **b.** Vortex Fragmentation Buffer. Verify there is no precipitate.
- c. Prepare Fragmentation Mix on ice. Pipette mix and centrifuge briefly.

Fragmentation Mix Add reagents in the order listed	PN	1Χ (μl)	4X + 10% (μl)	8X + 10% (μl)
Fragmentation Buffer	2000091	5	22	44
Fragmentation Enzyme	2000090/ 2000104	10	44	88
Total	-	15	66	132

d. Transfer ONLY 10 µl purified cDNA sample from Pellet Cleanup (step 2.3A-x) to a tube strip.

Note that only 10 µl (25%) cDNA sample transfer is sufficient for generating 3' Gene Expression library.

The remaining 30 µl (75%) cDNA sample can be stored at 4°C for up to 72 h or at **-20°C** for up to **4 weeks** for generating additional 3' Gene Expression libraries.

- e. Add 25 µl Buffer EB to each sample.
- f. Add 15 µl Fragmentation Mix to each sample.
- g. Pipette mix 15x (pipette set to 35 µl) on ice. Centrifuge briefly.
- h. Transfer into the pre-cooled thermal cycler (4°C) and press "SKIP" to initiate the protocol.

3.2 Post Fragmentation, End Repair & A-tailing **Double Sided Size** Selection – SPRIselect

- a. Vortex to resuspend SPRIselect reagent. Add 30 µl SPRIselect (0.6X) reagent to each sample. Pipette mix 15x (pipette set to 75 µl).
- b. Incubate 5 min at room temperature.
- c. Place on the magnet•High until the solution clears. DO NOT discard supernatant.





- **d.** Transfer **75** μ**l** supernatant to a new tube strip.
- e. Vortex to resuspend SPRIselect reagent. Add 10 µl SPRIselect reagent (0.8X) to each sample. Pipette mix 15x (pipette set to 80 µl).
- f. Incubate 5 min at room temperature.
- g. Place on the magnet. High until the solution clears.





- h. Remove 80 µl supernatant. DO NOT discard any beads.
- i. Add 125 µl 80% ethanol to the pellet. Wait 30 sec.
- i. Remove the ethanol.
- k. Repeat steps i and j for a total of 2 washes.
- l. Centrifuge briefly. Place on the magnet. Low until the solution clears. Remove remaining ethanol. DO NOT over dry to ensure maximum elution efficiency.
- m. Remove from the magnet. Add 50.5 µl Buffer EB to each sample. Pipette mix 15x.
- n. Incubate 2 min at room temperature.
- o. Place on the magnet•High until the solution clears.
- **p.** Transfer **50 μl** sample to a new tube strip.

3.3 **Adaptor Ligation**

a. Prepare Adaptor Ligation Mix. Pipette mix and centrifuge briefly.

Adaptor Ligation Mix Add reagents in the order listed	PN	1Χ (μl)	4X + 10% (μl)	8X + 10% (μl)
Ligation Buffer	2000092	20	88	176
ONA Ligase	220110/ 220131	10	44	88
Adaptor Oligos	2000094	20	88	176
Total	-	50	220	440

- **b.** Add **50 μl** Adaptor Ligation Mix to **50 μl** sample. Pipette mix 15x (pipette set to 90 μl). Centrifuge briefly.
- c. Incubate in a thermal cycler with the following protocol.

Lid Temperature	Reaction Volume	Run Time
30°C	100 μl	15 min
Step	Temperature	Time
1	20°C	00:15:00
2	4°C	Hold

3.4 Post Ligation Cleanup – SPRIselect

- a. Vortex to resuspend SPRIselect Reagent. Add 80 μ l SPRIselect Reagent (0.8X) to each sample. Pipette mix 15x (pipette set to 150 μ l).
- b. Incubate 5 min at room temperature.
- c. Place on the magnet•High until the solution clears.
- d. Remove the supernatant.
- e. Add 200 μl 80% ethanol to the pellet. Wait 30 sec.
- f. Remove the ethanol.
- g. Repeat steps e and f for a total of 2 washes.
- h. Centrifuge briefly. Place on the magnet•Low.
- i. Remove any remaining ethanol. Air dry for 2 min.
 DO NOT exceed 2 min as this will decrease elution efficiency.
- j. Remove from the magnet. Add 30.5 µl Buffer EB. Pipette mix 15x.
- k. Incubate 2 min at room temperature.
- I. Place on the magnet•Low until the solution clears.
- $\boldsymbol{m}.$ Transfer 30 μl sample to a new tube strip.

3.5 Sample Index PCR



- a. Choose the appropriate sample index sets to ensure that no sample indices overlap in a multiplexed sequencing run. Record the 10x sample index name (PN-2000240 Single Index Plate T Set A well ID) used.
- b. Prepare Sample Index PCR Mix.

Sample Index PCR Mix Add reagents in the order listed	PN	1Χ (μl)	4X + 10% (μl)	8X + 10% (µl)
○ AmpMix	2000047/ 2000103	50	220	440
SI Primer	2000095	10	44	88
Total	-	60	264	528

- c. Add 60 µl Sample Index PCR Mix to 30 µl sample.
- d. Add 10 μ l of an individual Single Index to each well and record the well ID used. Pipette mix 5x (pipette set to 90 μ l). Centrifuge briefly.
- e. Incubate in a thermal cycler with the following protocol.

Lid Temperature	Reaction Volume	Run Time
105°C	100 μl	~25-40 min
Step	Temperature	Time
1	98°C	00:00:45
2	98°C	00:00:20
3	54°C	00:00:30
4	72°C	00:00:20
5	Go to step 2, see be	low for # of cycles
6	72°C	00:01:00
7	4°C	Hold



The total cycles should be optimized based on 25% carry forward cDNA yield/input calculated during cDNA QC & Quantification (step 2.4)

 cDNA Input
 Total Cycles

 0.25-25 ng
 14-16

 25-150 ng
 12-14

 150-500 ng
 10-12

 500-1,000 ng
 8-10

 1,000-1,500 ng
 6-8

5

Recommended cycle numbers

>1500 ng



 f. Store at 4°C for up to 72 h or proceed to the next step.

3.6 Post Sample Index PCR Double Sided Size Selection – SPRIselect


- a. Vortex to resuspend the SPRIselect reagent. Add 60 µl SPRIselect Reagent (0.6X) to each sample. Pipette mix 15x (pipette set to 150 µl).
- b. Incubate 5 min at room temperature.
- c. Place the magnet•High until the solution clears. DO NOT discard supernatant.
- **d.** Transfer **150 μl** supernatant to a new tube strip.
- **e.** Vortex to resuspend the SPRIselect reagent. Add **20 µl** SPRIselect Reagent **(0.8X)** to each sample. Pipette mix 15x (pipette set to 150 µl).
- f. Incubate 5 min at room temperature.
- g. Place the magnet•High until the solution clears.
- h. Remove 165 μl supernatant. DO NOT discard any beads.
- i. With the tube still in the magnet, add $200 \,\mu l$ 80% ethanol to the pellet. Wait $30 \, sec.$
- j. Remove the ethanol.
- k. Repeat steps i and j for a total of 2 washes.
- l. Centrifuge briefly. Place on the magnet•Low. Remove remaining ethanol.
- m. Remove from the magnet. Add 35.5 µl Buffer EB. Pipette mix 15x.
- n. Incubate 2 min at room temperature.
- o. Place on the magnet Low until the solution clears.
- **p.** Transfer $35 \mu l$ to a new tube strip.



q. Store at 4°C for up to 72 h or at -20°C for long-term storage.

3.7 **Post Library Construction** QC

Run 1 µl sample at 1:10 dilution on an Agilent Bioanalyzer High Sensitivity chip.

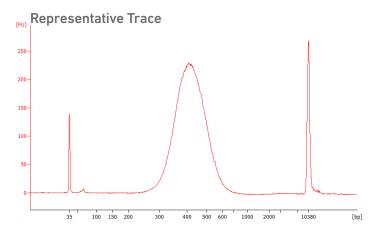

Determine the average fragment size from the Bioanalyzer trace. This will be used as the insert size for library quantification.

Alternate QC Method:

Agilent TapeStation. See Appendix for representative traces

See Appendix for Post Library Construction Quantification

Step 4

CRISPR Screening Library Construction

- 4.1 Guide RNA cDNA Cleanup— SPRIselect
- 4.2 Feature PCR
- 4.3 Post Feature PCR Cleanup – SPRIselect
- 4.4 Sample Index PCR
- 4.5 Post Sample Index PCR Double Sided Size Selection – SPRIselect
- 4.6 Post Library Construction QC

4.0 **CRISPR Screening Library Construction**



GET START	ED!			
Action	Item	10x PN	Preparation & Handling	Storage
Equilibrate to Room Temperature	Feature SI Primers 1	2000098	-	-20°C
	Feature SI Primers 2	2000099	-	-20°C
	Single Index Plate T Set A	2000240	-	-20°C
	Beckman Coulter SPRIselect Reagent	-	Manufacturer's recommendations.	-
	Agilent TapeStation Screen Tape and Reagents If used for QC		Manufacturer's recommendations.	-
	Agilent Bioanalyzer High Sensitivity kit If used for QC	-	Manufacturer's recommendations.	-
Place on Ice	○ Amp Mix	2000047	Centrifuge briefly.	-20°C
	KAPA Library Quantification Kit for Illumina Platforms	-	Manufacturer's recommendations.	-
Obtain	Qiagen Buffer EB	-	-	Ambient
	10x Magnetic Separator	230003	See Tips & Best Practices.	Ambient
	Prepare 80% Ethanol Prepare 20 ml for 8 reactions	-	Prepare fresh.	Ambient

4.1 Guide RNA cDNA Cleanup - SPRIselect

- a. Vortex to resuspend the SPRIselect reagent. Add 50 µl SPRIselect reagent (1.0X) to 50 μl Transfered Supernatant Cleanup (step 2.3B-xiv) and pipette mix 15x (pipette set to 150 µl).
- b. Incubate 5 min at room temperature.
- c. Place on the magnet•High until the solution clears.
- d. Remove supernatant.
- e. Add 200 µl 80% ethanol to the supernatant. Wait 30 sec.
- f. Remove the ethanol.
- g. Repeat steps e and f for a total of 2 washes.
- h. Centrifuge briefly and place on the magnet•Low.



- i. Remove any remaining ethanol. Air dry for 2 min. DO NOT exceed 2 min as this will decrease elution efficiency.
- j. Remove from the magnet. Add 30.5 µl Buffer EB. Pipette mix 15x.
- k. Incubate 2 min at room temperature.
- l. Place the tube strip on the magnet•High until the solution clears.
- m. Transfer 30 µl sample to a new tube strip.



n. Store at 4°C for up to 72 h or at -20°C for up to a week, or proceed to the next step.

4.2 Feature PCR

a. Prepare Feature PCR Mix on ice. Vortex and centrifuge briefly.

Feature PCR Mix Add reagents in the order listed	PN	1Χ (μl)	4X + 10% (μl)	8X + 10% (µl)
○ Amp Mix	2000047	50	220	440
Feature SI Primers 1	2000098	45	198	396
Total	-	95	418	836

- **b.** Transfer **ONLY 5** μ**l** from Guide RNA cDNA Cleanup (step 4.1m) to a new tube strip. Note that only $\mathbf{5}~\mu\mathbf{l}$ of the DNA sample transfer is sufficient for generating CRISPR Screening library. The remaining 25 µl sample can be stored at 4°C for up to 72 h or at -20°C for up to 4 weeks, for generating additional CRISPR Screening libraries.
- c. Add 95 µl Feature PCR Mix to 5 µl sample.
- **d.** Pipette mix 15x (pipette set to 90 μl). Centrifuge briefly.
- e. Incubate in a thermal cycler with the following protocol.

Lid Temperature	Reaction Volume	Run Time
105°C	100 μl	~30-45 min
Step	Temperature	Time
1	98°C	00:00:45
2	98°C	00:00:20
3	60°C	00:00:05
4	72°C	00:00:05
5	Go to Step 2, repeat	14X for a total of 15 cycles
6	72°C	00:01:00
7	4°C	Hold

4.3 Post Feature PCR Cleanup – SPRIselect

- a. Vortex to resuspend SPRIselect Reagent. Add 100 µl SPRIselect Reagent (1.0X) to each sample. Pipette mix 15x (pipette set to 150 µl).
- b. Incubate 5 min at room temperature.
- c. Place on the magnet•High until the solution clears.
- **d.** Remove the supernatant.
- e. Add 300 µl 80% ethanol to the pellet. Wait 30 sec.
- f. Remove the ethanol.
- g. Repeat steps e and f for a total of 2 washes.
- h. Centrifuge briefly. Place on the magnet•Low.
- i. Remove any remaining ethanol. Air dry for 1 min. DO NOT exceed 1 min as this will decrease elution efficiency.
- j. Remove from the magnet. Add 30.5 µl Buffer EB. Pipette mix 15x.
- k. Incubate 2 min at room temperature.
- I. Place on the magnet•Low until the solution clears.
- m. Transfer 30 µl sample to a new tube strip.

4.4 Sample Index PCR



- a. Choose the appropriate sample index sets to ensure that no sample indices overlap in a multiplexed sequencing run. Avoid using the same sample index sets for both 3' Gene Expression and CRISPR Screening libraries constructed from the same sample. Record the 10x sample index name (PN-2000240 Single Index Plate well ID) used.
- b. Prepare Sample Index PCR Mix.

Sample Index PCR Mix Add reagents in the order listed	PN	1Χ (μl)	4Χ + 10% (μl)	8X + 10% (μl)
○ AmpMix	2000047	50	220	440
Feature SI Primers 2	2000099	35	154	308
Total	-	85	374	748

- c. Transfer ONLY 5 µl Post Feature PCR Cleanup sample (step 4.3m) to a new tube strip. Note that only **5** µl sample transfer is sufficient for generating CRISPR Screening library. The remaining 25 µl sample can be stored at 4°C for up to 72 h or at -20°C for up to 4 weeks, for generating additional CRISPR Screening libraries.
- d. Add 85 μl Sample Index PCR Mix to 5 μl sample (Post Feature PCR Cleanup).
- e. Add 10 μl of an individual Single Index to each well and record their assignment. Pipette mix 5x (pipette set to 90 µl). Centrifuge briefly.
- f. Incubate in a thermal cycler with the following protocol.

Reaction Volume	Run Time
100 μl	~25-40 min
Temperature	Time
98°C	00:00:45
98°C	00:00:20
54°C	00:00:30
72°C	00:00:20
Go to step 2, repeat 4X	for a total of 5 cycles
72°C	00:01:00
4°C	Hold
	100 μl Temperature 98°C 98°C 54°C 72°C Go to step 2, repeat 4X

4.5 Post Sample Index **PCR Double Sided Size** Selection - SPRIselect


- a. Vortex to resuspend the SPRIselect reagent. Add 70 µl SPRIselect Reagent (0.7X) to each sample. Pipette mix 15x (pipette set to 150 µl).
- b. Incubate 5 min at room temperature.
- c. Place the magnet•High until the solution clears. DO NOT discard supernatant.
- **d.** Transfer 160 μl supernatant to a new tube strip.
- e. Vortex to resuspend the SPRIselect reagent. Add 30 µl SPRIselect Reagent (1.0X) to each sample. Pipette mix 15x (pipette set to 150 µl).
- f. Incubate 5 min at room temperature.
- g. Place the magnet. High until the solution clears.
- h. Remove the supernatant.
- i. Add 300 µl 80% ethanol to the pellet. Wait 30 sec.
- i. Remove the ethanol.
- k. Add 200 µl 80% ethanol to the pellet. Wait 30 sec.
- I. Remove the ethanol.
- m. Centrifuge briefly. Place on the magnet•Low.
- n. Remove remaining ethanol. Air dry for 1 min.
- o. Remove from the magnet. Add 30.5 µl Buffer EB. Pipette mix 15x.
- p. Incubate 2 min at room temperature.
- **q.** Place on the magnet•Low until the solution clears.
- r. Transfer 30 μ l to a new tube strip.



s. Store at 4°C for up to 72 h or at -20°C for long-term storage.

4.6 **Post Library Construction** QC

Run 1 µl sample at 1:25 dilution on an Agilent Bioanalyzer High Sensitivity chip.

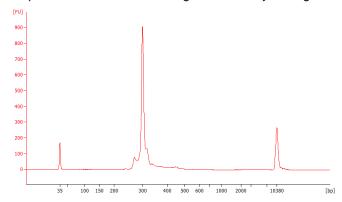

Determine the average fragment size from the Bioanalyzer trace. This will be used as the insert size for library quantification.

Alternate QC Method:

Agilent TapeStation. See Appendix for representative trace

See Appendix for Post Library Construction Quantification

Sequencing




Step 5 Sequencing


Sequencing Libraries

Chromium Single Cell 3' Gene Expression and CRISPR Screening libraries comprise standard Illumina paired-end constructs which begin with P5 and end with P7. These libraries include 16 bp 10x Barcodes at the start of TruSeq Read 1 and Nextera Read 1 (Read 1N) respectively while 8 bp sample index sequences are incorporated as the sample index read. TruSeq Read 1 and TruSeq Read 2 are standard Illumina sequencing primer sites used in paired-end sequencing of Single Cell 3' Gene Expression libraries. Nextera Read 1 (Read 1N) and TruSeg Read 2 are used for paired-end sequencing of Single Cell 3' CRISPR Screening libraries. Sequencing these libraries produce a standard Illumina BCL data output folder.

Chromium Single Cell 3' Gene Expression Library

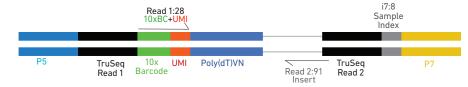

Chromium Single Cell 3' CRISPR Screening Library

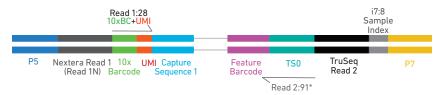

*Minimum required Read 2 length for CRISPR Screening libraries is 70 bp

Illumina Sequencer Compatibility

The compatibility of the listed sequencers has been verified by 10x Genomics. Some variation in assay performance is expected based on sequencer choice. For more information about performance variation, visit the 10x Genomics Support website.

- MiSeq
- NextSeg 500/550*
- HiSeg 2500 (Rapid Run)
- HiSeq 3000/4000
- NovaSeq

*Sequencing Chromium Single Cell libraries on the NextSeq 500/550 platform may yield reduced sequence quality and sensitivity relative to the MiSeq, HiSeq, and NovaSeq platforms. Refer to the 10x Genomics Support website for more information.

Sample Indices

Each sample index in the Single Index Kit T Set A (PN-1000213) is a mix of 4 different sequences to balance across all 4 nucleotides. If multiple samples are pooled in a sequence lane, the sample index name (i.e. the Single Index plate well ID, SI-GA-) is needed in the sample sheet used for generating FASTQs with "cellranger mkfastq". Samples utilizing the same sample index should not be pooled together, or run on the same flow cell lane, as this would not enable correct sample demultiplexing.

Step 5 Sequencing

21.0	Sequencing Depth	Minimum 20,000 read pairs per cell
3' Gene Expression Library Sequencing Depth	Sequencing Type	Paired-end, single indexing
& Run Parameters	Sequencing Read	Recommended Number of Cycles
	Read 1 i7 Index i5 Index Read 2	28 cycles 8 cycles 0 cycles 91 cycles
CDICDD Companies	Sequencing Depth	Minimum 5,000 read pairs per cell
CRISPR Screening Library Sequencing Depth	Sequencing Type	Paired-end, single indexing
& Run Parameters	Sequencing Read	Recommended Number of Cycles
	Read 1 i7 Index i5 Index Read 2	28 cycles 8 cycles 0 cycles 91 cycles

Library Loading

Once quantified and normalized, the 3' Gene Expression and CRISPR Screening libraries should be denatured and diluted as recommended for Illumina sequencing platforms. Refer to Illumina documentation for denaturing and diluting libraries. Refer to the 10x Genomics Support website, for more information.

Instrument	3' Gene Expression libr or 3' Gene Expression Screening librar	+ CRISPR	CRISPR Screening libr	aries only
	Loading Concentration (pM)	PhiX (%)	Loading Concentration (pM)	PhiX (%)
MiSeq	11	1	11	1
NextSeq 500/550	1.8	1	1.8	20**
HiSeq 2500 (RR)	11	1	11	1
HiSeq 4000	240	1	240	1
NovaSeq	150*/300	1	150*/300	1

^{*} Use 150pM loading concentration for Illumina XP workflow.

Library Pooling

The 3' Gene Expression and CRISPR Screening libraries maybe pooled for sequencing, taking into account the differences in cell number and per-cell read depth requirements between each library. Samples utilizing the same sample index should not be pooled together, or run on the same flow cell lane, as this would not enable correct sample demultiplexing.

Library Pooling Example:

Libraries	Sequencing Depth (read pairs per cell)	Library Pooling Ratio
3' Gene Expression library	20,000	4
CRISPR Screening library	5,000	1

^{**} When sequencing CRISPR screening libraries alone to create more diverse clusters for 30 cycles of Read 2.

Troubleshooting



GEMs

STEP

NORMAL

REAGENT CLOGS & WETTING FAILURES

1.4 d After Chip G is removed from the Controller and the wells are exposed



All 8 recovery wells are similar in volume and opacity.



Recovery well G indicates a reagent clog. Recovery well C and E indicate a wetting failure. Recovery wells B, D, and F are normal. Wells A and H contain 50% Glycerol Solution.

1.4 f Transfer GEMs from Chip G Row Labeled 3



All liquid levels are similar in volume and opacity without air trapped in the pipette tips.




Pipette tips C and E indicate a wetting failure. Pipette tip C contains partially emulsified GEMs. Emulsion is absent in pipette tip E. Pipette tip G indicates a reagent clog.

STEP


NORMAL

REAGENT CLOGS & WETTING FAILURES

2.1 a After transfer of the GEMs + **Recovery Agent**

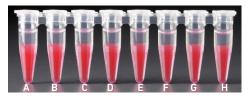

All liquid levels are similar in the aqueous sample volume (clear) and Recovery Agent/Partitioning Oil (pink).

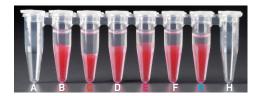

Tube G indicates a reagent clog has occurred. There is a decreased volume of aqueous layer (clear).

Tube C and E indicate a wetting failure has occurred. There is an abnormal volume of Recovery Agent/Partitioning Oil (pink).

2.1 b After aspiration of Recovery Agent/ Partitioning Oil




All liquid volumes are similar in the aqueous sample volume (clear) and residual Recovery Agent/Partitioning Oil (pink).



Tube G indicates a reagent clog has occurred. There is a decreased volume of aqueous layer (clear). There is also a greater residual volume of Recovery Agent/Partitioning Oil (pink).

Tube C and E indicate a wetting failure has occurred. There is an abnormal residual volume of Recovery Agent/Partitioning Oil (pink).

2.1 d After addition of **Dynabeads** Cleanup Mix

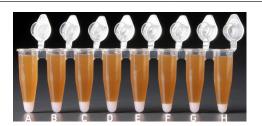

All liquid volumes are similar after addition of the Dynabeads Cleanup Mix.



Tube G indicates a reagent clog has occurred. There is an abnormal ratio of Dynabeads Cleanup Mix (brown) to Recovery Agent/Partitioning Oil (appears white).

Tube C and E indicate a wetting failure has occurred. There is an abnormal ratio of Dynabeads Cleanup Mix (brown) to Recovery Agent/Partitioning Oil (appears white).

If a channel clogs or wetting failure occurs during GEM generation, it is recommended that the sample be remade. If any of the listed issues occur, take a picture and send it to support@10xgenomics.com for further assistance.

Chromium Controller **Errors**

- If the Chromium Controller or the Chromium Single Cell Controller fails to start, an error tone will sound and one of the following error messages will be displayed:
- a. Chip not read Try again: Eject the tray, remove and/or reposition the Chromium Next GEM Secondary Holder assembly and try again. If the error message is still received after trying this more than twice, contact support@10xgenomics.com for further assistance.
- b. Check gasket: Eject the tray by pressing the eject button to check that the 10x Gasket is correctly installed on the Chromium Next GEM Chip. If the error message persists, contact support@10xgenomics.com for further assistance.
- c. Error Detected: Row _ Pressure:
 - i. If this message is received within a few seconds of starting a run, eject the tray by pressing the eject button and check for dirt or deposits on the 10x Gasket. If dirt is observed, replace with a new 10x Gasket and try again. If the error message is still received after trying this more than twice, contact support@10xgenomics.com for further assistance.
 - ii. If this message is received after a few minutes into the run, the Chromium Next GEM Chip must be discarded. Do not try running this Chromium Next GEM Chip again as this may damage the Chromium Controller.
- d. Invalid Chip CRC Value: This indicates that a Chromium Next GEM Chip has been used with an older firmware version. The chip must be discarded. Contact support@10xgenomics.com for further assistance.
- e. Chip Holder Not Present: Open the controller drawer and check if chip holder is present. Insert chip properly into chip holder and retry.
- f. Unauthorized Chip: This indicates that an incompatible non-Next GEM chip has been used with an instrument that only can run Next GEM assays. Use only Chromium Controller (PN-120223;120246) or Chromium Single Cell Controller (PN-120263;120212) to run that chip or chip must be discarded. Contact support@10xgenomics.com for further assistance.
- g. Endpoint Reached Early: If this message is received, contact support@10xgenomics.com for further assistance.

Appendix

Post Library Construction Quantification Agilent TapeStation Traces Compatible sgRNA Specifications Oligonucleotide Sequences

Post Library Construction Quantification

- a. Thaw KAPA Library Quantification Kit for Illumina Platforms.
- b. Dilute 2 µl sample with deionized water to appropriate dilutions that fall within the linear detection range of the KAPA Library Quantification Kit for Illumina Platforms. (For more accurate quantification, make the dilution(s) in duplicate).
- c. Make enough Quantification Master Mix for the DNA dilutions per sample and the DNA Standards (plus 10% excess) using the guidance for 1 reaction volume below.

Quantification Master Mix	1X (μl)
SYBR Fast Master Mix + Primer	12
Water	4
Total	16

- d. Dispense 16 µl Quantification Master Mix for sample dilutions and DNA Standards into a 96 well PCR plate.
- e. Add 4 μl sample dilutions and 4 μl DNA Standards to appropriate wells. Centrifuge briefly.
- f. Incubate in a thermal cycler with the following protocol.

Step	Temperature	Run Time
1	95°C	00:03:00
2	95°C	00:00:05
3	67°C	00:00:30
4	Go to Step 2, 29X (Total 30 cycles)	


g. Follow the manufacturer's recommendations for qPCR-based quantification. For library quantification for sequencer clustering, determine the concentration based on insert size derived from the Bioanalyzer/TapeStation trace.

Agilent TapeStation Traces


Agilent TapeStation Traces

Agilent Tape Station High Sensitivity D5000 ScreenTape was used. Protocol steps correspond to the Chromium Next GEM Single Cell 3 v3.1 Reagent Kits User Guide with Feature Barcoding technology for CRISPR Screening (CG000205)


Protocol Step 2.4 - cDNA QC & Quantification

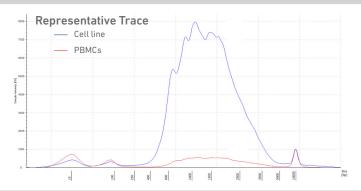

Protocol Step 3.7 - Post Library Construction QC

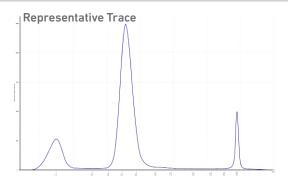

Protocol Step 4.6 - Post Library Construction QC

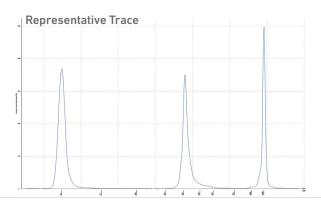

Alternate QC Method:

Qubit Fluorometer and Qubit dsDNA HS Assay Kit

Multiply the cDNA concentration reported via the Qubit Fluorometer by the elution volume (40 µl) to obtain the total cDNA yield in ng. To determine the equivalent range using the Agilent 2100 Expert Software, select the region encompassing 35-10,000 bp.

Compatible sgRNA **Specifications**


Integration of Capture Sequence 1 and Capture Sequence 2 in sgRNA

A representative sgRNA sequence along with the specific capture sequences integrated in two different locations in the sgRNA are shown

Capture Sequence 1


Capture Sequence 1 on Gel Bead: 5'-TTGCTAGGACCGGCCTTAAAGC-3'

Capture Sequence 1 integrated in sgRNA hairpin

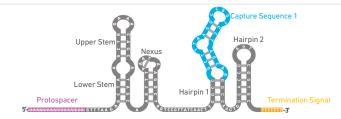

5-NNNNNNNNNNNNNNNNNNNNNNNNTTTAAGAGCTAAGCTGGAAACAGCATAGCAAGTTTAAATAAGGCTAGTCCGTTATCAACTTggccGCTTTTAAGGCCGGTCCTAGCAAggccAAGTGGCACCGAGTCCGTGTCTTTTTT-3'

Capture Sequence 1 integrated in sgRNA 3'-end

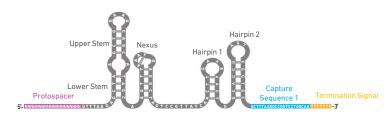


5-NNNNNNNNNNNNNNNNNNNNGTTTAAGAGCTAAGCTGGAAACAGCATAGCAAGTTTTAAATAAGGCTAGTCCGTTATCAACTTgaaaAAGTGGCACCGAGTCCGGTGCGCTTTAAGGCCCGGTCCTAGCAAGTTTTTTT-3*

Capture Sequence 2


Capture Sequence 2 on Gel Bead oligonucleotide strand: 5'-CCTTAGCCGCTAATAGGTGAGC-3'

Capture Sequence 2 integrated in sgRNA hairpin

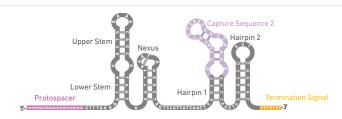

5-NNNNNNNNNNNNNNNNNNNNTTTAAGAGCTAAGCTGGAAACAGCATAGCAAGTTTAAATAAGGCTAGTCCGTTATCAACTTggccGCTCACCTATTAGCGGCCTAAGGggccAAGTGGCACCGAGTCGGTGCTTTTTTT-3'

Capture Sequence 2 integrated in sgRNA 3'-end

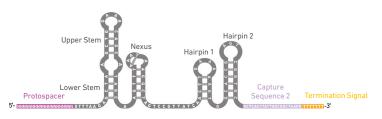


5- NINNINNINNINNINNINNINNINNTTTAAGAGCTAAGCTGGAAACAGCATAGCAAGTTTAAATAAGGCTAGTCCGTTATCAACTTgaaaAAGTGGCACCGAGTCGGTGCGCTCACCTATTAGCGGCTAAGGTTTTTTT -3'

Oligonucleotide Sequences


Protocol steps correspond to the Chromium Next GEM Single Cell 3 v3.1 Reagent Kits User Guide with Feature Barcoding technology for CRISPR Screening (CG000205)

Protocol Step 1.5 - GEM-RT Incubation



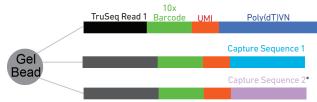

*Capture Sequence 2 is not illustrated in the protocol steps below

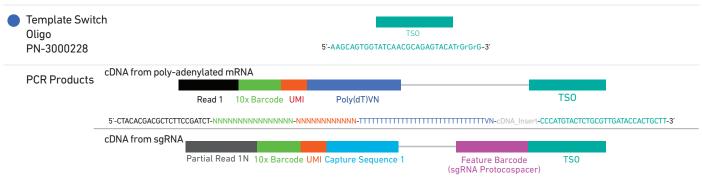

Protocol Step 2.2 - cDNA Amplification

Primers 1 PN-2000096

Amplifies cDNA from poly-adenylated mRNA

Forward Primer:

Partial Read 1
5'-CTACACGACGTCTTCCGATCT-3'

Reverse Primer:

Partial TSO
5'-AAGCAGTGGTATCAACGCAGAG-3'

Amplifies cDNA from sgRNA

Forward Primer: Partial Read 1N 5'-GCAGCGTCAGATGTGTATAAGAGACAG-3'

Reverse primer: Partial TS0


5'-AAGCAGTGGTATCAACGCAGAG-3'

Amplification Ar Products

Amplified cDNA from poly-adenylated mRNA

Read 1 10x Barcode UMI Poly(dT)VN TS0



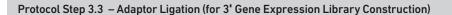

Adaptor Oligos PN -2000094




5'- GATCGGAAGAGCACACGTCTGAACTCCAGTCA-3' 3'-TCTAGCCTTCTCG-5'


Ligation Product 10x Barcode UMI $\mathsf{Poly}(\mathsf{dT})\mathsf{VN}$ Read 1

Protocol Step 3.5 - Sample Index PCR (for 3' Gene Expression Library Construction) SI Primer PN-2000095 Partial Read 1 Sample Partial Read 2 P5 Single Index T 5'-AATGATACGGCGACCACCGAGATCT-ACACTCTTTCCCTACACGACGCTC-3' 5'-CAAGCAGAAGACGGCATACGAGAT-NNNNNNNN GTGACTGGAGTTCAGACGTGT-3' Set A PN-2000240 Sample Index **PCR Product** Read 1 10x Barcode UMI Poly(dT)VN Sample Read 2 Index

5-AATGATACGGCGACCACCGAGATCT-ACACTCTTTCCCTACACGACGCTCTTCCCGATCT-NNNNNNN-ATCTCGTATGCCGTCTTCTGCTTG-3'

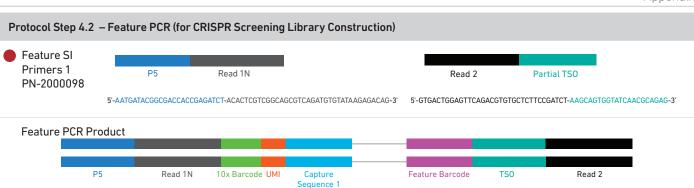

5-AATGATTACGGCGACCACCGGGATCT-ACACTCGTCGGCAGCGTCAGATGTGTATAAGAGACAG-16N-12N-TTGCTAGGACCGCCCTTAAAGC-sgRNA_Feature_Barcode-CCCATGTACTCTGCGTTGATACCACTGCTT-AGATCGGAAGAGCACACGTCTGAACTCCAGTCAC-3'

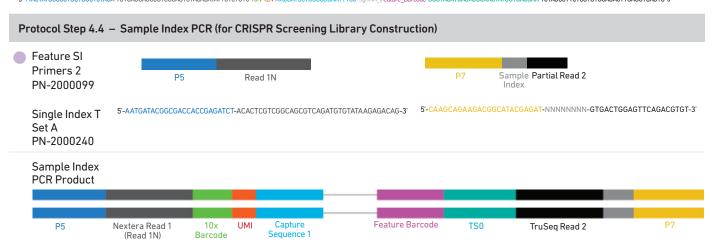

5-AATGATACGGCGACCACCGAGATCT-ACACTCGTCCGCAGCGCTCAGATGTGTATAAGAGACAG-Inn-12N-TTGCTAGGACCGGCCTTAAAGC-sglivin_reature_Barcode-CCCATGTACTCTGCGTTGATACCACTGCTT-AGATCGGAAGAGCACACGTCTGAACTCCAGTCAC-8N-ATCTCGTATGCCGTCTTCTGCTTG-3 3-TTACTATGCCGCTGGTGGCTCTAGA-TGTGAGCAGCCGTCCACAGTTCTACACATATTCTCTGTC-16N-12N-AACGATCTGGCCGGAATTTCG-3gNA_Feature_Barcode-G6GTACATGAGACCCAACTATGGTGACGAA-TCTAGCCTTCTCGTGTGCAGACTTGAGGTCAGTG-8N-TAGAGCATAGGGCAGAACACGAAC-5