TECHNICAL NOTE

Targeted Gene Expression Custom Panel Performance Metrics and Best Practices

Introduction

The Targeted Gene Expression product is a modular enrichment kit designed to enrich libraries for relevant genes, while decreasing sequencing requirements by up to 90%. With this product, panels of baits are used to target a selection of genes. When used to enrich whole transcriptome libraries, these panels enable significantly reduced sequencing costs while accurately reflecting gene expression information from the corresponding whole transcriptome analysis (WTA) parent library for targeted genes. 10x Genomics provides pre-designed panels of baits. For customers interested in creating custom panels, the 10x Genomics Custom Panel Designer enables two customization options: adding genes to pre-designed panels (addon panels) and designing fully custom panels. Panels can be a combination of human genes and exogenous sequences. This Technical Note describes the technical performance for each customization option and outlines best practices for panel design and sequencing.

Method

Each customization option was tested by choosing sets of genes arbitrarily to use as either add-on panels or as fully custom panels. Each custom bait was designed using the Custom Panel Designer available on the 10x Genomics website. Pre-designed panels used in this Technical Note were the Human Pan-Cancer Panel (PN-1000247), the Human Immunology Panel (PN-1000246), and the Human Gene Signature Panel (PN-1000245).

Targeted libraries were generated using the reagents and protocol described in the Targeted Gene Expression Reagent Kits User Guide (Document CG000293) using single cell gene expression libraries previously sequenced for WTA as shown in Figure 1. Add-on panels were tested with Chromium Single Cell 3' Gene Expression (v3.1) and 5' Gene Expression (v1.1, v2). Fully custom panels were tested with Chromium Single Cell 3' Gene Expression (v3, v3.1) and 5' Gene Expression (v1, v1.1, v2) libraries. Panels were tested across 53 human samples. A list of sample types is provided in the Appendix.

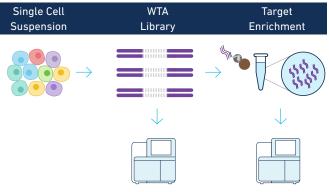


Figure 1. Targeted Gene Expression workflow overview. Whole Transcriptome Analysis (WTA) libraries are generated from single cell suspensions. WTA libraries are incubated with biotinylated baits from Targeted Gene Expression panels to generate targeted libraries. WTA and targeted libraries are both sequenced. These data are used for metrics comparisons.

Add-on Panel Experiment

Pre-designed panels customized with 10 or 200 additional genes (add-on panel) were generated to test the number of genes in the add-on panel as well as relative expression level as shown in Table 1. Bait sequences for each add-on panel were obtained from the Custom Panel Designer and ordered from Integrated DNA Technologies (IDT) as NGS Discovery Pools. A list of genes for each add-on panel is provided in the Appendix.

Add-on Panel	Number of Genes	Expected Relative Expression Level	Number of Samples
Cancer-High	10	High	10
Cancer-Low	10	Low	9
Immune-High	10	High	4
Immune-Low	10	Low	9
Immune-200	200	Even Distribution	24

Table 1. Summary of add-on panels tested with pre-designed panels. Genes in Cancer and Immune add-on panels appear in the Human Pan-Cancer Panel and Human Immunology Panel respectively. Immune-200 genes span a range of expected expression levels. Samples were run in pulldown reactions with 1, 4, or 8 input libraries.



Fully Custom Panel Experiment

Fully custom panels were created by selecting a 10, 25, 50, 100, or 200 gene subset from the Human Immunology Panel. Genes were chosen at random; however, genes with very high or very low levels of expression were omitted. Fully custom panels were ordered from IDT as NGS Discovery Pools. A list of genes for fully custom panels is provided in the Appendix.

Fully Custom Panel	Expected Panel Expression Level	Number of Samples
Immune-10	0.10%	12
Immune-25	0.10%	16
Immune-25	0.25 - 1%	4
Immune-50	0.25 - 1%	11
Immune-50	2 - 4%	4
Immune-100	0.25 - 1%	18
Immune-200	0.25 - 1%	17
Immune-200	2 - 4%	9

Table 2. Summary of conditions tested for fully custom panels. Samples were run in pulldown reactions with 1, 2, 4, or 8 input libraries.

Different combinations of samples and fully custom panels were chosen to create reactions that represented panels with a range of expression levels (Table 2). The targeted-depth subcommand in Cell Ranger was used to calculate the "Fraction of Reads from Targeted Genes" in WTA libraries, and this value is used as a proxy for the expression level of the targeted genes in that sample (Expected Panel Expression Level). For more information, consult the Targeted Depth section of the 10x Genomics Support website.

Sequencing Depth

To calculate parent-targeted comparison metrics, both libraries were computationally downsampled and compared at the following depths:

Parent WTA Libraries: 20,000 or 50,000 read pairs per cell.

Targeted Gene Expression Libraries: two-fold the number of on target panel reads in the parent WTA library.

Example: Parent WTA library with 6% on target reads=6,000 read pairs per cell ($0.06 \times 50,000 \times 2$) for the associated target library. Sequencing libraries were computationally downsampled and compared at the read numbers described in the calculation above.

<u>Data Analysis</u>

Data from Targeted Gene Expression libraries were analyzed by comparing parent and targeted data on a matched (parent derived) barcode set, and metrics related to UMI recovery that compare parent WTA and targeted libraries were calculated over the cell-associated barcodes. In cases where add-on gene panels were used, metrics related to UMI recovery (targeted total UMI recovery, fraction targeted genes with \geq 80% UMI recovery, and pseudo-bulk targeted UMI R²) can be calculated over the add-on genes or pre-designed panel genes separately.

Metrics Analyzed

Targeting Metrics:

- Reads mapped confidently to the targeted transcriptome: Fraction of reads that mapped to a unique and targeted gene in the transcriptome. The read must be consistent with annotated splice junctions. These reads are considered for UMI counting. This metric can be calculated for the parent WTA library, the targeted library, or both, and does not require a matched WTA library control.
- Targeted total UMI recovery: Fraction of cell-associated, on target UMIs recovered (pseudo-bulk) at the sequencing depth described above. Requires a matched WTA library control. This metric is calculated on matched cell barcodes across WTA and targeted libraries.
- Fraction targeted genes with ≥80% UMI recovery: Fraction of observed (defined as ≥ 10 UMI in WTA library) panel genes in a WTA library control for which ≥80% of UMIs are recovered in the matched targeted library at the sequencing depth described above. Requires a matched WTA library control.
- Pseudo-bulk targeted UMI R²: Pearson correlation coefficient of cell-associated panel UMIs across WTA and targeted libraries. Requires a matched WTA library control.

Cell calling concordance: A comparison of the number of cells detected, calculated as the number of cells detected in the targeted library divided by the number detected in the WTA library.

Results

Targeting Metrics - Add-on Panels:

Targeting metrics were assessed using similar criteria and thresholds for optimal performance used to demonstrate the performance of pre-designed panels as shown in Tables 3 and 4. Consult the Targeted Gene Expression Pre-designed Panel Performance Metric Technical Note (Document CG000345) for more information. All Single Cell 3' Gene Expression libraries passed the reads mapped confidently to the targeted transcriptome metric, and most (23/26) Single Cell 5' Gene Expression libraries passed the same metric as shown in Figure 2. Single Cell 5' Gene Expression libraries typically have a decreased fraction of reads mapped to the targeted transcriptome due to slightly higher levels of antisense reads present in these libraries compared to Single Cell 3' Gene Expression libraries.

Targeted total UMI recovery is a measure of how many mRNA molecules (i.e. UMIs for genes of interest) in the parent WTA library are recovered in the targeted library. For all libraries tested, this ratio is above 0.8 for entire panel (Figure 3A) and when analyzing pre-designed panel genes alone (Figure 3B) or add-on panel genes alone (Figure 3C).

UMI correlations between parent and targeted libraries typically exceeded the threshold of $R^2 > 0.9$ for the entire panel, pre-

Pre-designed Panels Customized with Add-on Panels

designed panel genes alone, or add-on panel genes alone (Figure 4A-C) for all samples tested. In a few instances, this correlation for the add-on panel genes is lower than 0.9. This is often due to calculating the correlation coefficient over very few genes.

Examining UMI recovery on a gene-by-gene basis is another way to assess the evenness of enrichment for targeted genes. The distribution of UMI recovery values of a representative example is shown for each of the pre-designed/add-on panel combinations (Figure 5). The majority of targeted genes, whether from the pre-designed or add-on panel genes, have greater than 80% UMI recovery (dashed line). The narrow distribution of UMI recovery values across genes in the pre-designed and add-on panels demonstrates the even enrichment of the targeted genes. Genes with low expression show additional variability, primarily because 20,000 or 50,000 read pairs per cell is not sufficient to read all UMIs for these genes in the parent WTA library. The UMI recovery metric is less variable for genes with higher expression because the number of UMIs detected in the parent WTA library with a set number of read pairs is a better reflection of the true number of UMIs for those genes in the parent WTA library, resulting in UMI recovery values near 1. For genes with low expression, the high degree of enrichment provided by the targeted approach results in UMI recovery >1.

Targeting Metric	Performance Threshold	Cancer-High % Pass	Cancer-Low % Pass	Immune-High % Pass	Immune-Low % Pass	Immune-200 % Pass
Reads Mapped Confidently to the Targeted Transcriptome	80% 70%	100%	100%	100%	89%	92%
Fraction Targeted Genes with ≥80% UMI Recovery	0.9	100%	100%	100%	100%	88%
Targeted Total UMI Recovery	0.8	100%	100%	100%	100%	100%
Pseudo-bulk Targeted UMI R ²	0.9	100%	100%	100%	100%	100%

Table 3. Percentage of pre-designed panels with add-on panels that meet optimal performance thresholds. The 80% performance threshold (red) applies to Single Cell 3' Gene Expression libraries, while the 70% performance threshold (blue) applies to Single Cell 5' Gene Expression libraries. Some Single Cell 5' Gene Expression samples are marginally below threshold due to elevated antisense reads. Pre-designed panels customized with add-on panels may require more sequencing depth for full UMI recovery.

Add-on Panels Alone

Targeting Metric	Performance Threshold	Cancer-High % Pass	Cancer-Low % Pass	Immune-High % Pass	Immune-Low % Pass	Immune-200 % Pass
Fraction Targeted Genes with ≥80% UMI Recovery	0.9	100%	67%	100%	100%	100%
Targeted Total UMI Recovery	0.8	100%	100%	100%	100%	100%
Pseudo-bulk Targeted UMI R ²	0.9	90%	78%	100%	100%	100%

Table 4. Percentage of samples that meet optimal performance thresholds based on add-on panels alone. For add-on panels with few genes, suboptimal performance of a single gene can prevent the sample from meeting performance thresholds. R² noise in small panels is due to the lower number of genes.

Targeting Metrics for Add-on Panels

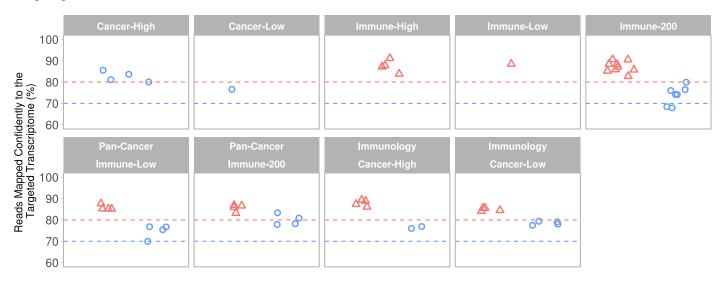


Figure 2. Reads mapped confidently to the targeted transcriptome for add-on panels. The names of the pre-designed panel and add-on panel used are indicated above each plot. Dashed line indicates threshold for optimal performance. Red triangles indicate Single Cell 3' Gene Expression libraries, blue circles indicate Single Cell 5' Gene Expression libraries. Thresholds differ for Single Cell 3' and 5' libraries for reasons described in the Results.

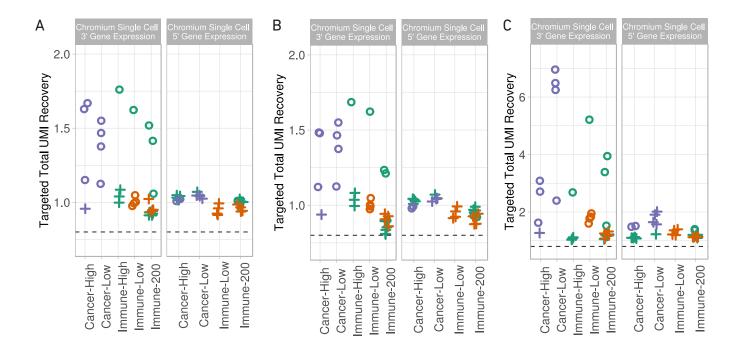


Figure 3. Targeted total UMI recovery for add-on panels, pre-designed panels or both. Each data point represents a single Targeted Gene Expression library for pre-designed panels customized with an add-on panel (A). The same data were calculated for the pre-designed panel alone (B) or the add-on panel alone (C). Data points are colored according to the pre-designed panel used: Human Gene Signature Panel (green), Human Pan-Cancer Panel (orange), Human Immunology Panel (purple). The targeted parent comparison was made using either 20,000 (circle) or 50,000 (plus sign) read pairs per cell from the parent WTA library. Dashed line indicates threshold for optimal performance.

Targeting Metrics for Add-on Panels

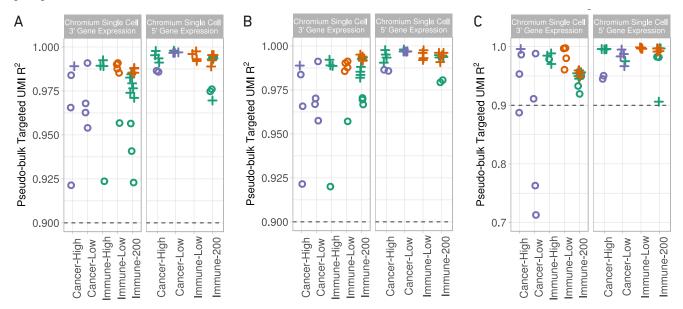


Figure 4. Pseudo-bulk targeted UMI R². Each data point represents a single Targeted Gene Expression library for pre-designed panels customized with an add-on panel (A). The same data were calculated for the pre-designed panel alone (B) or the add-on panel alone (C). Data points are colored according to the pre-designed panel used: Human Gene Signature Panel (green), Human Pan-Cancer Panel (orange), Human Immunology Panel (purple). The targeted parent comparison was made using either 20,000 (circle) or 50,000 (plus sign) read pairs per cell from the parent WTA library. In (C), the data point for one sample (A375 cell line sample with Gene Signature and Immune-Low panels) was omitted because only one gene of the add-on panel was observed in the parent WTA library; thus, pseudo-bulk targeted UMI R² cannot be calculated. Dashed line indicates threshold for optimal performance.

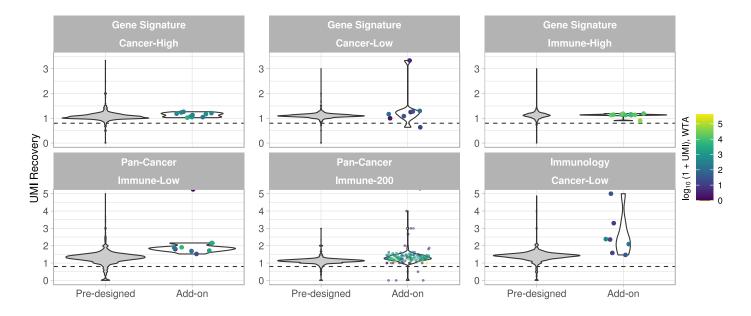


Figure 5. UMI recovery for 10 and 200 add-on panels. Representative samples for add-on panels tested. The names of the pre-designed panel and addon panel used are indicated above each plot. UMI recovery distributions for pre-designed panel genes are shown separately from the add-on genes. Each data point represents the UMI recovery value for an individual gene and data points are color coded by expression level (i.e. the number of UMIs in the parent WTA library). Data points are omitted when no UMIs are detected for a targeted gene in both the WTA and the targeted library. Data points appear clipped at the y-axis when no UMIs are detected for a targeted gene in the WTA library, but have a non-zero UMI count in the targeted library. Dashed line indicates threshold for optimal performance.

Targeting Metrics - Fully Custom Panels

Targeting metrics were assessed using similar criteria and thresholds for optimal performance used to demonstrate the performance of pre-designed panels. Consult the Targeted Gene Expression Pre-designed Panel Performance Metric Technical Note (Document CG000345) for more information. Fully custom panels perform well based on these metrics (Table 5). The expression level of fully custom panels is determined by both the sample and the genes on the panel. Figure 6 demonstrates that the same panel of genes will have different expression levels in different samples; thus, the fraction of reads in a WTA library for that panel of genes will vary and is only weakly correlated with the number of genes on the panel.

Reads confidently mapped to the targeted transcriptome values are generally on par with the pre-designed panels, but can be variable when the expression level of panel genes is low (Figure 7A). This can result from the enrichment of related, but nontargeted, genes and is related to the selection of genes in the fully custom panel (refer to Discussion). Despite a higher percentage of reads mapped to non-targeted genes, fully custom panels perform well as shown by targeting metrics that measure how accurately the targeted library reflects gene expression information from the corresponding WTA library (Figure 7B-D).

Samples that did not meet the optimal performance threshold for fraction targeted genes with \geq 80% UMI recovery (Figure 7D) often had low expression of panel genes. However, these samples were able to achieve the optimal performance threshold with additional sequencing of the targeted sample. Figure 7E also shows fraction targeted genes with \geq 80% UMI recovery, but after discarding non-panel reads from the targeted library and downsampling the targeted library at two-fold the number of on target panel reads in the parent WTA library. By considering only the on target panel reads in the targeted library, the fraction targeted genes with \geq 80% UMI recovery metric is decoupled from reads mapped confidently to the targeted transcriptome.

The amount of additional sequencing used to make the comparison shown in Figure 7E often represents, at most, a few hundred additional reads per cell. For example, the values in Figure 7D assess the targeted library at two-fold the number of on target panel reads in the parent WTA library as described in the Methods. This equates to 100 read pairs per cell for a fully custom panel where expression of panel genes in a WTA library at 50,000 read pairs per cell is 0.1% (0.001 x 50,000 x 2 = 100). For the sample with the lowest reads mapped to the targeted transcriptome in this data set (36%), the parent-targeted comparison is made considering only reads that map to panel genes from the targeted library (Figure 7E).

The targeted library is assessed at 278 read pairs per cell (100/0.36 = 277.7) and now achieves expected performance.

Targeting Metric	Performance Threshold	% Pass	Mean ± stdev
Reads Mapped Confidently to the Targeted Transcriptome	80% 70%	77%	79±12%
Fraction Targeted Genes with ≥80% UMI Recovery	0.9	91%	0.95±0.05
Targeted Total UMI Recovery	0.8	100%	1.12±0.19
Pseudo-bulk targeted UMI R²	0.9	100%	0.98±0.03
Cell calling concordance	± 25% of parent WTA library detected	88%	0.91±0.14

Table 5. Percentage of fully custom panels that meet optimal performance thresholds. Sample statistics are computed over all sample-panel combinations (N=91). Samples were combined into pulldown reactions to maintain similar expression levels for targeted genes based on the percentage of library reads in the parent WTA library mapped to the respective panel. The 80% performance threshold (red) applies to Single Cell 3' Gene Expression libraries, while the 70% performance threshold (blue) applies to Single Cell 5' Gene Expression libraries.


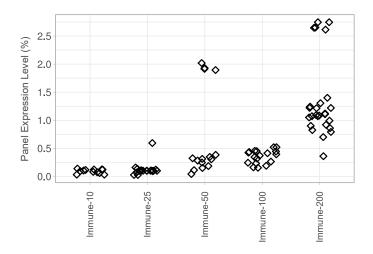


Figure 6. Expression level of a fully custom panel. Each data point represents a single WTA library. Panel Expression Level is the sum of the fraction of reads in the parent WTA library across targeted genes in a given panel.

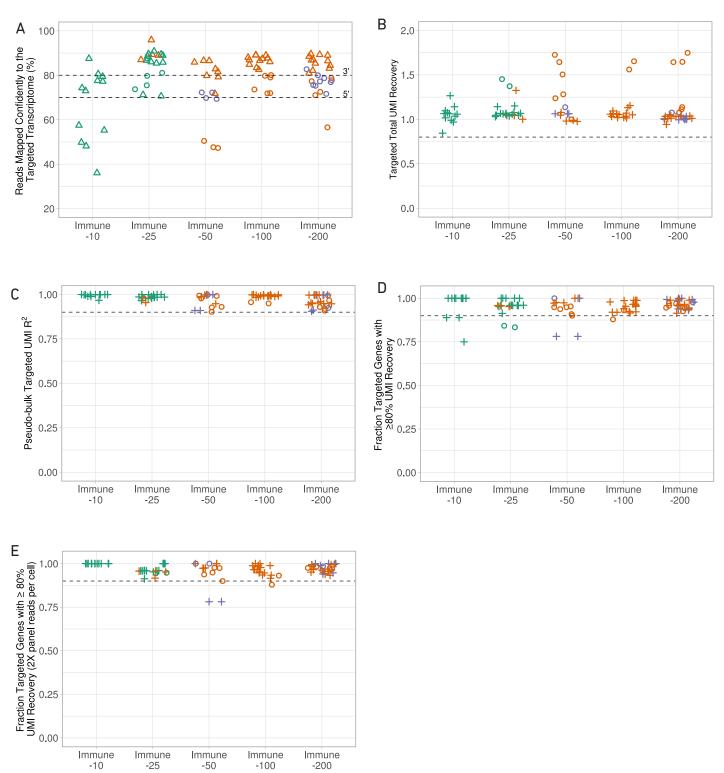

Targeting Metrics for Fully Custom Panels

Figure 7. Targeting metrics for fully custom panels. Each data point represents a single Targeted Gene Expression library that is colored for expression level: 0.10% (green), 0.25-1% (orange), and 2-4% (purple). For panels B-E, the targeted parent comparison was made using either 20,000 (circle) or 50,000 (plus sign) read pairs per cell from the parent WTA library. A. Reads mapped confidently to the targeted transcriptome. Sample data are grouped by the fully custom panel used. Dotted lines indicate threshold for optimal performance. Triangles indicate Chromium Single Cell 3' Gene Expression libraries and circles indicate Chromium Single Cell 5' Gene Expression libraries. B. Targeted Total UMI recovery. C. Pseudo-bulk targeted UMI R². D. Fraction of genes with UMI recovery \geq 80%. E. Same as (D) but assessed with a greater sequencing depth in the targeted sample to compensate for lower Reads Mapped Confidently to the Targeted Transcriptome. Refer to Results text for more information.

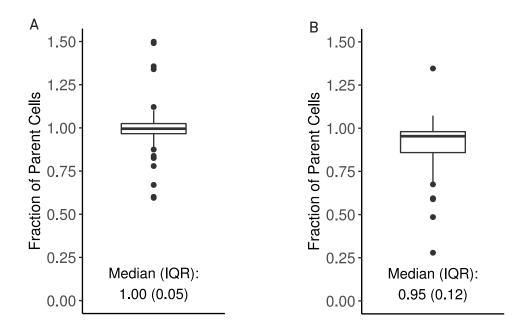


Figure 8. Cell calling concordance for add-on panels (A) and fully custom panels (B). Median and interquartile range (IQR) are indicated.

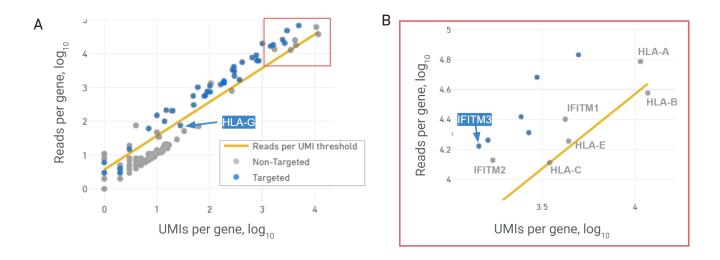


Figure 9. Target Enrichment plot from the Cell Ranger Web Summary for the Immune-50 custom panel. Panel (B) is a zoomed in view of the red highlighted area in (A).

Discussion

10x Genomics pre-designed panels supply a comprehensive set of genes relevant to several research areas. If additional genes are required, the Custom Panel Designer allows the design of add-on panels to supplement pre-designed content. Supplementing a pre-designed panel with add-on genes is a robust and cost-effective way to perform a targeted experiment.

Fully custom panels may vary in their number of genes, fraction of reads from targeted genes in the parent WTA library, and the expression of panel genes in cell subpopulations in a sample. Fully custom panels may be required when there is minimal overlap between the desired gene set and the genes in a pre-designed panel, or when a smaller panel (less than 1,000 genes) is desired.

Whether designing add-on panels or fully custom panels, consider the following:

Expression level of genes: Inclusion of highly expressed genes that are present in many cells in the sample will result in greater sequencing read pair requirements for the targeted sample. The Cell Ranger targeted-depth subcommand can help estimate the sequencing requirement for a custom panel. For instructions on running this script, refer to the Targeted Depth section of the 10x Genomics Support website. 10x Genomics recommends that the panel genes account for at least 0.1% of the WTA reads for optimal performance.

Differentially expressed genes that define a cell type: Choose differentially expressed genes that define cell types of interest to ensure that these cell types can be detected by the cell calling algorithm and differentiated during analysis. Cell calling should be robust, as shown in Figure 8 for this data set.

Presence of high-expression homologs: The Targeted Gene Expression workflow is sensitive to the sequence similarity of gene families. The inclusion of one gene in a homologous family can result in high non-targeted gene enrichment due to bait hybridization to family members. High expression homologs can consume baits and result in a lower relative enrichment of the desired gene. Conversely, the inclusion of baits for related members of a gene family that are not expressed may result in higher than expected UMI recovery for the gene family. Figure 9 shows an example of this phenomenon occurring in the Immune-50 fully custom panel. The panel gene IFITM3 shows good enrichment, but the related genes IFITM1 and IFITM2 that are also expressed in this cell are enriched despite not being included on the panel. This issue is also observed with the panel gene HLA-G and its related (but non-panel) genes HLA-A, HLA-B, HLA-C, and HLA-E. Presence of high-expression homologs decreases the fraction of reads mapping to the targeted transcriptome, and competition for baits may result in a lower relative enrichment of the desired gene.

Panel bait format (applicable to both fully custom and addon panels): Bait sequences obtained from the 10x Genomics Custom Panel Designer are synthesized through IDT or another oligo provider. Selecting the appropriate panel for the application is important for best results and for minimizing cost. 10x Genomics has tested two different panel formats from IDT: xGen Lockdown Probes and NGS Discovery Pools. xGen Lockdown Probes are individually synthesized and undergo per-bait quality control, while NGS Discovery Pools do not undergo the same level of quality control but are offered at a reduced cost. Although NGS Discovery Pools are appropriate for many applications, some applications require panels that have undergone per-bait quality control. In this data set, NGS Discovery Pools were compatible with both add-on panels and fully custom panels, but exhibited some bait dropout that was not detected when using xGen Lockdown Probes. Variability in the actual concentration of NGS Discovery Pool baits ocassionally resulted in over-representation of reads mapping to the add-on genes in the targeted library. This is illustrated in Figure 5, where a small concentration difference between add-on baits and pre-designed panel baits for the Pan-Cancer panel with Immune-Low add-on panel resulted in a higher proportion of reads mapping to add-on genes. This also led to an increased recovery of UMIs from add-on genes as compared to pre-designed panel genes.

IDT has discontinued these panel formats and now offers three different panel formats: xGen Custom Hyb Panels - Accel, xGen Custom Hyb Panels, and xGen Custom Hyb Panels - Production. All three formats are compatible with the Targeted Gene Expression Assay. xGen Custom Hyb Panel - Accel offers faster turnaround time and lower cost. The xGen Custom Hyb Panel includes an xGen NGS functional test report. The xGen Custom Hyb Panel - Production further includes quality documentation.

Conclusion

As shown by the metrics evaluated in this Technical Note, the 10x Genomics Custom Panel Designer may be used to generate add-on panels to supplement pre-designed panels or generate fully custom panels. These custom panels deliver a high recovery of unique molecules and efficient targeting of panel genes over a variety of sample and tissue types.

References

- Targeted Gene Expression Reagent Kits User Guide (CG000293)
- Targeted Gene Expression Pre-designed Panel Performance Metrics (CG000345)

Document Number	CG000349	

Document Revision Summary

Document Number	0000347
Title	Targted Gene Expression Custom Panel Performance Metrics and Best Practices
Revision	Rev B to Rev C
Revision Date	November 2022
General Changes	Updated for general minor consistency of language and terms throughout

Appendix

Panel							Genes						
Cancer-High	CALR	EIF	4A2	H3F3A	HNRNPK	MARCKSL	1 MCL	.1	RPL5	SF1	ZFP36L	.1 ZFF	P36L2
Cancer-Low	AR	CACN	A2D1	ERBB4	FZD9	HIST2H3	D LAM	C3 L	EFTY2	MPO	NKX2-	1 RPS	66KA6
Immune-High	ANP32B	AN	XA1	CD63	CD99	DUSP1	HMG	B1 L(GALS1	NAMPT	S100A1	0 SL0	25A3
Immune-Low	CES1	CLE	C4D	CXCR6	IGHD	KIR2DL3	KLR	F1	LCN2	MASP2	S100A	7 ZN	F683
Immune-10	BIN2	CC	L3	CXCL5	GNLY	HLA-DME	IGHO	52 N	BAN3	RORA	TOLLIF	, TR	EM2
Immune-25	BCL11B FCGR2B	BTLA IFIT2	CD160 INPP5D	CD33 ITGA1	CD36 ITGA5	CD37 LAMP3	CD46 MZB1	CD5 PBK	CD55 SH2D1B	CHI3L1 SLAMF6	CLEC4E SLC25A37	CX3CR1 WDHD1	DOCK8
Immune-50	ABCG2 CFP IGHM PYCARD	ADGRG3 CLEC6A IL12RB1 RUNX3	ADORA2A CR2 IL32 S100A12	AIRE CX3CR1 JAM3 SELL	C1QA CXCL10 LAMP2 SLC7A7	C4B CYBB LAMP3 SMPD3	CCL27 ENTPD1 MAF TAL1	CCL4 FCN1 MAGEA1 TNFSF11	CD180 FCRL2 MASP2 TNFSF18	CD2 GNLY MME VMO1	CD28 HLA-DMB MSH4 ZBED2	CD69 HLA-G NCF4	CD9 IFITM3 PI3
Immune-100	AICDA CAMP CD72 DOCK8 IL18 LILRB2 PSEN2 TLR10	ALCAM CASP5 CD9 EGR2 IL1RL2 LRRC32 RIPOR2 TLR5	ANP32B CCL18 CFI ENG IL27 LY9 RPN2 TLR7	AOC3 CCL20 CFP ENTPD1 IRGM MAF S100A8 TLR9	AQP9 CCL23 CHST10 ERAL1 ISG15 MAPK11 S100B TNFRSF11B	ATF2 CCL26 CLEC10A FAM30A ITCH MS4A2 SBN02 USP9Y	ATG5 CD1A CLEC4C FCGR3B ITGAX MSH5 SIGIRR VEGFB	ATG7 CD1B CN0T2 FUT7 ITGB1 NCAM1 SPANXB1 VM01	AZU1 CD247 CR2 GIMAP5 JAM3 NCR1 SPN YTHDF2	BIN2 CD46 CTRC HLA-DQA1 KLRC1 PNOC SPOCK2	BLK CD47 CXCL14 IGBP1 KLRD1 PPBP TFEB	BORA CD52 CXCL16 IGHM LCN2 PRF1 TICAM2	C1S CD55 CXCL IL17R LILRA PSEN TIRAI
Immune-200	A2M B3GAT1 CCL4 CD37 CLEC10A EBF1 FICD IFITM1 IL21R KLRG1 MARC0 PIK3AP1 S100A10 SLC25A3 TLR10 USP9Y	ABCB1 BCL11B CCR10 CD3E CMTM2 ECSIT FRYL IFNL1 IL2RG LILRA1 MCAM PNOC S100A8 SPA17 TLR5 VAV1	ABCB11 BTN3A1 CCR4 CD3G CSF1 EGR2 FUT4 IGHA1 IL9R LILRA4 MGST1 POU2AF1 S1PR1 S1PR1 SPINK5 TMEFF2 VSIG4	ABCC2 C1QA CCR9 CD40LG CSF2RB F13A1 GNLY IGHG3 IRF7 LILRB1 MNX1 PPBP SBN02 SP0CK2 TNFRSF11A	ADA C1QB CD160 CD46 CTSH F2RL1 GTF3C1 IGHG4 ITGA4 LILRB2 MSH4 PRAME SELL ST6GAL1 TNFRSF12A	ADGRG3 C7 CD180 CD5 CXCL1 F5 HAMP IGHM ITGB1 LRBA MZB1 PSMB7 SEMA7A STAT2 TNFRSF138	ANP32B CAMP CD1B CD52 CXCL10 FCAR HDC IGKC JAM3 LTB NAMPT REPS1 SH2B2 STAT6 TNFRSF13	IKZF2 JAML LTBR NCF4 RGS1 SH2D1B SYT17	CASP5 CD207 CD8B CXCL9 FCER1G I HLA-DQA1 IL17B KIR2DL3 LTK NCR3 RIPK2 SIGIRR TAP1	IL18R1 KIR3DL1 LY9 NOD1 RNASE2 SIGLEC1 TFEB	IL1RL1 KIR3DL2 MAGEA12 NOD2 RORA SKAP1 TGFBI	DOCK8 FCN1 IFIH1 IL1RL2 KLRC3 MAP3K7 NUP107 RRAD SLAMF1 THBD	ATG7 CCL3L1 CD36 CHIT1 DUSP1 FCRL2 IFIT2 IL2 KLRC4 MAPK14 PDIA6 RUNX3 SLAMF6 TICAM2 TRDC

Table 6. Genes used in add-on panels and fully custom panels.

Sample Sample A375 Cell Lines Jurkat/Raji Mixture Peripheral Blood Mononuclear Cells/Jurkat/GM12878 Mixture Gastroesophageal Glioblastoma Hodgkin's Lymphoma Lymphadenopathy Mucosa-associated Lymphoid Tissue Lymphoma Ovarian Cancer Peripheral Blood Mononuclear Cells		
Cell Lines Jurkat/Raji Mixture Peripheral Blood Mononuclear Cells/Jurkat/GM12878 Mixture Gastroesophageal Glioblastoma Image: Cell Science Scien		Sample
Peripheral Blood Mononuclear Cells/Jurkat/GM12878 Mixture Gastroesophageal Glioblastoma Glioblastoma Hodgkin's Lymphoma Lymphadenopathy Glioblastoma Mucosa-associated Lymphoid Tissue Lymphoma Glioblastoma		A375
Primary Cells Gastroesophageal Glioblastoma Upphadenopathy Aucosa-associated Lymphoid Tissue Lymphoma Aucosa-associated Lymphoma Upphoma Build Autophoma Upphoma Autophoma Ovarian Cancer	Cell Lines	Jurkat/Raji Mixture
Glioblastoma Hodgkin's Lymphoma Lymphadenopathy Mucosa-associated Lymphoid Tissue Lymphoma Multiple Myeloma Ovarian Cancer		Peripheral Blood Mononuclear Cells/Jurkat/GM12878 Mixture
Hodgkin's Lymphoma Lymphadenopathy Mucosa-associated Lymphoid Tissue Lymphoma Multiple Myeloma Ovarian Cancer		Gastroesophageal
Lymphadenopathy Mucosa-associated Lymphoid Tissue Lymphoma Multiple Myeloma Ovarian Cancer		Glioblastoma
Primary Cells A Constraint of Constraint o		Hodgkin's Lymphoma
Primary Cells Lymphoma Multiple Myeloma Ovarian Cancer		Lymphadenopathy
Primary Cells Multiple Myeloma Ovarian Cancer		Mucosa-associated Lymphoid Tissue
Multiple Myeloma Ovarian Cancer	Drimony Colle	Lymphoma
	Primary Cells	Multiple Myeloma
Peripheral Blood Mononuclear Cells		Ovarian Cancer
		Peripheral Blood Mononuclear Cells
Peripheral Blood Mononuclear Cells (Lupus)		Peripheral Blood Mononuclear Cells (Lupus)
Peripheral Blood Mononuclear Cells (Multiple Myeloma)		Peripheral Blood Mononuclear Cells (Multiple Myeloma)
Primary Bone Marrow Mononuclear Cells		Primary Bone Marrow Mononuclear Cells

Table 7. Human samples used for add-on and fully custom panel experiments.

© 2022 10x Genomics, Inc. (10x Genomics). All rights reserved. Duplication and/or reproduction of all or any portion of this document without the express written consent of 10x Genomics, is strictly forbidden. Nothing contained herein shall constitute any warranty, express or implied, as to the performance of any products described herein. Any and all warranties applicable to any products are set forth in the applicable terms and conditions of sale accompanying the purchase of such product. 10x Genomics provides no warranty and hereby disclaims any and all warranties as to the use of any third-party products or protocols described herein. The use of products described herein is subject to certain restrictions as set forth in the applicable terms and conditions of sale accompanying the purchase of such product. Non-exhaustive list of 10x Genomics' marks, many of which are registered in the United States and other countries can be viewed at: www.10xgenomics. com/trademarks. 10x Genomics may refer to the products or services offered by other companies by their brand name or company name solely for clarity, and does not claim any rights in those third-party marks or names. 10x Genomics products may be covered by one or more of the patents as indicated at: www.10xgenomics. com/patents. All products and services described herein are intended FOR RESEARCH USE ONLY and NOT FOR USE IN DIAGNOSTIC PROCEDURES.

The use of 10x Genomics products in practicing the methods set forth herein has not been validated by 10x Genomics, and such non-validated use is NOT COVERED BY 10X GENOMICS STANDARD WARRANTY, AND 10X GENOMICS HEREBY DISCLAIMS ANY AND ALL WARRANTIES FOR SUCH USE. Nothing in this document should be construed as altering, waiving or amending in any manner 10x Genomics terms and conditions of sale for the Chromium Controller or the Chromium Single Cell Controller, consumables or software, including without limitation such terms and conditions relating to certain use restrictions, limited license, warranty and limitation of liability, and nothing in this document shall be deemed to be Documentation, as that term is set forth in such terms and conditions of sale. Nothing in this document shall be construed as any representation by 10x Genomics that it currently or will at any time in the future offer or in any way support any application set forth herein. Contact: support@10xgenomics.com 10x Genomics 6230 Stoneridge Mall Road Pleasanton, CA 94588 USA

