

CG000352 Rev N

Chromium Connect Software Release Note

FOR USE WITH

Chromium Connect, PN-1000171

Chromium Connect, PN-1000180

Notices

Document Number

CG000352 • Rev N

Legal Notices

© 2023 10x Genomics, Inc. (10x Genomics). All rights reserved. Duplication and/or reproduction of all or any portion of this document without the express written consent of 10x Genomics, is strictly forbidden. Nothing contained herein shall constitute any warranty, express or implied, as to the performance of any products described herein. Any and all warranties applicable to any products are set forth in the applicable terms and conditions of sale accompanying the purchase of such product. 10x Genomics provides no warranty and hereby disclaims any and all warranties as to the use of any third-party products or protocols described herein. The use of products described herein is subject to certain restrictions as set forth in the applicable terms and conditions of sale accompanying the purchase of such product. A non-exhaustive list of 10x Genomics' marks, many of which are registered in the United States and other countries can be viewed at: www.10xgenomics.com/trademarks. 10x Genomics may refer to the products or services offered by other companies by their brand name or company name solely for clarity, and does not claim any rights in those third party marks or names. 10x Genomics products may be covered by one or more of the patents as indicated at: www.10xgenomics.com/patents. All products and services described herein are intended FOR RESEARCH USE ONLY and NOT FOR USE IN DIAGNOSTIC PROCEDURES.

Instrument & Licensed Software Updates Warranties

Updates to existing Instruments and Licensed Software may be required to enable customers to use new or existing products. In the event of an Instrument failure resulting from an update, such failed Instrument will be replaced or repaired in accordance with the 10x Genomics Limited Warranty, Assurance Plan or service agreement, only if such Instrument is covered by any of the foregoing at the time of such failure. Instruments not covered under a current 10x Genomics Limited Warranty, Assurance Plan or service agreement will not be replaced or repaired.

Support

Email: support@10xgenomics.com

10x Genomics

6230 Stoneridge Mall Road

Pleasanton, CA 94588 USA

Chromium Connect Software Release Note

Item	Description
For Use With	PN-1000171 Chromium Connect PN-1000180 Chromium Connect
Software Version	V2022.12.30
Software Release Date	December 2022
Compatibility	Chromium Connect Instrument
Features	<ol style="list-style-type: none">1. New manual barcode data entry popup screen2. All assay workflows now dispense a fixed amount of library volume for all channels3. New popup warning screen when selecting "Expert" or "Advanced" instruction levels4. New assay setup loading data from previously cancelled run during loading phase5. New remote support activation/deactivation screen
Bug Fixes & Changes	<ol style="list-style-type: none">6. HEPA fans activated during all maintenance runs7. Improved device connection error handling defects8. Improved assay run screen displays timers based upon variations in number of samples and thermal cycling

Previous Software Releases

Item	Description
Previous Software Version V2022.09.30	
Features	<ol style="list-style-type: none">1. New assay start notification to insert all carriers on the deck2. Network Connection Screen allows for network adapter display and option to disconnect from Wi-Fi
Bug Fixes & Changes	<ol style="list-style-type: none">1. User instructions to unload the deck incorrectly indicate to unload the chip from a carrier2. Error message incorrect for "Oil tube" Deck Check3. Import of CSV file defects

Previous Software Releases

Item	Description
	<p>Previous Software Version V2022.06.29</p> <p>Additional specifics:</p> <ul style="list-style-type: none"> Chromium Automated Controller V4.1.0 (FW 4.0.1, TEC FW 1.5.0) Chromium Connect User Interface V1.31.5.6 Chromium Connect Automated Assays: Build 669 Hamilton Nimbus8 Software Version 2.1.4 rev B SP2 Hamilton Nimbus8 Firmware Version 2.1.3 SP1 <p>Software Released: June 2022</p> <p>Features</p> <ol style="list-style-type: none"> 1. User Guides available on-instrument via 'Help' menu 2. User may select previous run setup via search function, use variables from previous run design 3. Unavailable options grayed out on user interface if component is not present (i.e. HEPA fan not installed) 4. cDNA input volumes user-controlled in library preparation setup 5. Add local drives other than C:\ <p>Bug Fixes & Changes</p> <ol style="list-style-type: none"> 1. Import of CSV files do not transfer volume information to Automated Single Cell 5' GEX Library Constructions setup
	<p>Previous Software Version V2022.03.30</p> <p>Additional Specifics:</p> <ul style="list-style-type: none"> Chromium Automated Controller V4.1.0 (FW 4.0.1, TEC FW 1.5.0) Chromium Connect User Interface V1.30.3.6 Chromium Connect Automated Assays: Build 606 Hamilton Nimbus8 Software Version 2.1.4 rev B SP2 Hamilton Nimbus8 Firmware Version 2.1.3 SP1 <p>Software Released: March 2022</p> <p>Features</p> <ol style="list-style-type: none"> 1. New Single Cell 5' Feature Barcode Library Construction Kit 2. Add "Resuspend Dynabeads" information to Load Sample Plate screen(s) 3. Disable/enable Assay Kits 4. Enhancements to instrument user interface Run screen information display 5. User interface Run Report list updated to include color codes to specify status (Red - ERROR; Yellow - CANCEL/ABORT; Green - SUCCESS) 6. New user interface screen image (10x Genomics Logo) viewer and image exporter <p>Bug Fixes & Changes</p> <ol style="list-style-type: none"> 1. Power cycling instrument failure fix to allow user to run assays 2. Automated Single Cell 3' and 5' Library Construction assay failure to run immediately after software installation 3. User interface invalid numeric input data does not properly display text and error indicator
	<p>Previous Software Version V2022.01.14</p> <p>Software Released: January 2022</p> <p>Bug Fixes & Changes</p> <ol style="list-style-type: none"> 1. Power cycling instrument failure fix to allow user to run assays 2. Automated Single Cell 3' and 5' Library Construction assay failure to run immediately after software installation 3. User interface invalid numeric input data does not properly display text and error indicator

Previous Software Releases

Previous Software Version V2021.12.14

Released: December 2021

Features

1. New Chromium Next GEM Automated Single Cell 3' v3.1 & 5' v2 cDNA and Library Construction Kit and __workflow compatibility
2. Automated universal Library Construction workflow
3. Increase in library pooling capability (8 to 32) and the number of pools to populate (1 to 6)
4. Comfort Lid quantity reduction for assays

Bug Fixes & Changes

1. Email notification improvements
2. Additional error recovery
3. LIMS workbook update for duplicate Sample Index entries

Previous Software Version V2021.09.28

Released: September 2021

Features

1. Chromium Next GEM Single Cell 3' v3.1 Gene Expression Automated workflow supports Dual Indexing
2. Assay run Pause button disables during critical sections of the workflow
3. Instrument controlled HEPA operation availability
4. HEPA Fans turn ON after Setup phase of workflow
5. New status bar indicators for network and remote server connection

Bug Fixes & Changes

1. Improvements in user interactive error recovery situations

Previous Software Version V2021.05.27

Released: May 2021

Features

1. Updated for Chromium Next GEM Single Cell 5' v2 assay
2. Prevents selection of inaccessible Gel Bead wells
3. New Run Setup Confirmation screen

Bug Fixes & Changes

1. Invalidates improperly formatted email address entered by user

Previous Software Releases

Previous Software Version V2021.03.31 Released: April 2021

Features

1. Instrument login status
2. Pause Button warning to indicate that the run may abort if instrument is paused
3. Updated LIMS import to assay template files

Bug Fixes & Changes

1. Anti-droplet Control (ADC) channel failure terminates assay
2. Troubleshooting package failure to include all data
3. Data transfer handling of Software Upgrade program from a USB
4. Control Panel failure when executing Thermal Cycler "Run protocol"
5. Incorrect recovery dialog appears when a Carrier "break" sensor error is detected
6. If a run is cancelled, screen prompts the user to remove Chip from the Automated Chromium Controller
7. Messaging update to Tip pickup failure error handling

Previous Software Version V2020.10.07 Released: October 2020

Features

1. Loading of pipette tip racks scales with number of samples
2. LIMS Output to include a run result field (i.e. SUCCESS | FAILURE | ABORT)
3. Screen display to indicate status of Chromium Connect Software update

Bug Fixes & Changes

1. HEPA light remains ON during a run
2. cDNA no longer aliquoted to TS-1 when mid-run QC is skipped
3. Updated maintenance screens
4. Pre-run deck check ensures Waste Bar is present
5. Repetitive prompt to refill pipette tips removed
6. Wraparound selection of SI wells now compatible with the assay
7. Secondary Assay transition corrected to proceed with beginning or end of run inputs

Previous Software Releases

Previous Software Version V2020.07.31
Released: August 2020

Features

1. Chromium Automated Controller Firmware update (v4.1.0) integrated into Chromium Connect Software update
2. Software integration for HEPA CAP hood functionality
3. Capability to export barcode information into LIMS and run audit logs
4. Added LIMS compatibility for qPCR and Pooling

Bug Fixes & Changes

1. Updated User Interface screens for general consistency of language and formatting
 2. Added sound notification for GEM Check Timer
 3. Minor bug fixes to user interface and barcode scanning
 4. Added ability to remove network drives in menu settings
 5. Enabled data purging off storage databases
 6. Easier access to waste bin during maintenance
 7. Pooling assay accepts sample input
-