

User Guide | CG000416 | Rev D

Chromium Next GEM Single Cell 3' HT Reagent Kits v3.1 (Dual Index)

For use with:

Chromium Next GEM Single Cell 3' HT Kit v3.1 48 rxns PN-1000348 | 8 rxns PN-1000370

Chromium Next GEM Chip M Single Cell Kit^{*} 80 rxns PN-1000349 | 16 rxns PN-1000371 (*Included with Chromium Next GEM Single Cell 3' HT Kit v3.1; 16 rxn kit can also be ordered separately)

Dual Index Kit TT Set A 96 rxns PN-1000215

Notices

Document Number

CG000416 | Rev D

Legal Notices

© 2022 10x Genomics, Inc. (10x Genomics). All rights reserved. Duplication and/or reproduction of all or any portion of this document without the express written consent of 10x Genomics, is strictly forbidden. Nothing contained herein shall constitute any warranty, express or implied, as to the performance of anyproducts described herein. Any and all warranties applicable to any products are set forth in the applicable terms and conditions of sale accompanying the purchase of such product. 10x Genomics provides no warranty and hereby disclaims any and all warranties as to the use of any third-party products or protocols described herein. The use of products described herein is subject to certain restrictions as set forth in the applicable terms and conditions of sale accompanying the purchase of such product. A non-exhaustive list of 10x Genomics' marks, many of which are registered in the United States and other countries can be viewed at: www.10xgenomics.com/trademarks. 10x Genomics may refer to the products or services offered by other companies by their brand name or company name solely for clarity, and does not claim any rights in those third-party marks or names. 10x Genomics products may be covered by one or more of the patents as indicated at: www.10xgenomics.com/patents. The use of products described herein is subject to 10x Genomics Terms and Conditions of Sale, available at www.10xgenomics.com/legal-notices, or such other terms that have been agreed to in writing between 10x Genomics and user. All products and services described herein are intended FOR RESEARCH USE ONLY and NOT FOR USE IN DIAGNOSTIC PROCEDURES.

Instrument & Licensed Software Updates Warranties

Updates to existing Instruments and Licensed Software may be required to enable customers to use new or existing products. In the event of an Instrument failure resulting from an update, such failed Instrument will be replaced or repaired in accordance with the 10x Limited Warranty, Assurance Plan or service agreement, only if such Instrument is covered by any of the foregoing at the time of such failure. Instruments not covered under a current 10x Limited Warranty, Assurance Plan or service agreement will not be replaced or repaired.

Support

Email: support@10xgenomics.com 10x Genomics 6230 Stoneridge Mall Road Pleasanton, CA

Document Revision Summary

Document Number

CG000416 | Rev D

Title

Chromium Next GEM Single Cell 3' HT Reagent Kits v3.1 (Dual Index) DOCUMENT SUBTITLE

Revision

Rev C to Rev D

Revision Date

September 26, 2022

Specific Changes

- Updated Thermal Cycler recommendations.
- Updated to include correct storage time (for up to 4 weeks) at step 2.3n.

General Changes

Updated for general minor consistency of language and terms throughout.

Table of Contents

Introduction

| Chromium Next GEM Single Cell 3' HT Reagent Kits v3.1 | 7 |
|---|----|
| 10x Genomics Accessories | 11 |
| Recommended Thermal Cyclers | 11 |
| Additional Kits, Reagents & Equipment | 12 |
| Protocol Steps & Timing | 16 |
| Stepwise Objectives | 17 |

Tips & Best Practices

Step 1: GEM Generation and Barcoding

| 1.0 Get Started | 37 |
|-----------------------------------|----|
| 1.1 Prepare Master Mix | 39 |
| Assemble Chromium Next GEM Chip M | 40 |
| Sample Loading Guidelines | 41 |
| 1.2 Load Chromium Next GEM Chip M | 43 |
| 1.3 Run Chromium X | 45 |
| GEM Transfer Overview | 46 |
| 1.4 Transfer GEMs | 47 |
| 1.5 GEM-RT Incubation | 49 |
| | |

Step 2: Post GEM-RT Cleanup & cDNA Amplification

| 2.0 Get Started | 51 |
|---|----|
| 2.1 Post GEM-RT Cleanup – Dynabeads | 53 |
| 2.2 cDNA Amplification | 55 |
| 2.3 cDNA Cleanup – SPRIselect | 57 |
| 2.4 Post cDNA Amplification QC & Quantification | 58 |

Step 3: 3' Gene Expression Library Construction

| 3.0 Get Started | 61 |
|--|----|
| Step Overview (Step 3.1d) | 63 |
| 3.1 Fragmentation, End Repair & A-tailing | 64 |
| 3.2 Post Fragmentation, End Repair & A-tailing Double Sided – SPRIselect | 65 |
| 3.3 Adaptor Ligation | 66 |
| 3.4 Post Ligation Cleanup – SPRIselect | 67 |

Table of Contents

| 3.5 Sample Index PCR | 68 |
|--|----|
| 3.6 Post Sample Index PCR Double Sided Size Selection – SPRIselect | 69 |
| 3.7 Post Library Construction QC | 70 |
| | |
| Step 4: Sequencing | |
| Sequencing Libraries | 72 |
| Illumina Sequencer Compatibility | 73 |
| Sample Indices | 73 |
| 3' Gene Expression Library Sequencing Parameters | 74 |
| Library Loading | 75 |
| Library Pooling | 75 |
| Data Analysis and Visualization | 76 |
| | |
| Troubleshooting | |
| GEMs | 78 |
| Chromium X Series Errors | 81 |
| Appendix | |
| Post Library Construction Quantification | 83 |
| | 84 |
| Agilent TapeStation Traces | 85 |
| LabChip Traces | |
| Oligonucleotide Sequences | 86 |
| | |



Introduction

| Chromium Next GEM Single Cell 3' HT Reagent Kits v3.1 | 7 |
|---|----|
| 10x Genomics Accessories | 11 |
| Recommended Thermal Cyclers | 11 |
| Additional Kits, Reagents & Equipment | 12 |
| Protocol Steps & Timing | 16 |
| Stepwise Objectives | 17 |

Chromium Next GEM Single Cell 3' HT Reagent Kits v3.1

Refer to SDS for handling and disposal information

Chromium Next GEM Single Cell 3' HT Kit v3.1, 48 rxns PN-1000348

| Sing 48 r | omium Next GEM gle Cell 3' HT GEM Kit xns, PN-1000351 'e at -20°C | v3.1 | | 48 r | ary Construction Kit xns, PN-1000352 re at -20°C | | |
|--------------|--|------|-----------------|------|--|---|---------|
| | | # | PN | | | # | PN |
| | RT Reagent B | 3 | 2000435 | | Fragmentation Enzyme | 3 | 2000090 |
| | RT Enzyme C | 3 | 2000436 | | Fragmentation Buffer | 3 | 2000091 |
| | Template Switch Oligo | 6 | 3000228 | | Ligation Buffer | 3 | 2000092 |
| 0 | Reducing Agent B | 3 | 2000087 | | DNA Ligase | 3 | 220110 |
| | Cleanup Buffer | 3 | 2000438 | | Adaptor Oligos | 3 | 2000094 |
| | cDNA Primers | 3 | 2000439 | 0 | Amp Mix | 3 | 2000047 |
| 0 | Amp Mix | 3 | 2000440 | | | | |
| | | | 10x genomics | | | | 10x |

| Chromium Next GEM Single Cell 3' HT Gel Bead Kit v3.1 48 rxns, PN-1000350 Store at -80°C | | | | | | |
|---|---|---------|--|--|--|--|
| | # | PN | | | | |
| Single Cell 3' HT Gel Beads v3.1 | 3 | 2000443 | | | | |
| | | 10× | | | | |

| Dynabeads [™] MyOne [™] SILANE PN-2000048 Store at 4°C | | | |
|--|---|---------|--|
| | # | PN | |
| Dynabeads MyOne SILANE | 6 | 2000048 | |
| | | | |


Chromium Next GEM Chip M Single Cell Kit, 80 rxns PN-1000349

| Part | Chromium Partitioning Oil Store at ambient temperature | | Rec | omium covery Agent re at ambient tem | iperat | ture | |
|------|---|---|--------|---|----------------|------|---------|
| | | # | PN | | | # | PN |
| | Partitioning Oil | 5 | 220088 | 0 | Recovery Agent | 5 | 2000434 |
| | | | | | | | |

| Chromium Next GEM Chip M & Gaskets Store at ambient temperature | | | |
|---|---|---------|-----|
| | # | PN | |
| Chromium Next GEM Chip M | 5 | 2000417 | |
| Chip Gasket, HT, 5-pack | 1 | 3000614 | |
| | | | 10× |

Chromium Next GEM Single Cell 3' HT Kit v3.1, 8 rxns PN-1000370

| Sing 8 rxr | omium Next GEM gle Cell 3' HT GEM Ki ns, PN-1000373 re at -20°C | t v3. 1 | I | | rary Construction Kit xns, PN-1000190 re at -20°C | | |
|---------------|--|----------------|---------|---|---|---|---------|
| | | # | PN | | | # | PN |
| | RT Reagent B | 1 | 2000165 | | Fragmentation Enzyme | 1 | 2000090 |
| | RT Enzyme C | 1 | 2000085 | | Fragmentation Buffer | 1 | 2000091 |
| | Template Switch Oligo | 1 | 3000228 | | Ligation Buffer | 1 | 2000092 |
| \bigcirc | Reducing Agent B | 1 | 2000087 | | DNA Ligase | 1 | 220110 |
| | Cleanup Buffer | 2 | 2000088 | | Adaptor Oligos | 1 | 2000094 |
| | cDNA Primers | 1 | 2000089 | 0 | Amp Mix | 1 | 2000047 |
| 0 | Amp Mix | 1 | 2000047 | | | | |
| | | | 10x | | | | 10x |

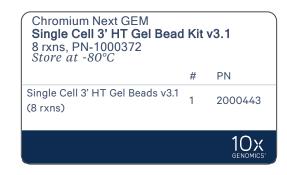

| Dynabeads[™] MyOne[™] SILANE PN-2000048 Store at 4°C | | | |
|---|---|---------|--|
| | # | PN | |
| Dynabeads MyOne SILANE | 2 | 2000048 | |

Chromium Next GEM Chip M Single Cell Kit, 16 rxns PN-1000371

| Chromium Partitioning Oil Store at ambient temperature | | | | Rec | omium covery Agent re at ambient tem | perat | ture | |
|---|------------------|---|--------|-----|---|----------------|------|---------|
| | | # | PN | | | | # | PN |
| | Partitioning Oil | 1 | 220088 | | 0 | Recovery Agent | 1 | 2000434 |
| | | | | J | | | | |

| Chromium Next GEM Chip M & Gaskets Store at ambient temperature | | | |
|---|---|---------|-----|
| | # | PN | |
| Chromium Next GEM Chip M | 1 | 2000417 | |
| Chip Gasket, HT, 2-pack | 1 | 3000656 | |
| | | | 10× |

Dual Index Kit TT Set A, 96 rxns PN-1000215

| Dual Index Kit TT Set A Store at -20°C | | Ň |
|--|---|---------|
| | # | PN |
| Dual Index Plate TT Set A | 1 | 3000431 |
| | | |

10x Genomics Accessories

| Product | Part Number | Part Number (Item) |
|---------------------------|-------------|--------------------|
| 10x Vortex Adapter | 120251 | 330002 |
| 10x Magnetic Separator HT | 1000394 | 2000431 |
| Chromium X Chip Holder | 1000393 | 3000598 |

Recommended Thermal Cyclers

The table below lists the thermal cyclers that have been validated by 10x Genomics. Thermal cyclers used must support uniform heating of 100 μl emulsion volumes

| Supplier | Description | Part Number |
|-----------------------------|---|---|
| BioRad | C1000 Touch Thermal Cycler with 96-Deep Well Reaction Module | 1851197 |
| Analytik Jena† | Biometra TAdvanced 96 SG | 846-x-070-241 (x=2 for 230 V; 4 for 115 V; 5 for 100 V, 50-60 Hz) |
| Eppendorf* | MasterCycler X50s | North America 6311000010 |
| | Mastercycler Pro (discontinued) | North America 950030010 International 6321 000.019 |
| Thermo Fisher Scientific | Veriti 96-Well Thermal Cycler | 4375786 |

Introduction

Additional Kits, Reagents & Equipment

The items in the table below have been validated by 10x Genomics and are highly recommended for 10x Genomics workflows, training, and system operations. Substituting materials may adversely affect system performance. This list does not include standard laboratory equipment, such as water baths, centrifuges, vortex mixers, pH meters, freezers, etc.

| Supplier | Description | Part Number (US) |
|------------------------------|---|------------------------|
| Plastics | | |
| Choose either Eppendorf, USA | A Scientific, or Thermo Fisher Scientific PCR 8-tube s | trips |
| Eppendorf | PCR Tubes 0.2 ml 8-tube strips | 951010022 |
| | DNA LoBind Tubes, 1.5 ml | 022431021 |
| | DNA LoBind Tubes, 2.0 ml | 022431048 |
| USA Scientific | TempAssure PCR 8-tube strip (alternate to Eppendorf or Thermo Fisher Scientific product) | 1402-4700 |
| Thermo Fisher Scientific | MicroAmp 8-Tube Strip, 0.2 ml (alternate to Eppendorf or USA Scientific product) | N8010580 |
| | MicroAmp 8 -Cap Strip, clear | N8010535 |
| Kits & Reagents | | |
| Thermo Fisher Scientific | Nuclease-free Water Low TE Buffer (10 mM Tris-HCl pH 8.0, 0.1 mM EDTA) | AM9937 12090-015 |
| Millipore Sigma | Ethanol, Pure (200 Proof, anhydrous) | E7023-500ML |
| Beckman Coulter | SPRIselect Reagent Kit | B23318 |
| Bio-Rad | 10% Tween 20 | 1662404 |
| Ricca Chemical Company | Glycerin (glycerol), 50% (v/v) Aqueous Solution | 3290-32 |
| Qiagen | Qiagen Buffer EB | 19086 |
| Equipment | | |
| VWR | Vortex Mixer Divided Polystyrene Reservoirs | 10153-838 41428-958 |
| Thermo Fisher Scientific | MYFUGE 12 Mini Centrifuge (alternatively, use any equivalent mini centrifuge) | C1012 |
| | lsotemp Drybath Incubators (when using 48 rxn reagent kits) (alternatively, use any equivalent drybath or waterbath) | 11-718-22Q |

| Supplier | Description | Part Number (US) |
|--------------------------------|---|-----------------------------------|
| Eppendorf | Eppendorf ThermoMixer C (when using 8 rxn reagent kits) | 5382000023 |
| | Eppendorf SmartBlock 1.5 ml, Thermoblock for 24 reaction vessel (alternatively, use a temperature-controlled Heat Block) | 5360000038 |
| Quantification & Quality Contr | ol | |
| Choose either Bioanalyzer, 2 | TapeStation, Fragment Analyzer, Labchip, or Qubit bas | sed on availability & preference. |
| Agilent | 2100 Bioanalyzer Laptop Bundle (discontinued) (Replacement 2100 Bioanalyzer Instrument/2100 Expert Laptop Bundle) | G2943CA G2939BA/G2953CA |
| | High Sensitivity DNA Kit | 5067-4626 |
| | 4200 TapeStation | G2991AA |
| | High Sensitivity D1000 ScreenTape/Reagents | 5067-5592/ 5067-5593 |
| | High Sensitivity D5000 ScreenTape/Reagents | 5067-5584/ 5067-5585 |
| Thermo Fisher Scientific | Qubit 4.0 Flourometer Qubit dsDNA HS Assay Kit | Q33226 Q32854 |
| Advanced Analytical | Fragment Analyzer Automated CE System - 12 cap | FSv2-CE2F |
| | Fragment Analyzer Automated CE System - 48/96 cap | FSv2-CE10F |
| | High Sensitivity NGS Fragment Analysis Kit | DNF-474 |
| PerkinElmer | LabChip GX Touch HT Nucleic Acid Analyzer DNA High Sensitivity Reagent Kit | CLS137031 CLS760672 |
| KAPA Biosystems | KAPA Library Quantification Kit for Illumina Platforms | KK4824 |

Recommended Pipette Tips

10x Genomics recommends using only validated emulsion-safe pipette tips for all Single Cell protocols. Rainin pipette tips have been extensively validated by 10x Genomics and are highly recommended for all single cell assays. If Rainin tips are unavailable, any of the listed alternate pipette tips validated by 10x Genomics may be used.

| Supplier | Description | Part Number (US) |
|---|--|------------------|
| Recommended Pipettes & Pipette tips | | |
| Rainin | Pipettes Pipet-Lite Multi Pipette L8-50XLS+ | 17013804 |
| | Pipet-Lite Multi Pipette L8-200XLS+ | 17013805 |
| | Pipet-Lite Multi Pipette L8-10XLS+ | 17013802 |
| | Pipet-Lite Multi Pipette L8-20XLS+ | 17013803 |
| | Pipet-Lite LTS Pipette L-2XLS+ | 17014393 |
| | Pipet-Lite LTS Pipette L-10XLS+ | 17014388 |
| | Pipet-Lite LTS Pipette L-20XLS+ | 17014392 |
| | Pipet-Lite LTS Pipette L-100XLS+ | 17014384 |
| | Pipet-Lite LTS Pipette L-200XLS+ | 17014391 |
| | Pipet-Lite LTS Pipette L-1000XLS+ | 17014382 |
| | Pipette Tips Tips LTS 200UL Filter RT-L200FLR | 30389240 |
| | Tips LTS 1ML Filter RT-L1000FLR | 30389213 |
| | Tips LTS 20UL Filter RT-L10FLR | 30389226 |
| Alternate Recommendations (If Rainin pipette tips are unavailable, | any of the listed pipette tips may be used) | |
| Eppendorf | Pipettes Eppendorf Research Plus, 8-channel, epT.I.P.S. Box, 0.5 – 10 μl | 3125000010 |
| | Eppendorf Research Plus, 8-channel, epT.I.P.S. Box, 10 – 100 µl | 3125000036 |
| | Eppendorf Research Plus, 8-channel, epT.I.P.S. Box, 30 – 300 µl | 3125000052 |
| | Eppendorf Research Plus, 1-channel, epT.I.P.S.® Box, 0.1 – 2.5 µl | 3123000012 |
| | Eppendorf Research Plus, 1-channel, epT.I.P.S.® Box, 0.5 – 10 μI | 3123000020 |
| | Eppendorf Research Plus, 1-channel, epT.I.P.S.® Box, 2 – 20 µl | 3123000039 |
| | Eppendorf Research Plus, 1-channel, epT.I.P.S.® Box, 2 – 200 μl | 3123000055 |
| | Eppendorf Research Plus, 1-channel, epT.I.P.S.® Box, 100 – 1000 µl | 3123000063 |

| Supplier | Description | Part Number (US) |
|----------|--|------------------|
| | Pipette Tips | |
| | (compatible with Eppendorf pipettes only) | |
| | ep Dualfilter T.I.P.S., 2-20 µl | 0030078535 |
| | ep Dualfilter T.I.P.S., 2-200 μl | 0030078551 |
| | ep Dualfilter T.I.P.S., 2-1,000 µl | 0030078578 |
| Labcon* | ZAP SLIK 20 μL Low Retention Aerosol Filter Pipet Tips for Rainin LTS | 4-1143-965-008 |
| | ZAP SLIK 200 μL Low Retention Aerosol Filter Pipet Tips for Rainin LTS | 4-1144-965-008 |
| | ZAP SLIK 1200 μL Low Retention Aerosol Filter Pipet Tips for Rainin LTS | 4-1145-965-008 |
| Biotix* | xTIP4 Racked Pipette Tips, Rainin LTS Pipette Compatible, 0.1-20 µl | 63300931 |
| | xTIP4 Racked Pipette Tips, Rainin LTS Pipette Compatible, 200 μI | 63300001 |
| | xTIP4 Racked Pipette Tips, Rainin LTS Pipette Compatible, 1200 µI | 63300004 |

*Compatible with Rainin pipettes

Protocol Steps & Timing

| Day | Steps | Timing | Stop & Store |
|-----|--|--|--|
| | Cell Preparation | | |
| 2 h | Dependent on Cell Type | ~1-1.5 h | |
| | Step 1 – GEM Generation & Barcoding | | |
| | Prepare Reaction Mix Load Chromium Next GEM Chip M Run the Chromium X | 20 min 10 min 18 min | |
| | 1.5Run the chromitan X1.4Transfer GEMs1.5GEM-RT Incubation | 3 min | 500 4°C ≤72 h or −20°C ≤1 week |
| 4 h | Step 2 – Post GEM-RT Cleanup & cDNA Amplification | | |
| | 2.1 Post GEM RT-Cleanup – Dynabead 2.2 cDNA Amplification 2.3 cDNA Cleanup – SPRIselect 2.4 cDNA QC & Quantification | 60 min 40 min 35 min 50 min | 510P 4°C ≤72 h or −20°C ≤1 week 510P 4°C ≤72 h −20°C ≤4 weeks |
| 6 h | Step 3 – 3' Gene Expression Library Construction | | |
| | 3.1 Fragmentation, End Repair & A-tailing 3.2 Post Fragmentation, End Repair & A-tailing Double Sided Size Selection – SPRIselect | 50 min 30 min | |
| 8 h | 3.3 Adaptor Ligation 3.4 Post Ligation Cleanup- SPRIselect 3.5 Sample Index PCR 3.6 Post Sample Index PCR Double Sided Size Selection- SPRIselect 3.7 Post Library Construction QC | 25 min 20 min 40 min 30 min 50 min | stop 4°C ≤72 h 5Top 4°C ≤72 h or -20°C long term |

😳 Stepwise Objectives


The Chromium Single Cell Gene Expression HT is a scalable microfluidic platform for 3' digital gene expression for profiling 2,000-20,000 individual cells per sample. A pool of ~3,500,000 10x Barcodes are sampled separately to index each cell's transcriptome. It is done by partitioning thousands of cells into nanoliter-scale Gel Beads-in-emulsion (GEMs), where all generated cDNA share a common 10x Barcode. Dual indexed libraries are generated and sequenced from the cDNA, and 10x Barcodes are used to associate individual reads back to the individual partitions.

This document outlines the high throughput (HT) protocol for generating Single Cell 3' Gene Expression dual index libraries from single cells.

Single Cell 3' HT v3.1 Gel Beads

In addition to the poly(dT) primer that enables the production of barcoded, full-length cDNA from poly-adenylated mRNA, the Single Cell 3' HT v3.1 Gel Beads also include two additional primer sequences (Capture Sequence 1 and Capture Sequence 2) that enable capture and priming of Feature Barcode technology compatible targets or analytes of interest.


Only the poly(dT) primers are used in this protocol for generating Single Cell 3' Gene Expression libraries.

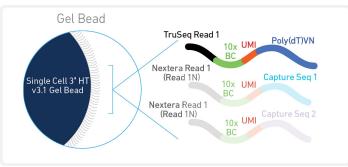




Step 1: GEM Generation & Barcoding

GEMs are generated by combining barcoded Gel Beads, a Master Mix containing cells, and Partitioning Oil onto Chromium Next GEM Chip M. To achieve single cell resolution, cells are delivered at a limiting dilution, such that the majority (~90-99%) of generated GEMs contain no cell, while the remainder largely contain a single cell.

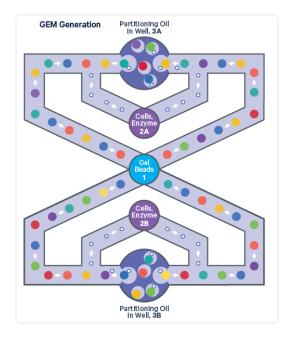

Immediately following GEM generation, the Gel Bead is dissolved, primers are released, and any copartitioned cell is lysed.

Primers containing an Illumina TruSeq Read 1 (read 1 sequencing primer), 16 nt 10x Barcode, 12 nt unique molecular identifier (UMI), and 30 nt poly(dT) sequence are mixed with both the cell lysate and a Master Mix containing reverse transcription (RT) reagents. Incubation of the GEMs produces barcoded, full-length cDNA from poly-adenylated mRNA.

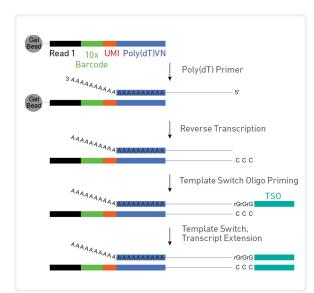


Inside Individual GEMs

Step 2: Post GEM-RT Cleanup & cDNA Amplification


After incubation, GEMs are broken and pooled fractions are recovered. Silane magnetic beads are used to purify the first-strand cDNA from the post GEM-RT reaction mixture, which includes leftover biochemical reagents and primers. Barcoded, full-length cDNA is amplified via PCR to generate sufficient mass for library.

Pooled cDNA Amplification

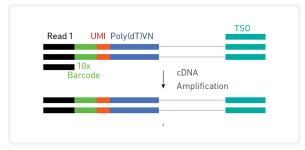

😳 Step 3: 3' Gene Expression Library Construction

Enzymatic fragmentation and size selection are used to optimize the cDNA amplicon size. P5 and P7 adapters, i7 and i5 sample indexes, and TruSeq Read 2 (read 2 primer sequence) are added via End Repair, A-tailing, Adaptor Ligation, and PCR. The final libraries contain the P5 and P7 primers used in Illumina amplification.

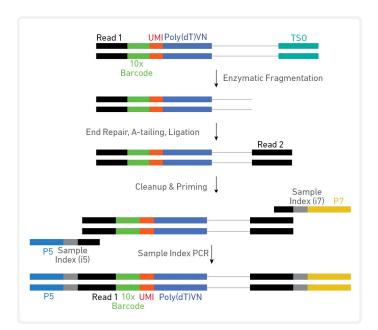


Amplified cDNA Processing (dual index)

Sequencing

The Single Cell 3' library comprises standard Illumina paired-end constructs which begin and end with P5 and P7 adapters. The 16 bp 10x Barcode and 12 bp UMI are encoded in Read 1, while Read 2 is used to sequence the cDNA fragment in 3' Gene Expression libraries. i7 and i5 index sequences are incorporated as the sample index reads. Standard Illumina sequencing primer sites TruSeq Read 1 and TruSeq Read 2 in the 3' Gene Expression libraries are used in paired-end sequencing.

Illumina sequencer compatibility, sample indices, library loading and pooling for sequencing are summarized in the sequencing chapter.

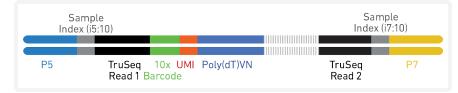

Chromium Single Cell 3' Gene Expression Dual Index Library

See Appendix for Oligonucleotide Sequences on page 86



Tips & Best Practices



DUAL

INDEX

Dual index

specific

protocol step

updates

lcons



Emulsion-safe Plastics

Use validated emulsion-safe plastic consumables when handling GEMs as some plastics can destabilize GEMs.

Cell Concentration

- Recommended starting point is to load ~3,000 cells per reaction, resulting in recovery of ~2,000 cells, and a multiplet rate of ~0.8%. The optimal input cell concentration is 700-1,200 cells/µl.
- The presence of dead cells and debris in the suspension may also reduce the recovery rate. Consult the 10x Genomics Single Cell Protocols Cell Preparation Guide and the Best Practices to Minimize Chromium Next GEM Chip Clogs and Wetting Failure (Documents CG00053 and CG000479 respectively) for more information on preparing cells.

| Multiplet Rate (%) | # of Cells Loaded | # of Cells Recovered |
|--------------------|-------------------|----------------------|
| ~0.8% | ~3,130 | ~2,000 |
| ~1.6% | ~6,320 | ~4,000 |
| ~2.4% | ~9,550 | ~6,000 |
| ~3.2% | ~12,800 | ~8,000 |
| ~4.0% | ~16,100 | ~10,000 |
| ~4.8% | ~19,500 | ~12,000 |
| ~5.6% | ~22,900 | ~14,000 |
| ~6.4% | ~26,300 | ~16,000 |
| ~7.2% | ~29,800 | ~18,000 |
| ~8.0% | ~33,300 | ~20,000 |

General Reagent Handling

- Fully thaw and thoroughly mix reagents before use.
- Keep all enzymes and Master Mixes on ice during setup and use. Promptly move reagents back to the recommended storage.
- Calculate reagent volumes with 10% excess of 1 reaction values.
- Cover Partitioning Oil tubes and reservoirs to minimize evaporation.
- If using multiple chips, use separate reagent reservoirs for each chip during loading.
- Thoroughly mix samples with the beads during bead-based cleanup steps.

50% Glycerol Solution

• Purchase 50% glycerol solution from Ricca Chemical Company, Glycerin (glycerol), 50% (v/v) Aqueous Solution, PN-3290-32.

OR

- Prepare 50% glycerol solution:
 - Mix an equal volume of water and ≥99% Glycerol, Molecular Biology Grade.
 - $\circ~$ Filter through a 0.2 μm filter.
 - Store at −**20**°**C** in 1-ml LoBind tubes. 50% glycerol solution should be equilibrated to room temperature before use.
- Adding glycerol to non-sample chip wells is essential to avoid chip failure.

Pipette Calibration

- Follow manufacturer's calibration and maintenance schedules.
- Pipette accuracy is particularly important when using SPRIselect reagents.

Chromium Next GEM Chip Handling

- Minimize exposure of reagents, chips, and gaskets to sources of particles and fibers, laboratory wipes, frequently opened flip-cap tubes, clothing that sheds fibers, and dusty surfaces.
- After removing the chip from the sealed bag, use in ≤24 h.
- Execute steps without pause or delay, unless indicated. When using multiple chips, load, run, and collect the content from one chip before loading the next.
- Only even number of reactions can be run on the chip. Refer to 1.2 Load Chromium Next GEM Chip M on page 43 for specific instructions.
- Fill all unused paired input wells on a chip with an appropriate volume of 50% glycerol solution before loading the used wells.
- Avoid contacting the bottom surface of the chip with gloved hands and other surfaces. Frictional charging can lead to inadequate priming of the channels, potentially leading to either clogs or wetting failures.
- Minimize the distance that a loaded chip is moved to reach the Chromium X.
- Keep the chip horizontal to prevent wetting the gasket with oil, which depletes the input volume and may adversely affect the quality of the resulting emulsion.

Chromium X Chip Holders

- Chromium X Chip Holders encase Chromium Next GEM Chips used for the HT (high throughput) assay.
- The holder lid flips over to become a stand, holding the chip at 45 degrees for optimal recovery well content removal.
- Squeeze the slider on the back side of the holder together to unlock the lid and return the holder to a flat position.



Chromium Next GEM Chip & Holder Assembly with Gasket

- Close the holder lid. Attach the gasket by holding the tongue (curved end, to the right) and hook the gasket on the left-hand tabs of the holder. Gently pull the gasket toward the right and hook it on the two right-hand tabs.
- DO NOT touch the smooth side of the gasket.
- Open the chip holder.
- Align notch on the chip (upper left corner) and the open holder with the gasket attached.
- Slide the chip to the left until the guide on the holder is inserted into the chip. Depress the right hand side of the chip until the spring-loaded clip engages.
- Keep the assembled unit with the attached gasket until ready for dispensing reagents into the wells.



Chromium Next GEM Chip Loading

- Place the assembled chip and holder flat (gasket attached) on the bench with the lid open.
- Dispense at the bottom of the wells without introducing bubbles.
- When dispensing Gel Beads into the chip, wait for the remainder to drain into the bottom of the pipette tips and dispense again to ensure complete transfer.
- Refer to 1.2 Load Chromium Next GEM Chip M on page 43 for specific instructions.

Gel Bead Handling

- Use one tube of Gel Beads per sample**pair**. DO NOT puncture the foil seals of tubes not used at the time.
- After removing the Gel Bead strip from the packaging, equilibrate the Gel Bead strip to **room temperature** for at least **30 min** before use.
- Store unused Gel Beads at -80°C and avoid more than 12 freeze-thaw cycles. DO NOT store Gel Beads at -20°C.
- Snap the tube strip holder with the Gel Bead strip into a 10x Vortex Adapter. Vortex **30 sec**.



- Centrifuge the Gel Bead strip for **~5 sec** after removing from the holder. Confirm there are no bubbles at the bottom of tubes and the liquid levels look even. Place Gel Bead strip back in the holder and secure the holder lid.
- Ensure that the gel bead strip is positioned with one tube in the left-most position (do not center the strip if using fewer than 8 tubes). Gently depress the lid until light resistance is met. DO NOT attempt to further depress the lid, even if it may be angled with respect to the strip holder.



• If the required volume of beads cannot be recovered, place the pipette tips against the sidewalls and slowly dispense the Gel Beads back into the tubes. DO NOT introduce bubbles into the tubes and verify that the pipette tips contain no leftover Gel Beads. Withdraw the full volume of beads again by pipetting slowly.

10x Gasket Attachment

- **Before reagents are loaded,** attach the gasket by holding the tongue (curved end, to the right) and hook the gasket on the left-hand tabs of the holder. Gently pull the gasket toward the right and hook it on the two right-hand tabs.
- DO NOT touch the smooth side of the gasket.
- Keep the assembled unit with the gasket attached until ready for dispensing reagents into the wells.
- After loading reagents, DO NOT press down on the top of the gasket. Keep the assembly horizontal to avoid wetting the gasket with Partitioning Oil.



10x Magnetic Separator HT

- Offers two positions of the magnets (high and low) relative to a tube, depending on its orientation. Flip the magnetic separator over to switch between high (magnet•**High**) or low (magnet•**Low**) positions.
- The 10x Magnetic Separator HT can accommodate four 8-Tube Strips
- If using MicroAmp 8-Tube Strips, use the high position (magnet•**High**) only throughout the protocol.


10x Magnetic Separator HT



Magnetic Bead Cleanup Steps


- During magnetic bead based cleanup steps that specify waiting "until the solution clears", visually confirm clearing of solution before proceeding to the next step. See panel below for an example.
- The time needed for the solution to clear may vary based on specific step, reagents, volume of reagents etc.
- When processing multiple tube strips, prevent bead over drying after ethanol removal by drying the beads only for the specified time. With the tube strips on the magnet, immediately add indicated buffer (EB or ES1) to all tube strips. Then, remove one tube strip at a time from the magnet and mix.

Visually Confirm Clearing of Magnetic Bead Solution

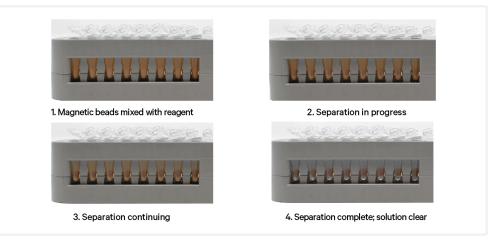

SPRIselect Cleanup & Size Selection

- After aspirating the desired volume of SPRIselect reagent, examine the pipette tips before dispensing to ensure the correct volume is transferred.
- Pipette mix thoroughly as insufficient mixing of sample and SPRIselect reagent will lead to inconsistent results.
- Use fresh preparations of 80% Ethanol.

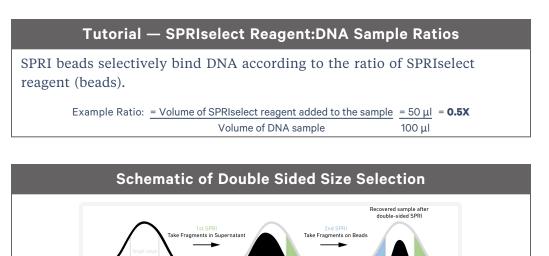

After the first SPRI, supernatant is transferred for a second SPRI while larger fragments are discarded (green). After the second SPRI, fragments on beads are eluted and kept while smaller fragments are discarded (blue). Final sample has a tight fragment size distribution with reduced overall amount (black).

Fragment Size

Tutorial — Double Sided Size Selection1. First SPRIselect: Add 50 µl SPRIselect reagent to 100 µl sample (0.5X).Ratio: = Volume of SPRIselect reagent added to the sample100 µl2. Second SPRIselect: Add 30 µl SPRIselect reagent to supernatant from step a (0.8X).Ratio: = Total Volume of reagent added to the sample (step a + b)= 50 µl + 30 µlOriginal Volume of DNA sample

Enzymatic Fragmentation

Ensure enzymatic fragmentation reactions are prepared on ice and then loaded into a thermal cycler pre-cooled to **4°C** prior to initiating the Fragmentation, End Repair, and A-tailing incubation steps.

Sample Indices in Sample Index PCR

- Choose the appropriate sample index sets to ensure that no sample indices overlap in a multiplexed sequencing run.
- Verify and use the specified index plate only. DO NOT use the plates interchangeably.
- Each well in the Dual Index Plate contains a unique i7 and a unique i5 oligonucleotide.

Index Hopping Mitigation

Index hopping can impact pooled samples sequenced on Illumina sequencing platforms that utilize patterned flow cells and exclusion amplification chemistry. To minimize index hopping, follow the guidelines listed below.


- Remove adapters during cleanup steps.
- Ensure no leftover primers and/or adapters are present when performing post-Library Construction QC.
- Store each library individually at 4°C for up to 72 h or at −20°C for long-term storage. DO NOT pool libraries during storage.
- Pool libraries prior to sequencing. An additional 0.8X SPRI may be performed for the pooled libraries to remove any free adapters before sequencing.
- Hopped indices can be computationally removed from the data generated from single cell dual index libraries.



Step 1:

GEM Generation and Barcoding

| 1.0 Get Started | 37 |
|-----------------------------------|----|
| 1.1 Prepare Master Mix | 39 |
| Assemble Chromium Next GEM Chip M | 40 |
| Sample Loading Guidelines | 41 |
| 1.2 Load Chromium Next GEM Chip M | 43 |
| 1.3 Run Chromium X | 45 |
| GEM Transfer Overview | 46 |
| 1.4 Transfer GEMs | 47 |
| 1.5 GEM-RT Incubation | 49 |

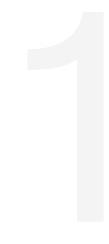

1.0 Get Started

| Overview GEM Generation 1 sample/chip well | Transfer GEM & GEM-RT | cDNA | ine Expression Library Constr. | |
|---|-------------------------------------|---------------------|--|---------|
| Action | ltem | 10x PN | Preparation & Handling | Storage |
| Equilibrate to Ro | oom Temperature | | | |
| | Single Cell 3' HT v3.1 Gel Beads | 2000443 | Equilibrate to room temperature 30 min before loading the chip. | -80°C |
| | RT Reagent B | 2000435/ 2000165 | Vortex, verify no precipitate, centrifuge briefly. | -20°C |
| | Template Switch Oligo | 3000228 | Centrifuge briefly, resuspend in 80 µl Low TE Buffer. Vortex 15 sec at maximum speed, centrifuge briefly, leave at room temperature for ≥ 30 min. After resuspension, store at -80°C. Thaw at room temperature for ≥ 30 min in subsequent uses. | -20°C |
| | Reducing Agent B | 2000087 | Vortex, verify no precipitate, centrifuge briefly. | -20°C |
| Place on Ice | | | | |
| | RT Enzyme C | 2000436/ 2000085 | Centrifuge briefly before adding to the mix. | -20°C |
| | Cell Suspension | _ | | — |
| Obtain | | | | |
| | Partitioning Oil | 220088 | _ | Ambient |
| | Chromium Next GEM Chip M | 2000417 | See Tips & Best Practices. | Ambient |
| | 10x Gasket, HT | 3000614/ 3000656 | See Tips & Best Practices. | Ambient |
| | 10x Vortex Adapter | 330002 | See Tips & Best Practices. | Ambient |

| Action | ltem | 10x PN | Preparation & Handling | Storage |
|--------|---|---------|----------------------------|---------|
| | 50% glycerol solution If using < 16 reactions | — | See Tips & Best Practices. | — |
| | Chromium X Chip Holder | 3000598 | See Tips & Best Practices. | Ambient |

1.1 Prepare Master Mix

a. Prepare Master Mix on ice. Pipette mix 15x and centrifuge briefly.

| Master Mix Add reagen | ts in the order listed | PN | 2X* + 10% (μl) | 8X + 10% (μl) | 16X +10% (μl) |
|---------------------------------|--------------------------|---------------------|----------------------|------------------|------------------|
| ٠ | RT Reagent B | 2000435/ 2000165 | 82.5 | 330.0 | 660.0 |
| | Template Switch Oligo | 3000228 | 10.3 | 41.4 | 82.7 |
| \bigcirc | Reducing Agent B | 2000087 | 8.5 | 34.3 | 68.6 |
| ٠ | RT Enzyme C | 2000085/ 2000436 | 38.5 | 154.0 | 308.0 |
| | Total | - | 139.8 | 559.7 | 1119.3 |
| *Volume fo | r 2 rxns; only even ni | umber of rxns | can be run | on the chip | |

a. Add 63.6 µl Master Mix into each tube of a PCR 8-tube strip on ice.

Assemble Chromium Next GEM Chip M




See Tips & Best Practices on page 23 for chip handling instructions.

- a. Close the holder lid. Attach the gasket by holding the tongue (curved end, to the right) and hook the gasket on the left-hand tabs of the holder. Gently pull the gasket toward the right and hook it on the two right-hand tabs.
- **b.** DO NOT touch the smooth side of the gasket.
- c. Open the chip holder.



- **d.** Remove the chip from the sealed bag. Use the chip within \leq 24 h.
- **e.** Align notch on the chip (upper left corner) and the open holder with the gasket attached.
- **f.** Slide the chip to the left until the guide on the holder is inserted into the chip. Depress the right hand side of the chip until the spring-loaded clip engages.
- **g.** Keep the assembled unit with the attached gasket open until ready for and while dispensing reagents into the wells.
- **h.** DO NOT touch the smooth side of the gasket.
- **i.** After loading reagents, close the chip holder. DO NOT press down on the top of the gasket.

For GEM generation, load the indicated reagents only in the specified rows, starting from row labeled 1, followed by rows labeled 2A & 2B and 3A & 3B.

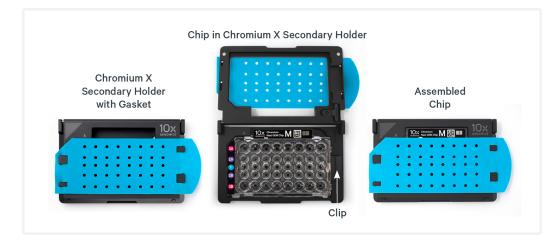

Sample Loading Guidelines

Read these guidelines before loading Chromium Next GEM Chip M

- Up to 16 independent samples can be run on the chip.
- Only even number of reactions can be run on the chip.
- Corresponding wells in rows 2A & 2B of the chip should both be either loaded with independent samples/mock-sample or with glycerol.
- For even number of samples:

Load independent samples (Master Mix + Cells) in pairs in rows 2A & 2B. See Example 1 below.

• For odd number of samples :

Load the unpaired sample in a well in row 2A and a mock-sample (Master Mix + Water) in the corresponding well in row 2B. Additionally, add Partitioning Oil to the corresponding well in row 3B. See Example 2 below.

• Follow the step-by-step chip loading instructions provided in step 1.2.



Sample loading configuration examples

Cell Suspension Volume Calculator

Volume of Cell Suspension Stock per reaction (µl) | Volume of Nuclease-free Water per reaction (µl)

DO NOT add nuclease-free water directly to single cell suspension. Add nuclease-free water to the Master Mix. Refer to step 1.2c.

| Cell Stock | Targeted Cell Recovery | | | | | | | | | | | |
|---------------------|------------------------------------|--------------|--------------|---------------|--------------------------------|---------------------|--------------|--------------|---|--------------|--|--|
| Conc. (Cells/µl) | 2000 | 4000 | 6000 | 8000 | 10000 | 12000 | 14000 | 16000 | 18000 | 20000 | | |
| 100 | 31.3 55.1 | 63.2 23.2 | n/a | n/a | n/a | n/a | n/a | n/a | n/a | n/a | | |
| 200 | 15.6 70.8 | 31.6 54.8 | 47.8 38.6 | 64.1 22.3 | 80.7 5.7 | n/a | n/a | n/a | n/a | n/a | | |
| 300 | 10.4 | 21.1 | 31.8 | 42.7 | 53.8 | 64.9 | 76.2 | n/a | n/a | n/a | | |
| | 76.0 | 65.3 15.8 | 54.6 23.9 | 43.7 32.1 | 32.6 40.3 | 21.5 48.7 | 10.2 57.2 | 65.8 | 74.4 | 83.2 | | |
| 400 | 78.6 | 70.6 | 62.5 | 54.3 | 46.1 | 37.7 | 29.2 | 20.6 | 12.0 | 3.2 | | |
| 500 | 6.3 80.1 | 12.6 73.8 | 19.1 67.3 | 25.6 60.8 | 32.3 54.1 | 39.0 47.4 | 45.7 40.7 | 52.6 33.8 | 59.6 26.8 | 66.6 19.8 | | |
| | 5.2 | 10.5 | 15.9 | 21.4 | 26.9 | 32.5 | 38.1 | 43.8 | 49.6 | 55.5 | | |
| 600 | 81.2 | 75.9 | 70.5 | 65.0 | 59.5 | 53.9 | 48.3 | 42.6 | 36.8 | 30.9 | | |
| 700 | 4.5 81.9 | 9.0 77.4 | 13.6 72.8 | 18.3 68.1 | 23.0 63.4 | 27.8 58.6 | 32.7 53.7 | 37.6 48.8 | 42.5 43.9 | 47.6 38.8 | | |
| | 3.9 | 77.4 | 11.9 | 16.0 | 20.2 | 24.4 | 28.6 | 32.9 | 37.2 | 41.6 | | |
| 800 | 82.5 | 78.5 | 74.5 | 70.4 | 66.2 | 62.0 | 57.8 | 53.5 | 49.2 | 44.8 | | |
| 900 | 3.5 82.9 | 7.0 79.4 | 10.6 75.8 | 14.2 72.2 | 17.9 68.5 | 21.6 64.8 | 25.4 61.0 | 29.2 57.2 | 33.1 53.3 | 37.0 49.4 | | |
| 1000 | 3.1 | 6.3 | 9.6 | 12.8 | 16.1 | 19.5 | 22.9 | 26.3 | 29.8 | 33.3 | | |
| 1000 | 83.3 | 80.1 | 76.8 | 73.6 | 70.3 | 66.9 | 63.5 | 60.1 | 56.6 | 53.1 | | |
| 1100 | 2.8 83.6 | 5.7 80.7 | 8.7 77.7 | 11.7 74.7 | 14.7 71.7 | 17.7 68.7 | 20.8 65.6 | 23.9 62.5 | 27.1 59.3 | 30.3 56.1 | | |
| 1200 | 2.6 | 5.3 | 8.0 | 10.7 | 13.4 | 16.2 | 19.1 | 21.9 | 24.8 | 27.7 | | |
| 1200 | 83.8 | 81.1 | 78.4 | 75.7 | 73.0 | 70.2 | 67.3 | 64.5 | 61.6 | 58.7 | | |
| 1300 | 2.4 84.0 | 4.9 81.5 | 7.3 79.1 | 9.9 76.5 | 12.4 74.0 | 15.0 71.4 | 17.6 68.8 | 20.2 66.2 | 22.9 63.5 | 25.6 60.8 | | |
| | 2.2 | 4.5 | 6.8 | 9.2 | 11.5 | 13.9 | 16.3 | 18.8 | 21.3 | 23.8 | | |
| 1400 | 84.2 | 81.9 | 79.6 | 77.2 | 74.9 | 72.5 | 70.1 | 67.6 | 65.1 | 62.6 | | |
| 1500 | 2.1 | 4.2 | 6.4 | 8.5 | 10.8 | 13.0 | 15.2 | 17.5 | 19.9 | 22.2 | | |
| | 84.3 2.0 | 82.2 4.0 | 80.0 6.0 | 77.9 8.0 | 75.6 | 73.4 12.2 | 71.2 | 68.9 16.4 | 66.5 18.6 | 64.2 20.8 | | |
| 1600 | 84.4 | 82.4 | 80.4 | 78.4 | 76.3 | 74.2 | 72.1 | 70.0 | 67.8 | 65.6 | | |
| 1800 | 1.8 | 3.7 | 5.6 | 7.5 | 9.5 | 11.5 | 13.5 | 15.5 | 17.5 | 19.6 | | |
| 1700 | 84.6 | 82.7 | 80.8 | 78.9 | 76.9 | 74.9 | 72.9 | 70.9 | 68.9 | 66.8 | | |
| 1800 | 1.7 | 3.5 | 5.3 | 7.1 | 9.0 | 10.8 | 12.7 | 14.6 | 16.5 | 18.5 | | |
| 1000 | 84.7 | 82.9 | 81.1 | 79.3 | 77.4 | 75.6 | 73.7 | 71.8 | 69.9 | 67.9 | | |
| 1900 | 1.6 | 3.3 | 5.0 | 6.7 | 8.5 | 10.3 | 12.0 | 13.8 | 15.7 | 17.5 | | |
| | 84.8 | 83.1 3.2 | 81.4 4.8 | 79.7 6.4 | 77.9 8.1 | 76.1 9.7 | 74.4 | 72.6 13.2 | 70.7 | 68.9 16.6 | | |
| 2000 | 84.8 | 83.2 | 81.6 | 80.0 | 78.3 | 76.7 | 75.0 | 73.2 | 71.5 | 69.8 | | |
| | G rey boxes: allowable v | olume | Optimal ce | l stock conc. | Blue boxes: for cell recove | ery target of 2 | ,000-20,000 | Low transf | Yellow boxes er volume tha er cell load va | t may resu | | |

1.2 Load Chromium Next GEM Chip M

- After removing chip from the sealed bag, use in ≤24 h.
 - Open the lid (gasket attached) of the assembled chip and lay flat for loading.
 - When loading the chip, raising and depressing the pipette plunger should each take **~5 sec**. When dispensing, raise the pipette tips at the same rate as the liquid is rising, keeping the tips slightly submerged.

a. Add 50% glycerol solution to each unused well

(if loading less than 16 samples/chip)

- 130 µl in each unused well in row labeled 1
- 140 µl in each unused well in rows labeled 2A & 2B
- 140 μl in each unused well in rows labeled 3A & 3B

DO NOT use any substitute for 50% glycerol solution. For odd number of samples, if a sample is loaded in a well in row 2A, load a mock-sample (Master Mix + 86.4 μ l water) and NOT glycerol in the corresponding well in row 2B.

b. Prepare Gel Beads

- Snap the tube strip holder with the Gel Bead strip into a 10x Vortex Adapter. Vortex **30 sec**.
- Remove the Gel Bead strip from the holder and centrifuge it for
 ~5 sec. Confirm there are no bubbles at the bottom of the tubes & the liquid levels are even.
- Place the Gel Bead strip back in the holder. Secure the holder lid.

c. Prepare Master Mix + Cell suspension

- Refer to the Cell Suspension Volume Calculator Table.
- Add the appropriate volume of nuclease-free water to Master Mix. Pipette mix 5X.
- Gently pipette mix the cell suspension and add corresponding volume of single cell suspension to Master Mix. Total of 150 μl in each tube.



 \forall







d. Load Row Labeled 1

- Puncture the foil seal of the Gel Bead tubes. Slowly aspirate 130 µl Gel Beads.
- Dispense into the wells in **row labeled 1** without introducing bubbles.
- Wait 30 sec.

e. Load Rows Labeled 2A, 2B

ONLY even number of reactions should be run on the chip. See Sample Loading Guidelines for more information and examples.

- Up to 16 independent samples can be run on the chip. Sample inputs for 2A should be equal to 2B (e.g. if processing only 8 samples, run 4 in 2A and 4 in 2B)
- First, process up to 8 samples: Gently pipette mix the Master Mix + Cell Suspension (prepared at step 1.2c) using a multichannel pipette. Using the same pipette tips, dispense 140 µl Master Mix + Cell Suspension into the bottom center of wells in row labeled 2A.
- Next, process up to 8 additional samples: Gently pipette mix the Master Mix + Cell Suspension (prepared at step 1.2c) using a multichannel pipette. Using the same pipette tips, dispense 140 µl Master Mix + Cell Suspension into the bottom center of wells in row labeled 2B.
- Wait 30 sec.

f. Load Rows Labeled 3A, 3B

• Dispense 140 μl Partitioning Oil into the wells in rows labeled **3A & 3B** from a reagent reservoir.

Failure to add Partitioning Oil to the rows labeled 3A and 3B will prevent GEM generation and can damage Chromium X.

g. Prepare for Run

• Close the lid (gasket already attached). DO NOT touch the smooth side of the gasket. DO NOT press down on the top of the gasket.

Run the chip in Chromium X **immediately** after loading the Partitioning Oil. ONLY even number of reactions should be run on the chip. See Sample Loading Guidelines on page 41 for more information & examples.



130 µl





1.3 Run Chromium X




Consult the Chromium X Series (X/iX) User Guide (CG000396) for detailed instrument operation instructions and follow the Chromium X touchscreen prompts for execution.

- a. Press the eject button on the Chromium X to eject the tray.If the eject button is not touched within 1 min, tray will close automatically. System requires a few seconds before the tray can be ejected again.
- **b.** Place the assembled chip with the gasket in the tray, ensuring that the chip stays horizontal. Press the button to retract the tray.
- **c.** Press the play button.




d. At completion of the run (~18 min), Chromium X will chime. **Immediately** proceed to the next step.

Chromium X

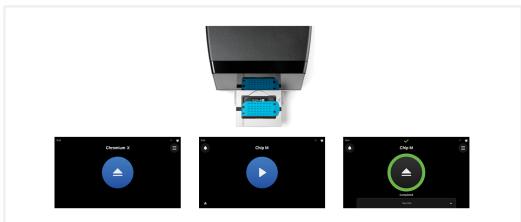


GEM Transfer Overview

For a sample loaded in a well in either row 2A or 2B of the chip, GEMs are retrieved from the corresponding well in row 3A and 3B and transferred to two tubes. The example below shows transfer of GEMs generated from 16 samples. GEMs from the chip are transferred to four tube strips, where the GEMs generated from each sample are transferred to 2 corresponding tubes in the indicated tube strips.

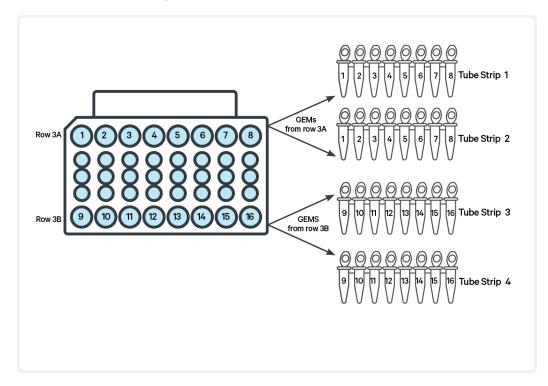

1.4 Transfer GEMs

- **a.** Label four tube strips and place on ice.
- **b.** Press the eject button of Chromium X and remove the chip.
- **c.** Discard the gasket. Open the chip holder. Fold the lid back until it clicks to expose the wells at 45 degrees.
- **d.** Visually check the volume in rows labeled 1, 2A, and 2B. Abnormally high volume relative to other wells indicates a clog. Significant volume of non-sample fluid is expected in rows 2A and 2B after a successful run and does not indicate a sample clog.
- **e.** Retrieve GEMs from row labeled 3A: Slowly aspirate **90 μl** GEMs from the lowest points of the recovery wells in the top row labeled 3A without creating a seal between the tips and the bottom of the wells.

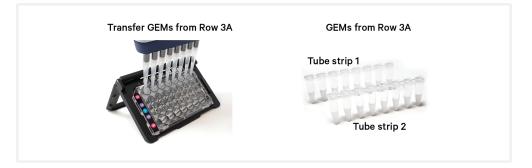

f. Withdraw pipette tips from the wells. GEMs should appear opaque and uniform across all channels. Excess Partitioning Oil (clear) in the pipette tips indicates a potential clog.



- **g.** Over the course of **~20 sec**, dispense GEMs into first tube strip on ice with the pipette tips against the sidewalls of the tubes.
- **h.** Using the same pipette tips, slowly aspirate remaining **90 μl** GEMs from the wells in the top row labeled 3A and dispense in second tube strip as

described above.

i. Retrieve from row labeled 3B: Slowly aspirate **90 μl** GEMs from the lowest points of the recovery wells in the bottom row labeled 3B without creating a seal between the tips and the bottom of the wells.



- **j.** Repeat steps f and g, dispensing GEMs into third tube strip on ice with the pipette tips against the sidewalls of the tubes.
- **k.** Using the same pipette tips, slowly aspirate remaining **90 μl** GEMs from the wells in the bottom row labeled 3B and dispense in fourth tube strip as described above.
- **1.** If multiple chips are run back-to-back, cap/cover the GEM-containing tube strips and place on ice for no more than **1 h**.

1.5 GEM-RT Incubation

Use a thermal cycler that can accommodate at least 100 μ l volume. A volume of 125 μ l is the preferred setting on Bio-Rad C1000 Touch. In alternate thermal cyclers, use highest reaction volume setting.

a. Incubate in a thermal cycler with the following protocol.

| Lid Temperature | Reaction Volume | Run Time |
|-----------------|------------------------|----------------------|
| 53°C | 125 µl | ~55 min |
| Step | Temperature | Time <i>hh:mm:ss</i> |
| 1 | 53°C | 00:45:00 |
| 2 | 85°C | 00:05:00 |
| 3 | 4°C | Hold |

b. Store at **4°C** for up to **72 h** or at −**20°C** for up to **a week**, or proceed to the next step.



Step 2:

Post GEM-RT Cleanup & cDNA Amplification

| 2.0 Get Started | 51 |
|---|----|
| 2.1 Post GEM-RT Cleanup - Dynabeads | 53 |
| 2.2 cDNA Amplification | 55 |
| 2.3 cDNA Cleanup - SPRIselect | 57 |
| 2.4 Post cDNA Amplification QC & Quantification | 58 |

2.0 Get Started

| Overview | | | | |
|------------------|--|--|--|---------|
| | Transfer GEM & GEM-RT | GEM-RT Cleanup cDNA Amp. & Cleanup | Combine cDNA (post Cleanup) | |
| Action | ltem | 10x PN | Preparation & Handling | Storage |
| Equilibrate to R | Room Temperature | | | |
| | Reducing Agent B | 2000087 | Thaw, vortex, verify no precipitate, centrifuge. | -20°C |
| | cDNA Primers | 2000089/ 2000439 | Thaw, vortex, centrifuge briefly. | -20°C |
| | Dynabeads MyOne SILANE | 2000048 | Vortex thoroughly (≥30 sec) immediately before adding to the mix. | 4°C |
| | Beckman Coulter SPRIselect Reagent | _ | Manufacturer's recommendations. | _ |
| | Agilent Bioanalyzer High Sensitivity Kit If used for QC & quantification | _ | Manufacturer's recommendations. | _ |
| | Agilent TapeStation ScreenTape & Reagents If used for QC & quantification | _ | Manufacturer's recommendations. | _ |
| | Qubit dsDNA HS Assay Kit If used for QC & quantification | _ | Manufacturer's recommendations. | _ |
| | DNA High Sensitivity Reagent Kit If used for QC | _ | Manufacturer's recommendations. | _ |
| Place on Ice | | | | |
| | Amp Mix Retrieve from | 2000440/ 2000047 | Vortex. Ensure all liquid is at the bottom of the tube. | -20°C |

User Guide | Chromium Next GEM Single Cell 3' HT Reagent Kits v3.1 (Dual Index)

| Action | ltem | 10x PN | Preparation & Handling | Storage |
|--------------|--|---------------------|---|---------|
| | Single Cell 3' HT GEM Kit | | | |
| Thaw at 65°C | | | | |
| | Cleanup Buffer | 2000088/ 2000438 | Thaw for 10 min at 65° C either in a water or bead bath, mixing every 5 min (for 16 rxn kit) or at max speed on a thermomixer (for 8 rxn kit). Verify no visible crystals. Cool to room temperature. | -20°C |
| Obtain | | | | |
| | Recovery Agent | 2000434 | _ | Ambient |
| | Qiagen Buffer EB | — | Manufacturer's recommendations. | — |
| | Bio-Rad 10% Tween 20 | — | Manufacturer's recommendations. | _ |
| | 10x Magnetic Separator HT | 2000431 | _ | Ambient |
| | Prepare 80% Ethanol Prepare 60 ml for 16 reactions. | _ | _ | _ |

2.1 Post GEM-RT Cleanup – Dynabeads



a. Add **125 μl** Recovery Agent to each sample at room temperature. DO NOT pipette mix or vortex the biphasic mixture. Wait **2 min**.

The resulting biphasic mixture contains Recovery Agent/Partitioning Oil (pink) and aqueous phase (clear), with no persisting emulsion (opaque).

If biphasic separation is incomplete:

- Firmly secure the cap on the tube strip, ensuring that no liquid is trapped between the cap and the tube rim.
- Mix by inverting the capped tube strip 5x, centrifuge briefly, and proceed to step b. DO NOT invert without firmly securing the caps.

A smaller aqueous phase volume indicates a clog during GEM generation.



- **b.** Slowly remove and discard **125 μl** Recovery Agent/Partitioning Oil (pink) from the bottom of the tube. DO NOT aspirate any aqueous sample.
- c. Prepare Dynabeads Cleanup Mix.

Before using Dynabeads MyOne SILANE to prepare the Dynabeads Cleanup Mix:



• Vortex the Dynabeads thoroughly (≥**30 sec**) **immediately** before adding to the mix.

• Aspirate full liquid volume in the Dynabead tube with a pipette tip to verify that beads have not settled in the bottom of the tube. If clumps are present, pipette mix to resuspend completely. DO NOT centrifuge before adding to the mix.

| Dynabeads Cleanup Mix Add reagents in the order listed | | PN | 2X + 10% (μl) | 16X + 10% (μl) | 32X + 10% (μl) | |
|--|------------------------|---------------------|---------------------|----------------------|----------------------|--|
| | Cleanup Buffer | 2000088/ 2000438 | 400.4 | 3203.2 | 6406.4 | |
| | Dynabeads MyOne SILANE | 2000048 | 17.6 | 140.8 | 281.6 | |
| \bigcirc | Reducing Agent B | 2000087 | 11 | 88 | 176 | |
| | Nuclease-free Water | _ | 11 | 88 | 176 | |
| | Total | | 440 | 3520 | 7040 | |

d. Vortex and add **200 \mul** to each tube. Pipette mix 10x (pipette set to 200 μ l).



- e. Incubate 10 min at room temperature. Pipette mix again at ~5 min after start of incubation to resuspend settled beads.
- **f.** Prepare Elution Solution I. Vortex and centrifuge briefly.

| Elution Soluti Add reagents i | on I in the order listed | PN | 1X (μl) | 40X (μl) |
|----------------------------------|-----------------------------|---------|---------|----------|
| | Buffer EB | — | 98 | 3920 |
| | 10% Tween 20 | _ | 1 | 40 |
| 0 | Reducing Agent B | 2000087 | 1 | 40 |
| | Total | | 100 | 4000 |

TIPS

g. At the end of 10 min incubation, place the tube strips in slots 1-4 of a 10x Magnetic Separator HT• High position (magnet•High) until the solution clears.

A white interface between the aqueous phase and Recovery Agent is normal.

- **h.** Remove the supernatant.
- i. Add **300 µl** 80% ethanol to the pellet while on the magnet. Wait **30 sec**.

- j. Remove the ethanol.
- k. Add 200 µl 80% ethanol to pellet. Wait 30 sec.
- **1.** Remove the ethanol.
- m. Centrifuge briefly. Place on the magnet- Low.
- n. Remove remaining ethanol. Air dry for 1 min.
- **ο.** To avoid over drying the pellet, **immediately** add **36 μl** Elution Solution I to the tube strips, cap the tubes, remove all tube strips from the magnet, and centrifuge briefly.
- **p.** Pipette mix (pipette set to 30 µl) without introducing bubbles.
- q. Incubate 2 min at room temperature.
- **r.** Place on the magnet• **Low** until the solution clears.
- s. Transfer 35 µl sample to a new tube strip.

2.2 cDNA Amplification

| a. | Prepare | cDNA | Amplification | Mix of | 1 ice. | Vortex | and | centrifuge | briefly. |
|----|---------|------|---------------|--------|--------|--------|-----|------------|----------|
|----|---------|------|---------------|--------|--------|--------|-----|------------|----------|

| Amplification Reaction Mix Add reagents in the order listed | | PN | 2X*+ 10% (μl) | 16X + 10% (μl) | 32X 10% (μl) |
|--|---|---------------------|---------------------|----------------------|--------------------|
| 0 | Amp Mix Retrieve from Single Cell 3' HT GEM kit | 2000047/ 2000440 | 110 | 880 | 176 |
| | cDNA Primers | 2000439/ 2000089 | 33 | 264 | 528 |
| | Total | | 143 | 1144 | 228 |

well.

- **b.** Add **65 µl** cDNA Amplification Reaction Mix to **35 µl** sample.
- **c.** Pipette mix 15x (pipette set to 90 µl). Centrifuge briefly.
- **d.** Incubate in a thermal cycler with the following protocol.

| Lid Temperature | Reaction Volume | Run Time |
|-----------------|----------------------------|---------------------------|
| 105°C | 100 µl | ~30-45 min |
| Step | Temperature | Time <i>hh:mm:ss</i> |
| 1 | 98°C | 00:03:00 |
| 2 | 98°C | 00:00:15 |
| 3 | 63°C | 00:00:20 |
| 4 | 72°C | 00:01:00 |
| 5 | Go to Step 2, see table be | low for total # of cycles |
| 6 | 72°C | 00:01:00 |
| 7 | 4°C | Hold |



Recommended starting point for cycle number optimization.The optimal cycle number is a trade-off between generating sufficient final mass for libraries & minimizing PCR amplification artifacts. Select PCR cycles based on the target recovery per GEM well and not on the estimated number of cells in the two tubes containing the split sample. The cDNA cycles should be reduced if sampling large numbers of cells.

| Targeted Cell Recovery (per GEM well) | Total Cycles |
|---------------------------------------|--------------|
| <12,000 | 12 |
| >12,000 | 11 |


STOP

e. Store at 4°C for up to 72 h or −20°C for ≤1 week, or proceed to the next step.

2.3 cDNA Cleanup – SPRIselect

- a. Vortex to resuspend the SPRIselect reagent. Add 60 µl SPRIselect reagent (0.6X) to each sample and pipette mix 15x (pipette set to 150 µl).
- **b.** Incubate **5 min** at **room temperature**.
- c. Place on the magnet-High until the solution clears.
- d. Remove the supernatant.
- e. Add 200 µl 80% ethanol to the pellet. Wait 30 sec.
- f. Remove the ethanol.
- g. Repeat steps e and f for a total of 2 washes.
- **h.** Centrifuge briefly and place on the magnet-Low.
- **i.** Remove any remaining ethanol. Air dry for **2 min**. DO NOT exceed **2 min** as this will decrease elution efficiency.
- j. Remove from the magnet. Add 41 µl Buffer EB. Pipette mix 15x.
- k. Incubate 2 min at room temperature.
- 1. Place the tube strip on the magnet-High until the solution clears.
- **m.** Transfer **40 µl** sample to a new tube strip.

For each sample, the amplified cDNA will be in two corresponding tubes. Transfer **15** μ I cDNA from each of the two sample tubes to a third tube and pipette mix. Use for downstream steps. Store remaining cDNA volume in tubes at **4°C** for up to **72** h or **-20°C** for up to **4 weeks** for generating additional libraries.

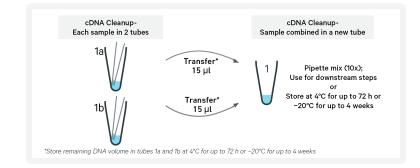


STOP

n. Store at 4°C for up to 72 h or −20°C for ≤4 weeks, or proceed to the next step.

2.4 Post cDNA Amplification QC & Quantification

- **a.** Run **1 μl** sample at **1:10 dilution** on an Agilent Bioanalyzer High Sensitivity chip.
- Samples of RNA-rich cells may require additional dilution in nuclease-free water. The number of distinct peaks may vary. Higher molecular weight product (2,000-9,000 bp) may be present. This does not affect sequencing.

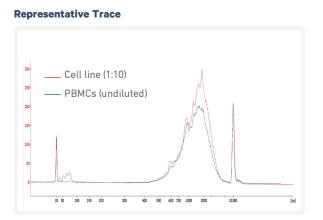

See example calculation in the following page.

Alternate Quantification Methods

Agilent TapeStation

LabChip


Qubit Fluoromter and Qubit dsDNA HS Assay Kit

See Appendix for

- Agilent TapeStation Traces on page 84
- LabChip Traces on page 85

Example Calculation

- i. Select Region: Under the "Electropherogram" view, choose the "Region Table." Manually select the region of ~200 ~9000 bp.
- **ii.** Note Concentration [pg/µl]

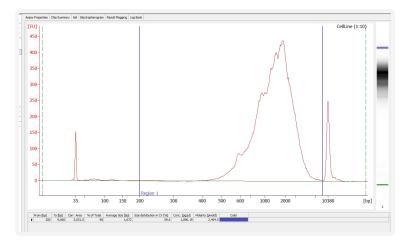

iii. Calculate: Multiply the cDNA concentration [pg/µl] reported via Agilent
 2100 Expert Software by elution volume and dilution factor and divide by
 1000 to obtain the total cDNA yield in ng. Carry forward ONLY 25% of
 total cDNA yield into 3' Gene Expression Library Construction.

Example Calculation of cDNA Total Yield

Concentration: 1890.19 pg/µl Elution volume: 80 µl; Dilution Factor: 10

Total cDNA Yield

= <u>Conc'n (pg/μl) x Elution Vol. (μl) x Dilution Factor</u> 1000 (pg/ng)

> = <u>1890.19 x 80 x 10</u> 1000 (pg/ng)

=1512.15 ng

Carrying Forward ONLY 25% of total cDNA yield for 3' GEX Library

=Total cDNA x 0.25 =1512.15 ng x 0.25

=378.03 ng

Refer to step 3.5 for appropriate number of Sample Index PCR cycles based on carry forward cDNA yield/input cDNA.



Step 3:

3' Gene Expression Library Construction

| 3.0 Get Started | 61 |
|--|----|
| Step Overview (Step 3.1d) | 63 |
| 3.1 Fragmentation, End Repair & A-tailing | 64 |
| 3.2 Post Fragmentation, End Repair & A-tailing Double Sided - SPRIselect | 65 |
| 3.3 Adaptor Ligation | 66 |
| 3.4 Post Ligation Cleanup - SPRIselect | 67 |
| 3.5 Sample Index PCR | 68 |
| 3.6 Post Sample Index PCR Double Sided Size Selection - SPRIselect | 69 |
| 3.7 Post Library Construction QC | 70 |



3.0 Get Started

| Overview | | |
|----------|---|--------------------|
| | Transfer GEM & GEM-RT Cleanup Combine cDNA Amp. & Cleanup (post Cleanup) | ssion y Constr. |

| Actio | n | ltem | 10x PN | Preparation & | Storage |
|-------|--------|--|---------|---|---------|
| | | Room Temperature | | Handling | |
| | ٠ | Fragmentation Buffer | 2000091 | Thaw, vortex, verify no precipitate, centrifuge. | -20°C |
| | • | Adaptor Oligos | 2000094 | Thaw, vortex, centrifuge briefly. | -20°C |
| | • | Ligation Buffer | 2000092 | Thaw, vortex thoroughly (≥30 sec) immediately before adding to the mix. | -20°C |
| | A | Dual Index Plate TT Set A Verify name & PN. Use indicated plate only | 3000431 | _ | -20°C |
| | | Beckman Coulter SPRIselect Reagent | _ | Manufacturer's recommendations. | _ |
| | | Agilent Bioanalyzer High Sensitivity Kit If used for QC | — | Manufacturer's recommendations. | _ |
| | | Agilent TapeStation ScreenTape & Reagents If used for QC | _ | Manufacturer's recommendations. | _ |
| | | DNA High Sensitivity Reagent Kit If used for QC | _ | Manufacturer's recommendations. | _ |
| Place | on Ice | | | | |
| | • | Fragmentation Enzyme Ensure that Fragmentation Buffer and Fragmentation Enzyme from the same kit are used together. Lots are matched for optimal performance | 2000090 | Centrifuge briefly. | -20°C |
| | | DNA Ligase | 220110 | Centrifuge briefly. | -20°C |
| | 0 | Amp Mix Retrieve from Library | 2000047 | Centrifuge briefly. | -20°C |


User Guide | Chromium Next GEM Single Cell 3' HT Reagent Kits v3.1 (Dual Index)

| Action | Item | 10x PN | Preparation & Handling | Storage |
|--------|--|---------|---------------------------------|---------|
| | Construction Kit. | | | |
| | KAPA Library Quantification Kit for Illumina Platforms | _ | Manufacturer's recommendations. | _ |
| Obtain | | | | |
| | Qiagen Buffer EB | — | Manufacturer's recommendations. | Ambient |
| | 10x Magnetic Separator HT | 2000431 | See Tips & Best Practices. | Ambient |
| | Prepare 80% Ethanol Prepare 20 ml for 8 reactions. | _ | Prepare fresh. | Ambient |

Step Overview (Step 3.1d)

Correlation between input & library complexity

A Single Cell 3' Gene Expression library is generated using a fixed proportion (20 µl, 25%) of the total cDNA (80 µl) obtained at step 2.3m. The complexity of this library will be comparable to one generated using a higher proportion (>25%) of the cDNA. The remaining proportion (60 µl, 75%) of the cDNA may be stored at **4°C** for up to **72 h** or at **-20°C** for longer-term storage (up to **4 weeks**).

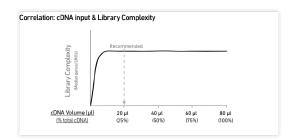

Note that irrespective of the total cDNA yield (ng), which may vary based on cell type, targeted cell recovery etc., this protocol has been optimized for a broad range of input mass (ng), as shown in the example below. The total number of SI PCR cycles (step 3.5d) should be optimized based on carrying forward a fixed proportion (20μ l, 25%) of the total cDNA yield calculated during Post cDNA Amplification QC & Quantification (step 2.4).

| | Torgo | Total cDNA Input into cDNA Fragmentation | | SI PCR Cycle | | |
|---------------------|-------|---|---------------|----------------|-----------|--------|
| Cell | | overy | Yield (ng) | Volume (µl) | Mass (ng) | Number |
| High RNA Content | Low | • | 500 ng | 20 µl | 125 ng | 12 |
| | High | | 3800 ng | 20 µl | 950 ng | 9 |
| Low RNA Content | Low | | 2 ng | 20 µl | 0.5 ng | 16 |
| 203 | High | | 400 ng | 20 µl | 100 ng | 13 |

Example: Library Construction Input Mass & SI PCR Cycles

3.1 Fragmentation, End Repair & A-tailing

a. Prepare a thermal cycler with the following incubation protocol.

| Lid Temperature | Reaction Volume | Run Time |
|---|------------------------|-------------------------|
| 65°C | 50 µl | ~35 min |
| Step | Temperature | Time hh:mm:ss |
| Pre-cool block Pre-cool block prior to preparing Fragmentation Mix | 4°C | Hold |
| Fragmentation | 32°C | 00:05:00 |
| End Repair & A-Tailing | 65°C | 00:30:00 |
| Hold | 4°C | Hold |

- **b.** Vortex Fragmentation Buffer. Verify there is no precipitate.
- c. Prepare Fragmentation Mix on ice. Pipette mix and centrifuge briefly.

| Fragmentation Mix Add reagents in the order listed | PN | 1Χ (μl) | 8X + 10% (μl) | 16X + 10% (μl) |
|--|---------|------------|------------------|-------------------|
| Fragmentation Buffer | 2000091 | 5 | 44 | 88 |
| Fragmentation Enzyme | 2000090 | 10 | 88 | 176 |
| Total | | 15 | 132 | 264 |

- **d.** Transfer **ONLY 20 μl** purified cDNA sample from Cleanup (step 2.3n) to a tube strip.
 - Note that only 20 μl (25%) cDNA sample is sufficient for generating 3' Gene Expression library.
 - The remaining cDNA sample can be stored at 4°C for up to 72 h or at -20°C for up to 4 weeks for generating additional 3' Gene Expression libraries.
- e. Add 15 µl Buffer EB to each sample.
- **f.** Add **15** μ l Fragmentation Mix to each sample.
- g. Pipette mix 15x (pipette set to 35 µl) on ice. Centrifuge briefly.
- **h.** Transfer into the pre-cooled thermal cycler (**4**°**C**) and press "SKIP" to initiate the protocol.

3.2 Post Fragmentation, End Repair & A-tailing Double Sided – SPRIselect

- **a.** Vortex to resuspend SPRIselect reagent. Add **30 μl** SPRIselect **(0.6X)** reagent to each sample. Pipette mix 15x (pipette set to 75 μl).
- b. Incubate 5 min at room temperature.
- **c.** Place on the magnet**·High** until the solution clears. DO NOT discard supernatant.
- **d.** Transfer **75** µl supernatant to a new tube strip.
- e. Vortex to resuspend SPRIselect reagent. Add 10 μl SPRIselect reagent (0.8X) to each sample. Pipette mix 15x (pipette set to 80 μl).
- f. Incubate 5 min at room temperature.
- g. Place on the magnet-High until the solution clears.
- h. Remove 80 µl supernatant. DO NOT discard any beads.
- i. Add 125 µl 80% ethanol to the pellet. Wait 30 sec.
- **j.** Remove the ethanol.
- **k.** Repeat steps i and j for a total of 2 washes.
- **I.** Centrifuge briefly. Place on the magnet**·Low** until the solution clears. Remove remaining ethanol pipetting slowly. DO NOT over dry to ensure maximum elution efficiency.
- m. Remove from the magnet. Add 51 μl Buffer EB to each sample. Pipette mix 15x.
- **n.** Incubate **2 min** at **room temperature**.
- **o.** Place on the magnet**·High** until the solution clears.
- **p.** Transfer **50 µl** sample to a new tube strip pipetting slowly.

3.3 Adaptor Ligation

a. Prepare Adaptor Ligation Mix. Pipette mix and centrifuge briefly.

| Adaptor Ligation Mix Add reagents in the order listed | PN | 1Χ (μl) | 8X + 10% (μl) | 16X + 10% (μl) |
|--|---------|------------|------------------|-------------------|
| Ligation Buffer | 2000092 | 20 | 176 | 352 |
| DNA Ligase | 220110 | 10 | 88 | 176 |
| Adaptor Oligos | 2000094 | 20 | 176 | 352 |
| Total | | 50 | 440 | 880 |

- **b.** Add **50 μl** Adaptor Ligation Mix to **50 μl** sample. Pipette mix 15x (pipette set to 90 μl). Centrifuge briefly.
- **c.** Incubate in a thermal cycler with the following protocol.

| Lid Temperature | Reaction Volume | Run Time |
|-----------------|------------------------|----------------------|
| 30°C | 100 µl | 15 min |
| Step | Temperature | Time <i>hh:mm:ss</i> |
| 1 | 20°C | 00:15:00 |
| 2 | 4°C | Hold |

3.4 Post Ligation Cleanup – SPRIselect

- a. Vortex to resuspend SPRIselect Reagent. Add 80 μl SPRIselect Reagent (0.8X) to each sample. Pipette mix 15x (pipette set to 150 μl).
- **b.** Incubate **5 min** at **room temperature**.
- c. Place on the magnet-High until the solution clears.
- d. Remove the supernatant.
- e. Add 200 µl 80% ethanol to the pellet. Wait 30 sec.
- **f.** Remove the ethanol.
- g. Repeat steps e and f for a total of 2 washes.
- h. Centrifuge briefly. Place on the magnet-Low.
- **i.** Remove any remaining ethanol. Air dry for **2 min**. DO NOT exceed **2 min** as this will decrease elution efficiency.
- j. Remove from the magnet. Add **31 µl** Buffer EB. Pipette mix 15x.
- **k.** Incubate **2 min** at **room temperature**.
- **I.** Place on the magnet-Low until the solution clears.
- **m.** Transfer **30 µl** sample to a new tube strip.

3.5 Sample Index PCR

- **a.** Choose the appropriate sample index sets to ensure that no sample indices overlap in a multiplexed sequencing run. Record the 10x sample index name (PN-3000431 Dual Index Plate TT Set A well ID) used.
- **b.** Add **50 μl** Amp Mix (PN-2000047) to **30 μl** sample.
- **c.** Add **20 μl** of an individual Dual Index TT Set A to each sample and record the well ID used. Pipette mix 5x (pipette set to 90 μl). Centrifuge briefly.

| Lid Temperature | Reaction Volume | Run Time |
|-----------------|---|----------------------|
| 105°C | 100 μl | ~25-40 min |
| Step | Temperature | Time <i>hh:mm:ss</i> |
| 1 | 98°C | 00:00:45 |
| 2 | 98°C | 00:00:20 |
| 3 | 54°C | 00:00:30 |
| 4 | 72°C | 00:00:20 |
| 5 | Go to step 2, see below for # of cycles | |
| 6 | 72°C | 00:01:00 |
| 7 | 4°C | Hold |

d. Incubate in a thermal cycler with the following protocol.



The total cycles should be optimized based on 25% carry forward cDNA yield/input calculated during Post cDNA Amplification QC & Quantification (step 2.4).

Recommended Cycle Numbers

| cDNA Input | Total Cycles |
|----------------|--------------|
| 0.25-25 ng | 14-16 |
| 25-150 ng | 12-14 |
| 150-500 ng | 10-12 |
| 500-1,000 ng | 8-10 |
| 1,000-1,500 ng | 6-8 |
| >1500 ng | 5 |

STOP

e. Store at 4°C for up to 72 h or proceed to the next step.


3.6 Post Sample Index PCR Double Sided Size Selection – SPRIselect

- a. Vortex to resuspend the SPRIselect reagent. Add 60 μl SPRIselect Reagent (0.6X) to each sample. Pipette mix 15x (pipette set to 150 μl).
- b. Incubate 5 min at room temperature.
- **c.** Place the magnet**·High** until the solution clears. DO NOT discard supernatant.
- **d.** Transfer **150** µl supernatant to a new tube strip.
- e. Vortex to resuspend the SPRIselect reagent. Add 20 μl SPRIselect Reagent (0.8X) to each transferred supernatant. Pipette mix 15x (pipette set to 150 μl).
- f. Incubate 5 min at room temperature.
- g. Place the magnet-High until the solution clears.
- h. Remove 165 µl supernatant. DO NOT discard any beads.
- With the tube still in the magnet, add 200 μl 80% ethanol to the pellet.
 Wait 30 sec.
- j. Remove the ethanol.
- **k.** Repeat steps i and j for a total of 2 washes.
- **l.** Centrifuge briefly. Place on the magnet**·Low**. Remove remaining ethanol.
- m. Remove from the magnet. Add 36 µl Buffer EB. Pipette mix 15x.
- n. Incubate 2 min at room temperature.
- o. Place on the magnet-Low until the solution clears.
- **p.** Transfer **35** µl to a new tube strip.
- **q.** Store at **4°C** for up to **72 h** or at **-20°C** for **long-term** storage.

3.7 Post Library Construction QC

Run 1 μl sample at 1:10 dilution on an Agilent Bioanalyzer High Sensitivity chip.

Representative Trace

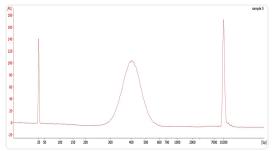

Determine the average fragment size from the Bioanalyzer trace. This will be used as the insert size for library quantification.

-ġ-

If peaks below 200 bp are more prominent, repeat step 3.6 Post Sample Index PCR Double Sided Size Selection - SPRIselect.

Alternate QC Method

Agilent TapeStation

LabChip

See Appendix for:

- Post Library Construction Quantification on page 83
- Agilent TapeStation Traces on page 84
- LabChip Traces on page 85



Step 4:


Sequencing

| Sequencing Libraries | 72 |
|--|----|
| Illumina Sequencer Compatibility | 73 |
| Sample Indices | 73 |
| 3' Gene Expression Library Sequencing Parameters | 74 |
| Library Loading | 75 |
| Library Pooling | 75 |
| Data Analysis and Visualization | 76 |

Sequencing Libraries

Chromium Single Cell 3' Gene Expression Dual Index libraries comprise standard Illumina paired-end constructs which begin with P5 and end with P7. These libraries include 16 bp 10x Barcodes at the start of TruSeq Read 1 while i7 and i5 sample index sequences are incorporated as the sample index read. TruSeq Read 1 and TruSeq Read 2 are standard Illumina sequencing primer sites used in paired-end sequencing of Single Cell 3' Gene Expression libraries. Sequencing these libraries produces a standard Illumina BCL data output folder.

Chromium Single Cell 3' Gene Expression Dual Index Library

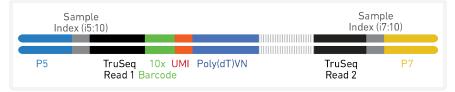

Illumina Sequencer Compatibility

The compatibility of the listed sequencers has been verified by 10x Genomics. Some variation in assay performance is expected based on sequencer choice. For more information about performance variation, visit the 10x Genomics Support website.

- MiSeq
- NextSeq 500/550
- NextSeq 1000/2000
- HiSeq 2500 (Rapid Run)
- HiSeq 3000/4000
- NovaSeq

Sample Indices

Each sample index in the Dual Index Kit TT Set A (PN-1000215) is a mix of one unique i7 and one unique i5 sample index. If multiple samples are pooled in a sequence lane, the sample index name (i.e. the Dual Index TT Set A plate well ID, SI-TT-) is needed in the sample sheet used for generating FASTQs. Samples utilizing the same sample index should not be pooled together, or run on the same flow cell lane, as this would not enable correct sample demultiplexing.

3' Gene Expression Library Sequencing Parameters

| Parameter | Description | |
|------------------|------------------------------------|--|
| Sequencing Depth | Minimum 20,000 read pairs per cell | |
| Sequencing Type | Paired-end, dual indexing | |
| Sequencing Read | Recommended Number of Cycles | |
| Read 1 | 28 cycles | |
| i7 Index | 10 cycles | |
| i5 Index | 10 cycles | |
| Read 2 | 90 cycles | |

Library Loading

Once quantified and normalized, the libraries should be denatured and diluted as recommended for Illumina sequencing platforms. Refer to Illumina documentation for denaturing and diluting libraries. Refer to the 10x Genomics Support website for more information.

3' Gene Expression libraries alone or in combination with other libraries

| Instrument | Loading Concentration (pM) | PhiX (%) |
|-------------------|----------------------------|----------|
| MiSeq | 11 | 1 |
| NextSeq 500/550 | 1.8 | 1 |
| NextSeq 1000/2000 | 650 | 1 |
| HiSeq 2500 (RR) | 11 | 1 |
| HiSeq 4000 | 240 | 1 |
| NovaSeq | 150*/300 | 1 |

Library Loading

Library Pooling

The 3' Gene Expression libraries maybe pooled for sequencing, taking into account the differences in cell number and per-cell read depth requirements between each library. Samples utilizing the same sample index should not be pooled together or run on the same flow cell lane, as this would not enable correct sample demultiplexing.

Library Pooling Example

Data Analysis and Visualization

Sequencing data may be analyzed using Cell Ranger or 10x Genomics Cloud Analysis and visualized using Loupe Browser. Key features for these tools are listed below. For detailed product-specific information, visit the 10x Genomics Support website.

Cell Ranger

Cell Ranger is a set of analysis pipelines that processes Chromium Single Gene Expression data to align reads, generate Feature Barcode matrices and perform clustering and gene expression analysis.

- Input: Base call (BCL) and FASTQ
- Output: BAM, MEX, CSV, HDF5, Web Summary, .cloupe/.loupe
- Operating System: Linux

Cloud Analysis

Cloud Analysis is currently only available for US customers.

Cloud Analysis allows users to run Cell Ranger analysis pipelines from a web browser while computation is handled in the cloud.

- Key features: scalable, highly secure, simple to set up and run
- Input: FASTQ
- Output: BAM, MEX, CSV, HDF5, Web Summary, .cloupe/.loupe

Loupe Browser

Loupe Browser is an interactive data visualization tool that requires no prior programming knowledge.

- Input: .cloupe
- Output: Data visualization, including t-SNE and UMAP projections, custom clusters, differentially expressed genes
- Operating System: MacOS, Windows



Troubleshooting



GEMs 78 Chromium X Series Errors 81

GEMs

Normal Step **Reagent Clogs & Wetting Failures** 1.4d After Chip M is removed from Chromium X and the wells are exposed N 0 М N 0 η. M All 16 recovery wells (rows 3A, 3B) (rows 3A, 3B) Recovery well G indicates a reagent clog. Recovery are similar in volume and opacity. well C and E indicate a wetting failure. Wells A, H, I & P contain 50% Glycerol Solution. Rest of the recovery wells have normal GEM generation.

1.4f Transfer GEMs



All liquid levels are similar in volume and opacity without air trapped in the pipette tips.




Pipette tips C and E indicate a wetting failure. Pipette tip C contains partially emulsified GEMs. Emulsion is absent in pipette tip E. Pipette tip G indicates a reagent clog.

2.1a After transfer of the GEMs + Recovery Agent



All liquid levels are similar in the aqueous sample volume (clear) and Recovery Agent/ Partitioning Oil (pink).

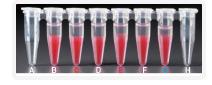

Tube G indicates a reagent clog has occurred. There is a decreased volume of aqueous layer (clear).

Tube C and E indicate a wetting failure has occurred. There is an abnormal volume of Recovery Agent/Partitioning Oil (pink).


Step

2.1 b After aspiration of Recovery Agent/ Partitioning Oil




All liquid volumes are similar in the aqueous sample volume (clear) and residual Recovery Agent/Partitioning Oil (pink).

Reagent Clogs & Wetting Failures

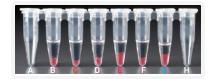

Tube G indicates a reagent clog has occurred. There is a decreased volume of aqueous layer (clear). There is also a greater residual volume of Recovery Agent/ Partitioning Oil (pink). Tube C and E indicate a wetting failure has occurred. There is an abnormal residual volume of Recovery Agent/Partitioning Oil (pink).

2.1 d After addition of Dynabeads Cleanup Mix

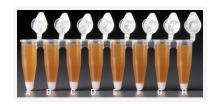


All liquid volumes are similar after addition of the Dynabeads Cleanup Mix.



Tube G indicates a reagent clog has occurred. There is an abnormal ratio of Dynabeads Cleanup Mix (brown) to Recovery Agent/Partitioning Oil (appears white). Tube C and E indicate a wetting failure has occurred. There is an abnormal ratio of Dynabeads Cleanup Mix (brown) to Recovery Agent/Partitioning Oil (appears white).

If a channel clogs or wetting failure occurs during GEM generation, it is recommended that the sample be remade. If any of the listed issues occur, take a picture and send it to support@10xgenomics.com for further assistance.

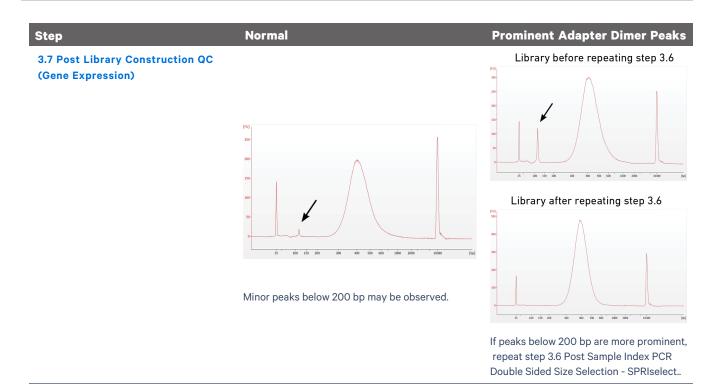

Chromium X Series Errors

The Chromium X touchscreen will guide the user through recoverable errors. If the error continues, or if the instrument has seen critical or intermediate errors, email support@10xgenomics.com with the displayed error code. Support will request a troubleshooting package. Upload pertinent logs to 10x Genomics by navigating to the Logs menu option on screen.

There are two types of errors:

Critical Errors – When the instrument has seen a critical error, the run will immediately abort. Do not proceed with any further runs. Contact support@10xgenomics.com with the error code.

- a. System Error
- **b.** Pressure Error
- c. Chip Error
- d. Run Error
- e. Temperature Error
- **f.** Software Error

User Recoverable Errors – Follow error handling instructions through the touchscreen and continue the run.

- a. Gasket Error
- **b.** Tray Error
- c. Chip Error
- d. Unsupported Chip Error
- e. Network Error
- **f.** Update Error

Consult the Chromium X Series (X/iX) User Guide (CG000396) for additional information and follow the Chromium X touchscreen prompts for execution. The Chromium X touchscreen will guide the user through recoverable errors.



Appendix

| Post Library Construction Quantification | 83 |
|--|----|
| Agilent TapeStation Traces | 84 |
| LabChip Traces | 85 |
| Oligonucleotide Sequences | 86 |

Post Library Construction Quantification

- a. Thaw KAPA Library Quantification Kit for Illumina Platforms.
- **b.** Dilute **2 μl** sample with deionized water to appropriate dilutions that fall within the linear detection range of the KAPA Library Quantification Kit for Illumina Platforms. (For more accurate quantification, make the dilution(s) in duplicate).
- **c.** Make enough Quantification Master Mix for the DNA dilutions per sample and the DNA Standards (plus 10% excess) using the guidance for 1 reaction volume below.

| Quantification Master Mix | 1Χ (μl) |
|-------------------------------|---------|
| SYBR Fast Master Mix + Primer | 12 |
| Water | 4 |
| Total | 16 |


- **d.** Dispense **16 μl** Quantification Master Mix for sample dilutions and DNA Standards into a 96 well PCR plate.
- **e.** Add **4 μl** sample dilutions and **4 μl** DNA Standards to appropriate wells. Centrifuge briefly.
- **f.** Incubate in a thermal cycler with the following protocol.

| Step | Temperature | Run Time |
|------|-------------------------------------|----------|
| 1 | 95°C | 00:03:00 |
| 2 | 95°C | 00:00:05 |
| 3 | 67°C Read Signal | 00:00:30 |
| 4 | Go to Step 2, 29X (Total 30 cycles) | |

g. Follow the manufacturer's recommendations for qPCR-based quantification. For library quantification for sequencer clustering, determine the concentration based on insert size derived from the Bioanalyzer/TapeStation trace.


Agilent TapeStation Traces

Agilent TapeStation High Sensitivity D5000 ScreenTape was used. Protocol steps correspond to the steps in this user guide.

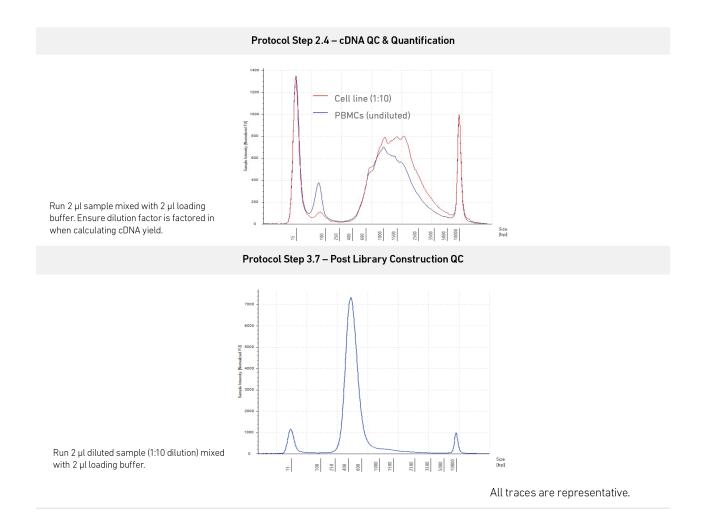

LabChip Traces

A DNA High Sensitivity Reagent Kit was used. Protocol steps correspond to the steps in this user guide.

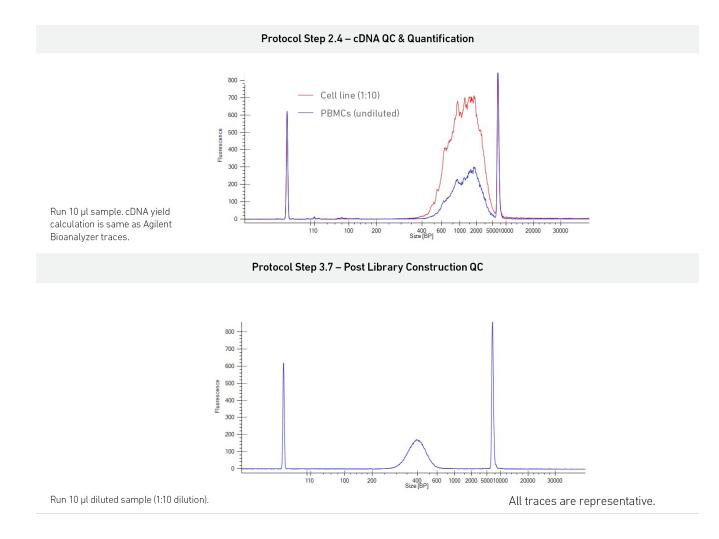

Alternate QC Method:

Qubit Fluorometer and Qubit dsDNA HS Assay Kit

Multiply the cDNA concentration reported via the Qubit Fluorometer by the elution volume (80 μ l) to obtain the total cDNA yield in ng. To determine the equivalent range using the Agilent 2100 Expert Software, select the region encompassing 35-10,000 bp.

Oligonucleotide Sequences

